

ED 147 192

SE 023 380

TITLE Effluent Monitoring Procedures: Metals Analyses.
Student Reference Manual.

INSTITUTION Environmental Protection Agency, Washington, D.C.
Office of Water Programs.

REPORT NO EPA-430-1-77-010

PUB DATE Sep 77

NOTE 319p.; For related documents, see SE 023 377-383

EDRS PRICE MF-\$0.83 HC-\$16.73 Plus Postage.

DESCRIPTORS Behavioral Objectives; Course Content; *Educational Programs; Environmental Education; *Instructional Materials; Laboratory Equipment; Laboratory Techniques; *Metals; *Pollution; Post Secondary Education; Skill Development; *Water Pollution Control

IDENTIFIERS *Waste Water Treatment

ABSTRACT

This is one of several short-term courses developed to assist in the training of waste water treatment plant operational personnel in the tests, measurements, and report preparation required for compliance with their NPDES Permits. The Student Reference Manual provides step-by-step procedures for laboratory application of equipment operating procedures for effluent monitoring. Each lesson outlines a specific objective, description of the analysis, and the applicability of the procedure. Parameters of this course include analyses of selected metals, including boron, copper, iron, mercury, sodium, and zinc. (CS)

 * Documents acquired by ERIC include many informal unpublished *
 * materials not available from other sources. ERIC makes every effort *
 * to obtain the best copy available. Nevertheless, items of marginal *
 * reproducibility are often encountered and this affects the quality *
 * of the microfiche and hardcopy reproductions ERIC makes available *
 * via the ERIC Document Reproduction Service (EDRS). EDRS is not *
 * responsible for the quality of the original document. Reproductions *
 * supplied by EDRS are the best that can be made from the original. *

EFFLUENT MONITORING PROCEDURES: METALS ANALYSES

FD147192

U.S. DEPARTMENT OF HEALTH
EDUCATION & WELFARE
NATIONAL INSTITUTE OF
EDUCATION

THIS DOCUMENT HAS BEEN REPRODUCED EXACTLY AS RECEIVED FROM THE PERSON OR ORGANIZATION ORIGINATING IT. POINTS OF VIEW OR OPINIONS STATED DO NOT NECESSARILY REPRESENT OFFICIAL NATIONAL INSTITUTE OF EDUCATION POSITION OR POLICY.

PERMISSION TO REPRODUCE THIS MATERIAL HAS BEEN GRANTED BY

Bernard Lukco

TO THE EDUCATIONAL RESOURCES INFORMATION CENTER (ERIC) AND USERS OF THE ERIC SYSTEM.



STUDENT REFERENCE MANUAL

U.S. ENVIRONMENTAL PROTECTION AGENCY
OFFICE OF WATER PROGRAM OPERATIONS

023 380

EFFLUENT MONITORING PROCEDURES: METALS ANALYSES

This course is designed for wastewater treatment plant technicians who will be responsible for performing the selected analyses.

Using methods approved by the U.S. Environmental Protection Agency for NPDES applications and reports, the student will perform selected metals analyses including: boron, copper, iron, mercury, sodium, and zinc.

Classroom instruction is limited to information about performing the selected analyses and reporting the results. Most of the time is given to laboratory experience for the trainee who uses detailed, stepwise procedures to analyze typical samples.

Those attending this course should be able to use laboratory glassware, and be able to do arithmetic calculations for formulas provided.

U.S. ENVIRONMENTAL PROTECTION AGENCY
Office of Water Program Operations
National Training and Operational Technology Center

DISCLAIMER

Reference to commercial products, trade names, or manufacturers is for purposes of example and illustration. Such references do not constitute endorsement by the Office of Water Program Operations, U. S. Environmental Protection Agency.

CONTENTS

<u>Title or Description</u>	<u>Outline Number</u>
Determination of Boron, Curcumin Method	1
Determination of Total Calcium (Volumetric Method)	2
Determination of Copper (Cu^{++}), Magnesium (Mg^{++}) Manganese (Mn^{++}), and Zinc (Zn^{++})	3
Determination of Lead by Atomic Absorption Using the Extraction Procedure	4
Determination of Mercury Using the Flameless Atomic Absorption (Cold Vapor) Technique	5
Determination of Potassium Using Flame Photometry	6
Determination of Sodium Using Flame Photometry	7

A PROTOTYPE FOR DEVELOPMENT OF
ROUTINE OPERATIONAL PROCEDURES

for the
DETERMINATION OF BORON, CURCUMIN, METHOD

as applied in
WASTEWATER TREATMENT FACILITIES
and in the
MONITORING OF EFFLUENT WASTEWATERS

Developed by the

National Training and Operational Technology Center,
Municipal Operations and Training Division
Office of Water Program Operations
U.S. Environmental Protection Agency

EFFLUENT MONITORING PROCEDURE: Determination of Boron, Curcumin Method

This operational procedure was developed by:

NAME : Audrey D. Kroner

ADDRESS EPA, OWPO, NTOTC, Cincinnati, Ohio 45268

POSITION Chemist-Instructor

EDUCATION AND TECHNICAL BACKGROUND

B.A. - Edgecliff College

1 year Industrial Research Chemist

8 years Secondary School Chemistry Instructor

4 years DHEW-DI, Water Quality Program Chemist

7-1/2 years DI-EPA, Chemist-Instructor

EFFLUENT MONITORING PROCEDURE: Determination of Boron, Curcumin Method

1. Objective:

To determine the mg/l concentration of boron in an unfiltered sample.

2. Brief Description of Analysis:

A sample of water containing boron is acidified with hydrochloric acid and evaporated in the presence of curcumin. The reaction forms a red-colored product called rosocyanine. The rosocyanine is taken up in ethyl alcohol and the red color in the solution is compared to that of standards in a spectrophotometer.

Calcium, magnesium and other cations interfere with the spectrophotometric measurements. These form salts which will not dissolve in ethyl alcohol, thereby contributing turbidity to the solution of rosocyanine. Such cationic interferences are removed by filtering the samples after rosocyanine is formed and dissolved in the ethyl alcohol.

3. Applicability of this Procedure:

a. Range of Concentration:

0.10 to 1.0 mg/liter boron (The range may be extended for samples by dilution.)

b. Pretreatment of Samples:

The Federal Register Guidelines do not specify any pretreatment (such as digestion) for determining total boron. A 0.45 micron filtration of samples is specified for determining dissolved boron. This procedure does not include directions for the filtration process.


c. Treatment of Interferences in Samples:

The Source of Procedure* cites turbidity caused by hardness cations and nitrate nitrogen concentrations in excess of 20 mg/liter as interferences. This procedure includes directions for removing turbidity in samples but does not include determination and removal of nitrate nitrogen. Consult the Source of Procedure* if the presence of excessive nitrate nitrogen is suspected in samples.

*Source of Procedure: Standard Methods for the Examination of Water and Wastewater, 14th ed., 1976, APHA, Washington, DC, p. 287.

EFFLUENT MONITORING PROCEDURE: Determination of Boron, Curcumin Method

FLOW SHEET

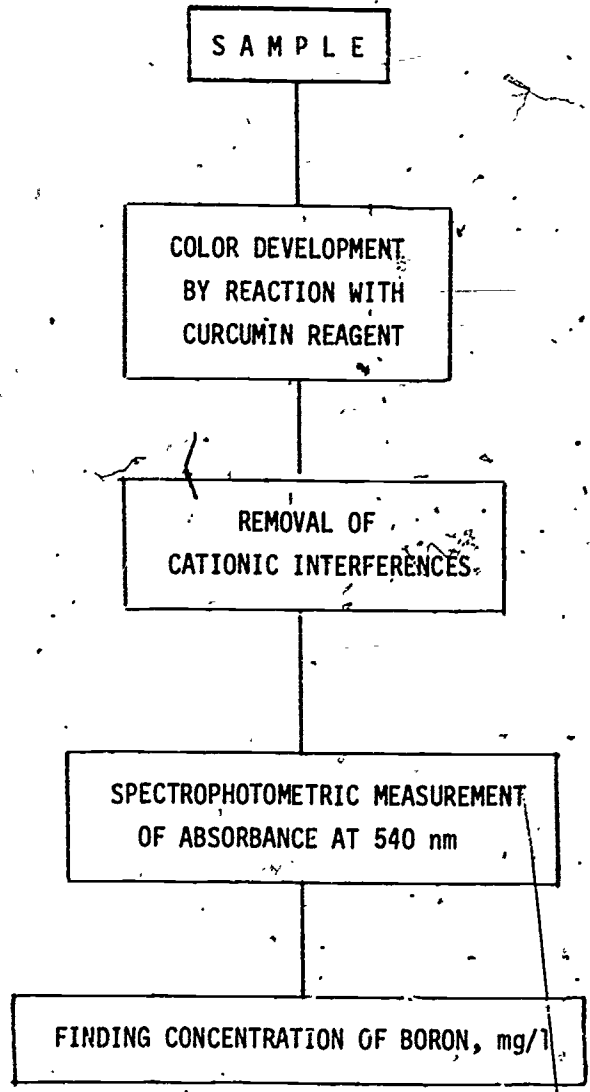

EFFLUENT MONITORING PROCEDURE: Determination of Boron, Curcumin Method

Equipment and Supply Requirements

A. Capital Equipment

1. Balance, analytical with a 0.1 milligram sensitivity.
2. Fume hood or equivalent.
3. Refrigerator capable of maintaining a 4°C temperature.
4. Spectrophotometer for use at 540 nm, with a minimum path length of 1 cm.
5. Water bath, constant temperature to maintain 55 ± 2°C. Size depends on number of samples to be processed along with 5 standards in 100 ml evaporating dishes. Each requires about 5 square inches of space.
6. Water still, metal or a mixed bed ion exchange resin cartridge to prepare boron-free distilled water.

B. Reusable Supplies

NOTE: Glassware made of Corning alkali-resistant glass and storage vessels made of ordinary soft glass have been satisfactory as "boron-free" glassware.

1. 1 Beaker, 50 ml
2. 1 Bottle, 500 ml, boron-free* glass with stopper.
3. 1 Bottle, 1000 ml, polyethylene with screw cap or boron-free* glass with stopper.
4. 1 Bottle, 5000 ml, polyethylene with screw cap (for distilled water).
5. 1 Cylinder, graduated, 250 ml.
6. Dishes, evaporating, porcelain, 100 ml, 5 plus 1 for each sample to be run.
7. 1 Dropper, 1 to 2 ml with bulb.
8. Flasks, volumetric, 25 ml, with stoppers. 5 plus 1 for each sample.
9. Flasks, volumetric, 100 ml, with stoppers. 6 plus 1 for each sample to be run.
10. 1 Flask, volumetric, 1000 ml, with stopper.
11. Funnels, fluted, 600, 40 mm diameter, 50 mm stem length. 1 for each sample.
12. 1 Glass rod, short.
13. 1 Mortar and pestle, 70 ml, porcelain (needed only if curcumin is not in finely powdered form).
14. 1 Pipet, measuring, Mohr, 10 ml.
15. Pipets, volumetric, 1.0 ml, 5 plus 1 for each sample.
16. 1 Pipet, volumetric, 2.0 ml.
17. 1 Pipet, volumetric, 4.0 ml.
18. 1 Pipet, volumetric, 5.0 ml.
19. 1 Pipet, volumetric, 8.0 ml. Alternatively, use a 10 ml measuring (Mohr) pipet.
20. 1 Pipet, volumetric, 10.0 ml.
21. 1 Pipet, volumetric, 100.0 ml.
22. 1 Stirring rod (policeman), polyethylene, about 6 inch length.
23. 1 Propipet or pipet bulb.
24. 1 Support, ring stand with small ring (for filtration)
25. 1 Thermometer, degrees C. (To check water bath temperature.)
26. 1 Triangle (for filter funnel).
27. 2 Wash bottles, polyethylene squeeze type.
28. 1 Laboratory apron.
29. 1 Pair safety glasses.

EFFLUENT MONITORING PROCEDURE: Determination of Boron, Curcumin Method

C. Consumable Supplies*

1. 5 Weighing boats (disposable plastic is suitable).
2. Graph paper: 8 1/2 inch by 11 inch with 10 major divisions along the 11 inch side is suitable.
3. Wax pencil.
4. Tissues, lint-free for wiping spectrophotometer cells.
5. Water, distilled, boron-free 2.5 - 3 liters.
6. Boric acid (H_3BO_3), ACS Reagent, anhydrous powder (1 lb unit).
7. Curcumin [$(2-CH_3OC_6H_3-1-OH-4-CH:GHCO)_2CH_2$] Eastman No. 1179 or equivalent (25 g unit).
8. 95% Ethyl alcohol (CH_3CH_2OH), Denatured is acceptable (1 gal. or 3 kg unit).
9. Hydrochloric acid (HCl), ACS Reagent, (5 pt. or 7 lb. unit).
10. Oxalic acid ($H_2C_2O_4$), ACS Reagent (1 lb. unit).
11. 1 Box filter paper, Whatman No. 30 or equivalent with diameter to fit filter funnel. (A diameter of 7 cm will fit the filter funnel described in B.)

*NOTE: Use polyethylene or boron-free glass bottles for storage of reagents after preparation.

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOAL 'SPECIFICATIONS	TRAINING GUIDE NOTES
<p>A. Preparing to Test the Sample.</p>	<ol style="list-style-type: none"> 1. Assemble all equipment to be used. 2. Check the water bath for maintenance of the $55 \pm 2^\circ\text{C}$ temperature required for the test. 3. If necessary, clean all labware to be used for the test. 4. Prepare the reagents for the test. 5. Record the identification information for the sample(s). 6. Record the date and time you are beginning the analysis. 7. Begin the analysis with Procedure E, Preparation of Standards: 	<ol style="list-style-type: none"> 1a. Equipment list is on pp. 6 and 7. 2a. Use Procedure B, Checking the Water Bath. 3a. Use Procedure C, Cleaning Labware. 3b. It is most convenient to clean labware immediately after performing the test so it is dry and ready for the next use. 4a. Use Procedure D, Reagent Preparation. 5a. The sample(s) should be at hand before continuing with the test. 5b. Use a laboratory notebook with space for information similar to the "Example Laboratory Data Sheet." 5c. In one column for each sample record: "Identification Code," "Type" (grab or composite), "Date and Time Collected," and the name of the "Sample Collector." 6a. Use the "Example Laboratory Data Sheet." 6b. In the column(s) of information you recorded for the sample(s). 	<p>IX Sheet 1 (p. 33)</p> <p>IX.A. (p. 33)</p> <p>IX.A.6. (p. 33)</p>

EFFLUENT MONITORING PROCEDURE: Determination of Boron, Curcumin Method

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>B. Checking the Water Bath</p>	<ol style="list-style-type: none"> 1. Fill the water bath to a level where the evaporating dishes for the test can float. 2. Place a Centigrade thermometer into the water bath. 3. Set the temperature regulator to 55°C. 4. Turn the bath heating element on. 5. Allow the bath to heat to 55°C. 6. At least ten minutes after the bath reaches 55°C, check the temperature of the water bath. 7. Leave the bath turned on. 	<ol style="list-style-type: none"> 1a. The water bath should be in a fume hood. 2a. The thermometer should be reliably accurate. 5a. You could start Procedure D, Reagent Preparation, during this time. 6a. A longer time is all right. 6b. The temperature should be between 53 and 57°C. 6c. If the temperature is below 53°C, adjust the temperature regulator upward and repeat steps 5 and 6. 6d. If the temperature is above 57°C, adjust the temperature regulator downward and repeat steps 5 and 6. 	
<p>14</p>			<p>15</p>

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>C. Cleaning Labware</p>	<ol style="list-style-type: none"> 1. All bottles, evaporating dishes, flasks, graduates, pipets, stirring rods, filter funnels and spectrophotometry cells for this procedure should be cleaned using the following steps. 2. Rinse each item thoroughly with tap water. 3. Rinse each item with warm 1:1 hydrochloric acid (HCl) solution. 4. Rinse each item thoroughly with tap water. 5. Rinse each item with seven individual rinses of boron-free distilled water. 6. Place items in a clean area to drain dry on clean towels. 	<ol style="list-style-type: none"> 1a. New labware should be cleaned before using it for this test. 3a. In a fume hood. 3b. Wear rubber gloves. 3c. CAUTION: Hydrochloric acid and its fumes are hazardous. 3d. The steps to prepare 1:1 hydrochloric acid solution are in D., Reagent Preparation, Reagent 2. 3e. 1:1 nitric acid solution could be used instead. 5a. Several rinses with smaller volumes of water are more effective than one or two large-volume rinsings. 	
<p>16</p>		<p>17</p>	

EFFLUENT MONITORING PROCEDURE: Determination of Boron, Curcumin Method

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>D. Reagent Preparation</p> <p>1. Boron-free distilled water..</p> <p>2. 1:1 Hydrochloric acid solution</p>	<p>1. Prepare boron-free distilled water.</p> <p>1. Measure 250 ml boron-free distilled water.</p>	<p>1a. Use a metal still or a mixed bed ion exchange resin</p> <p>1b. About 5 liters are required for reagents, etc., plus about 1 ml for each standard and sample to be analyzed.</p> <p>1c. Store the water in polyethylene bottles or boron-free glass bottles.</p> <p>1a. Use a graduate to measure the water.</p>	
<p>18</p>	<p>2. Pour the water into a clean, boron-free 500 ml storage bottle.</p> <p>3. Measure 250 ml concentrated hydrochloric acid (HCl).</p> <p>4. Slowly pour the acid into the bottle of water, along the inside wall.</p> <p>5. Swirl the mixture.</p>	<p>2a. Use a polyethylene, screw-capped bottle or a boron-free glass bottle with stopper.</p> <p>2b. The bottle should be cleaned with 1:1 hydrochloric acid solution. To do this, follow these steps using 0.1 of the volumes given, then use Procedure C, Cleaning Labware.</p> <p>3a. Use a graduate.</p> <p>3b. In a fume hood.</p> <p>3c. CAUTION; Hydrochloric acid causes severe burns. Also, the fumes are hazardous and can cause poisoning if inhaled.</p> <p>3d. 1:1 nitric acid could be prepared using these steps if its use for cleaning labware is preferred.</p> <p>4a. To avoid spattering of the mixture.</p> <p>5a. To thoroughly mix the contents.</p>	<p>19</p>



OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
D. Reagent Preparation (continued)	6. Label the bottle.	6a. This is 1:1 hydrochloric acid. Also, write the date and your name. 6b. This solution is used to clean glassware used in the test procedure.	
3. Curcumin reagent	1. Add about 35 ml of 95% ethyl alcohol ($\text{CH}_3\text{CH}_2\text{OH}$) to a 100 ml volumetric flask. 2. Finely grind the curcumin [$(2\text{-CH}_3\text{OC}_6\text{H}_3\text{-1-OH-4-CH:CHCO})_2\text{CH}_2$] if it is not in this form. 3. Weigh out 0.040 g curcumin 4. Carefully transfer the curcumin to the 100 ml volumetric flask. 5. Swirl the mixture. 6. Weigh out 5.0 g oxalic acid ($\text{H}_2\text{C}_2\text{O}_4$). 7. Carefully transfer the oxalic acid to the same flask. 8. Swirl the mixture.	1a. You need 4.0 ml of this reagent for each standard and sample to be analyzed. 3a. Use a weighing boat. 3b. Use an analytical balance. 4a. Use 95% ethyl alcohol in a squeeze-type wash bottle. 6a. This oxalic acid intensifies the color of the reaction product and also prevents the formation of interfering colors produced by some metals, if present. 7a. Again, use 95% ethyl alcohol.	

EFFLUENT MONITORING PROCEDURE: Determination of Boron, Curcumin Method

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>D. Reagent Preparation (continued)</p> <p>4. 95% Ethyl alcohol</p> <p>5. Stock boron solution</p>	<p>9. Measure 4.2 ml concentrated hydrochloric acid (HCl).</p> <p>10. Add the acid to the same flask.</p> <p>11. Swirl the mixture.</p> <p>12. Dilute the mixture up to the 100.0 ml mark with 95% ethyl alcohol.</p> <p>13. Stopper the flask.</p> <p>14. Invert the flask several times.</p> <p>15. Label the flask.</p> <p>16. Store in a refrigerator.</p> <p>1. This needs no preparation but have a supply at hand to perform the procedure.</p> <p>1. Add about 500 ml boron-free distilled water to a 1000 ml volumetric flask.</p>	<p>9a. Use a graduated pipet.</p> <p>14a. To mix the contents.</p> <p>15a. This is curcumin reagent. Also write the date and your name.</p> <p>16a. This reagent will be stable up to a week when stored in a cool, dark place.</p> <p>1a. You need 25 ml for each standard and sample.</p> <p>1b. Some of the alcohol should be in a squeeze-type wash bottle.</p>	



OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
D. Reagent Preparation (continued)	<ol style="list-style-type: none"> 2. Weigh out 0.5716 g anhydrous boric acid (H_3BO_3). 3. Carefully transfer the boric acid to the 1000 ml flask. 4. Swirl the mixture. 5. Dilute the mixture up to the 1000.0 ml mark with boron-free distilled water. 6. Pour this stock solution into a 1000 ml boron-free reagent bottle. 7. Label the bottle. 8. Store in a refrigerator. 	<ol style="list-style-type: none"> 2a. Use a weighing boat. 2b. Use an analytical balance. 2c. Do not dry the boric acid in a 103°C oven; it loses weight. Use ACS reagent grade boric acid and keep the chemical bottle tightly stoppered. 6a. Use a polyethylene, screw-capped bottle or a boron-free, glass bottle with a stopper. 7a. This is the stock boron solution. 1.00 ml = 100 µg B. Also, write the date and your name on the label. 1a. The standard solution should be prepared fresh each day samples are to be analyzed. 1b. Although small quantities of this solution are required, make one liter to increase the accuracy of its concentration. 	
6. Standard boron solution	<ol style="list-style-type: none"> 1. Add about 500 ml boron-free distilled water to a 1000 ml volumetric flask. 		

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>D. Reagent Preparation (continued)</p>	<ol style="list-style-type: none"> 2. Measure 100.0 ml of the stock boron solution. 3. Deliver the stock solution to the volumetric flask. 4. Swirl the flask 5. Dilute the mixture up to the 1000.0 ml mark with boron-free distilled water. 6. Label the flask. 	<ol style="list-style-type: none"> 2a. The stock solution must be at room temperature before this measurement is done. 2b. Use a volumetric pipet. 6a. This is the standard boron solution. 1.00 ml = 10.0 µg B. Also, write the date and your name on the label. 	
<p>E. Preparation of Standards</p>	<ol style="list-style-type: none"> 1. If you are establishing a calibration curve, prepare standards using the following steps. If a calibration curve has already been established, prepare a blank and a check standard beginning below at Step 11. 	<ol style="list-style-type: none"> 1a. You should prepare a calibration curve each day you do this test. If several samples are to be analyzed, a blank and a standard to check the curve should be run with each group of samples run after the curve has been prepared. 	

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES																		
<p>E. Preparation of Standards (continued)</p>	<p>2. Mark five 100 ml volumetric flasks with the concentrations of boron, mg/liter which are shown in the second column in Table 2d.</p> <p>3. Add about 40 ml boron-free distilled water to each flask.</p> <p>4. Do not put any standard solution into the flask marked "0.00".</p> <p>5. Into each of the remaining marked flasks, pipet the corresponding volume of standard boron solution as shown in the third column in Table 2d.</p> <p>6. Do not add any more water to the flask marked "0.00".</p>	<p>2a. Flasks must be clean. 2b. Use a wax pencil. 2c. These concentrations are recorded on the "Example Laboratory Data Sheet". 2d. Table 2d:</p> <table border="1" data-bbox="968 371 1478 697"> <thead> <tr> <th>Flask</th> <th>Boron, mg per liter</th> <th>Boron, ml Standard</th> </tr> </thead> <tbody> <tr> <td>1</td> <td>0.00</td> <td>None</td> </tr> <tr> <td>2</td> <td>0.20</td> <td>2</td> </tr> <tr> <td>3</td> <td>0.50</td> <td>5</td> </tr> <tr> <td>4</td> <td>0.80</td> <td>8</td> </tr> <tr> <td>5</td> <td>1.00</td> <td>10</td> </tr> </tbody> </table> <p>4a. This flask contains the blank.</p> <p>5a. The concentration of the standard boron solution is: 1:00 ml = 10.0 μg B. 5b. Use volumetric pipets.</p> <p>6a. A volumetric measurement is not necessary for the blank at this time.</p>	Flask	Boron, mg per liter	Boron, ml Standard	1	0.00	None	2	0.20	2	3	0.50	5	4	0.80	8	5	1.00	10	<p>IX.E.2. (p. 33)</p> <p>29</p>
Flask	Boron, mg per liter	Boron, ml Standard																			
1	0.00	None																			
2	0.20	2																			
3	0.50	5																			
4	0.80	8																			
5	1.00	10																			

EFFLUENT MONITORING PROCEDURE: Determination of Boron, Curcumin Method

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>E. Preparation of Standards. (continued)</p>	<p>7. To each of the other 4 flasks, carefully add boron-free distilled water up to the 100.0 ml mark.</p> <p>8. Put a stopper in each flask.</p> <p>9. Gently invert each flask several times.</p> <p>10. Go to the next Procedure, E. Color Development.</p> <p>11. If a calibration curve has been established, prepare a blank and a check standard using the following steps.</p> <p>12. Mark one 100 ml volumetric flask "0.00" mg/liter B.</p> <p>13. Add about 40 ml boron-free distilled water.</p> <p>14. Mark a second 100 ml volumetric flask "0.50" mg/liter B.</p>	<p>7a. Use a squeeze bottle for the final addition of water.</p> <p>9a. To ensure thorough mixing.</p> <p>11a. The blank is used to zero the spectrophotometer.</p> <p>11b. The check standard is compared to the standard curve to check reproducibility.</p> <p>12a. Use a wax pencil.</p> <p>13a. This flask contains the blank.</p> <p>13b. A volumetric measurement is not necessary for the blank at this time.</p> <p>14a. This concentration is recorded on the "Example Laboratory Data Sheet."</p>	<p>IX.E.14 (p. 33)</p>

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>E. Preparation of Standards (continued)</p>	<p>15. Into this flask, pipet 5.0 ml of the standard boron solution.</p> <p>16. Add boron-free distilled water up to the 100.0 ml mark on the flask.</p> <p>17. Put a stopper in the flask.</p> <p>18. Gently invert the flask several times.</p> <p>19. Go to the next Procedure, F. Color Development.</p>	<p>15a. The concentration of the standard boron solution is: 1.00 ml = 10.0 μg B.</p> <p>15b. Use a volumetric pipet.</p> <p>16a. Use a squeeze bottle for the final addition of water.</p> <p>18a. To ensure thorough mixing.</p>	
<p>F. Color Development of Standards and Sample</p>	<p>1. If you are establishing a calibration curve for the day, use the following steps. If you are using a blank and check standard, begin below at Step 5.</p> <p>2. Mark five 100 ml evaporating dishes with the concentrations of boron, mg/liter which are shown in Table 2e.</p>	<p>2a. Each evaporating dish should be clean, dry, and free from scratches and scoring.</p> <p>2b. The dishes must be identical in shape, size, and composition to ensure equal evaporation time.</p> <p>2c. Use wax pencil markings on the outside near the top.</p> <p>2d. These concentrations are recorded on the "Example Laboratory Data Sheet."</p>	

(continued)



EFFLUENT MONITORING PROCEDURE: Determination of Boron, Curcumin Method

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES																		
<p>F. Color Development of Standards and Sample (continued)</p>	<p>3. Pipet 1.0 ml of the 0.00 mg/l B (blank) solution and of each of the standards into the correspondingly marked evaporating dish.</p> <p>4. Skip Steps 5, 6 and 7 and continue at Step 8.</p> <p>5. Mark two 100 ml evaporating dishes with the concentrations of boron, mg/liter which are shown in Table 5e.</p>	<p>2e. Table 2e:</p> <table border="1" data-bbox="1207 333 1692 601"> <thead> <tr> <th>Dish</th> <th>Boron, mg per liter</th> </tr> </thead> <tbody> <tr> <td>1</td> <td>0.00</td> </tr> <tr> <td>2</td> <td>0.20</td> </tr> <tr> <td>3</td> <td>0.50</td> </tr> <tr> <td>4</td> <td>0.80</td> </tr> <tr> <td>5</td> <td>1.00</td> </tr> </tbody> </table> <p>3a. Use a clean, dry, volumetric pipet for each solution.</p> <p>3b. The standards were prepared in 100 ml flasks using E, Preparation of Standards.</p> <p>4a. These steps are used for a blank and check standard.</p> <p>5a. Each evaporating dish should be clean, dry, and free from scratches and scoring.</p> <p>5b. The dishes must be identical in shape, size, and composition to ensure equal evaporation time.</p> <p>5c. Use wax pencil markings on the outside near the top.</p> <p>5d. These concentrations are recorded on the "Example Laboratory Data Sheet."</p> <p>5e.</p> <table border="1" data-bbox="1207 1202 1735 1335"> <thead> <tr> <th>Dish</th> <th>Boron, mg per liter</th> </tr> </thead> <tbody> <tr> <td>1</td> <td>0.00</td> </tr> <tr> <td>2</td> <td>0.50</td> </tr> </tbody> </table>	Dish	Boron, mg per liter	1	0.00	2	0.20	3	0.50	4	0.80	5	1.00	Dish	Boron, mg per liter	1	0.00	2	0.50	
Dish	Boron, mg per liter																				
1	0.00																				
2	0.20																				
3	0.50																				
4	0.80																				
5	1.00																				
Dish	Boron, mg per liter																				
1	0.00																				
2	0.50																				

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>F. Color Development of Standards and Sample (continued)</p>	<p>6. Pipet 1.0 ml of the 0.00 mg/l B (blank) into the dish marked "0.00".</p> <p>7. Pipet 1.0 ml of the 0.50 mg per liter standard into the dish marked "0.50".</p> <p>8. Record the sample identification code for the sample.</p> <p>9. Mark the sample identification code on a 100 ml evaporating dish.</p> <p>10. Measure 1.0 ml of the sample (or of diluted sample) into the marked evaporating dish.</p>	<p>6a. Use a clean, dry, volumetric pipet.</p> <p>7a. Use a clean, dry, volumetric pipet.</p> <p>8a. On the "Example Laboratory Data Sheet" to the left of the blank where "Absorbance" will be recorded for the sample.</p> <p>9a. The dish should be clean, dry, and free from scratches and scoring.</p> <p>9b. The dish should be identical in shape, size, and composition to the dishes used for the standards.</p> <p>9c. Make a wax pencil mark on the outside near the top.</p> <p>10a. For samples containing 0.10 to 1.00 mg per liter boron, use undiluted sample.</p> <p>10b. For samples containing more than 1.00 mg per liter boron, make an appropriate dilution with boron-free distilled water so that a 1.00 ml portion contains approximately 0.50 mg boron (50 mg per liter).</p> <p>10c. If the boron concentration of the sample is unknown, you might use 1.00 ml of undiluted sample and also 1.00 ml of a 50% dilution (1:1) of the sample.</p>	<p>IX.F.8. (p. 33)</p>



EFFLUENT MONITORING PROCEDURE: Determination of Boron, Curcumin Method

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>F. Color Development of Standards and Sample (continued)</p>	<p>11. Repeat Steps 8, 9, and 10 for each sample to be analyzed.</p> <p>12. Add 4.0 ml curcumin reagent to each evaporating dish.</p> <p>13. Gently swirl each dish.</p> <p>14. Float the dishes on a water bath set at $55 \pm 20C$ and let them remain for 80 minutes to evaporate to dryness.</p> <p>15. Record the time after 80 minutes have elapsed.</p>	<p>11a. The number of samples that can be analyzed along with the standards depends on the size of the water bath.</p> <p>12a. Use a volumetric pipet for better control of this alcohol solution.</p> <p>12b. If you are establishing a standard curve, there are 5 dishes of standards plus the dish(es) of sample(s).</p> <p>12c. If a standard curve is established, there are 2 dishes of standards plus the dish(es) of sample(s).</p> <p>13a. To thoroughly mix contents.</p> <p>14a. The water bath should have been checked for maintenance of this temperature and turned on as described in B, Checking the Water Bath.</p> <p>14b. The temperature is critical to the test. Higher temperatures cause loss of color.</p> <p>14c. A red-colored product, rosocyanine, is formed.</p> <p>14d. The odor of hydrochloric acid (HCl) is gone when evaporation is complete.</p> <p>14e. Keep drying time constant for standards and samples.</p> <p>14f. The length of drying time critically affects the intensity of color.</p> <p>15a. This is the "Time Evaporation Ended" on the "Example Laboratory Data Sheet".</p> <p>15b. Spectrophotometric readings must be done within the next hour.</p>	<p>IX.F.15. (p. 33)</p>

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
F. Color Development of Standards and Sample (continued)	16. Turn the water bath heating element off.	16a. Unless more samples are to be determined.	
	17. Remove each dish from the water bath.	17a. Be careful that drops of water from the bottom of one dish do not drip into other dishes still in the water bath.	
	18. Carefully wipe the moisture from the bottom of each dish.		
	19. Allow the dishes to cool to room temperature.	19a. Five to 10 minutes.	
	20. Turn on the spectrophotometer.	20a. The instrument can warm up during the next steps.	
	21. Mark 25 ml volumetric flask with the concentrations of the standards you are running.	21a. If you are establishing a standard curve, mark 5 flasks as 0.00, 0.20, 0.50, 0.80, and 1.00 mg per liter, respectively. 21b. If you are running a blank and check standard, mark 2 flasks as 0.00 and 0.50 mg per liter, respectively.	
	22. Mark 25 ml volumetric flask(s) with the identification code(s) of the sample(s).	22a. If any flasks are not dry, rinse them with 95% ethyl alcohol.	
23. Add about 7 ml of 95% ethyl alcohol ($\text{CH}_3\text{CH}_2\text{OH}$) to the dish containing sample.	23a. Use a measuring pipet and very slowly direct the alcohol along the walls of the dish so any spattered solids are washed down to the bottom. 23b. Transfer samples first, then the standards.		

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>F. Color Development of Standards and Sample (continued)</p>	<p>24. Use a clean, dry polyethylene rod to gently stir the mixture until all the red-colored product is dissolved.</p> <p>25. Use a squeeze bottle of 95% ethyl alcohol to rinse the mixture from the tip of the rod into the dish.</p> <p>26. Using a dropper, transfer the solution from the dish into the corresponding 25 ml volumetric flask.</p> <p>27. Add a few ml of 95% ethyl alcohol to the dish.</p> <p>28. Use the same dropper to transfer this rinse to the corresponding flask.</p> <p>29. Repeat steps 27 and 28 two more times.</p> <p>30. Remove the bulb from the dropper and rinse 95% ethyl alcohol from the top through the tip and into the flask.</p>	<p>24a. Keep the mixture in the bottom part of the dish.</p> <p>25a. Keep in mind that 25 ml is the final volume allowed in the following steps.</p> <p>26a. Be careful not to scratch the surface of the evaporating dish with the tip of the glass dropper.</p> <p>26b. Be careful not to lose any of the solution during the transfer.</p> <p>26c. A very small funnel can be used for this transfer.</p> <p>27a. Use a squeeze-type plastic wash bottle for the alcohol.</p> <p>29a. Add just a few ml each time to keep within the 25 ml limit of the receiving volumetric flask.</p> <p>29b. The absence of yellow color from all areas of the dish indicates all original solution has been transferred.</p>	

OPERATING, PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
F. Color Development of Standards and Sample (continued)	<p>31. Rinse the outside of the tip of the dropper into the flask.</p> <p>32. Replace the bulb on the dropper.</p> <p>33. Add a few more ml of 95% ethyl alcohol to the dish:</p> <p>34. With the dropper, use this alcohol to bring the volume in the flask to the 25.0 ml mark.</p> <p>35. Put a stopper in the flask.</p> <p>36. Gently invert the flask several times.</p> <p>37. Repeat steps 23 through 36 for any other samples, then for the standards in evaporating dishes.</p> <p>38. Obtain all spectrophotometric readings within one hour of the time evaporation ended by using H, Spectrophotometric Measurements.</p>	<p>36a. To thoroughly mix the contents.</p> <p>36b. IMPORTANT: Any turbidity in the sample solution(s) at this point interferes with spectrophotometric readings. Use Procedure G to filter sample(s) now.</p> <p>37a. Check that each is transferred to the flask marked with the corresponding sample identification code or concentration.</p> <p>38a. The time evaporation ended is recorded on the "Example Laboratory Data Sheet".</p>	

EFFLUENT MONITORING PROCEDURE: Determination of Boron, Curcumin Method

OPERATING PROCEDURES	STEP, SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>G. Removal of Cationic Interferences from the Sample(s)</p>	<ol style="list-style-type: none"> 1. Mark a small receiving vessel with the sample identification code. 2. Place a small funnel into a clay triangle on a ring stand support. 3. Flute a piece of filter paper to fit into the funnel. 4. Place the paper in the funnel. 5. DO NOT "wet" the filter paper. 6. Using a small glass rod, transfer the sample from the flask onto the filter paper. 7. DO NOT rinse the original flask and DO NOT add alcohol to the filtered solution. 8. Remove the receiving vessel from under the funnel. 	<ol style="list-style-type: none"> 1a. Samples should be filtered to remove turbidity. 1b. One of a set of matched spectrophotometer tubes or a very small beaker would be suitable receiver. 2a. A funnel with a 40 mm diameter and a 50 mm stem works well with this volume. 3a. Use Whatman No. 30 or equivalent filter paper. 3b. A diameter of 7 cm fits the above funnel. 5a. You could change the concentration of the sample. 6a. While sample is being filtered, continue with transfers of other samples, then the standards from evaporation dishes to flasks at Step F.23 7a. Adding any alcohol will change the concentration of the sample. You only need about 15 ml of filtered sample to get spectrophotometric readings. 	

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>G. Removal of Cationic Interferences from the Sample(s) (continued)</p>	<p>9. Repeat Steps 1 through 8 for each sample being analyzed.</p> <p>10. Obtain spectrophotometric readings within one hour of the time evaporation ended by using H, Spectrophotometric Measurements.</p>	<p>9a. Rinse out the funnel and dry it between uses so no liquid is added to the next sample.</p> <p>10a. The time evaporation ended is recorded on the "Example Laboratory Data Sheet".</p>	
<p>H. Spectrophotometric Measurements</p>	<p>1. Check that the instrument is warmed up and ready to use.</p> <p>2. Set the wavelength at 540 nm.</p> <p>3. Set the instrument at infinite absorbance.</p> <p>4. Use the 0.00 mg per liter boron standard to set the instrument at zero absorbance.</p> <p>5. Remove the 0.00 mg per liter boron standard.</p> <p>6. Check that the scale reading returns to infinite absorbance.</p>	<p>1a. Allow a warm-up period of approximately 20 minutes (5 minutes minimum).</p> <p>1b. There is an EMP on "Use of a Spectrophotometer".</p> <p>4a. Use alcohol for any rinsings of the spectrophotometer tube.</p> <p>4b. Wipe the outside of the spectrophotometer tube to remove moisture and fingerprints before inserting the tube into the instrument.</p> <p>4c. Use this same spectrophotometer for standards and samples processed in the same-size evaporating dish as was used to process this blank.</p> <p>6a. If the reading is not at infinite absorbance, set the control knob until that reading occurs, reinsert the 0.00 mg per liter standard and reset the instrument at zero absorbance.</p> <p>6b. Then repeat steps 5 and 6.</p>	



EFFLUENT MONITORING PROCEDURE: Determination of Boron, Curcumin Method

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>H. Spectrophotometric Measurements (continued)</p>	<p>7. Record the time.</p> <p>8. Measure and record the absorbance of each standard.</p> <p>9. Measure and record the absorbance of each sample.</p> <p>10. Turn off the spectrophotometer.</p>	<p>7a. On the "Example Laboratory Data Sheet" as "Time Absorbances (were) Read".</p> <p>8a. If you are establishing a standard curve, there are 4 standards.</p> <p>8b. If a standard curve is established, there is 1 check standard, 0.50 mg per liter.</p> <p>8c. Keep flasks tightly stoppered when not in use because the alcohol solvent can evaporate, thereby changing the concentration.</p> <p>8d. Use the same cell for each solution, rinsing it once with alcohol, then two times with the solution to be measured and wiping off the outside before inserting it into the instrument. (You can also use matched cells.)</p> <p>8e. Record results on the "Example Laboratory Data Sheet" in the column next to Column E.2.</p> <p>9a. Use the same cell and techniques as in 8d above.</p> <p>9b. Record the results on the "Example Laboratory Data Sheet" in the column next to Column F.8, the "Sample Identification Code."</p>	<p>IX.H.7 (p. 33)</p> <p>IX.H.8 (p. 33)</p> <p>IX.H.9 (p. 33)</p>

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
I. Making a Calibration Curve	<ol style="list-style-type: none"> 1. If a calibration curve has been established, omit this Operating Procedure and do Operating Procedure J, Checking the Calibration Curve. If a calibration curve has not been established, do the following steps. 2. Plot the absorbance values for the four standards obtained in H, Step 8 vs. the concentrations of boron in the calibration standards in E, Step 2. 3. Draw the best straight line through all the points to produce a calibration curve. 4. This curve can be used for Procedure K, Finding the Concentration of Boron, mg/liter for the Samples. 	<ol style="list-style-type: none"> 1a. There is an EMP on "Preparation of Calibration Graphs". 2a. All this information should be available from the "Example Laboratory Data Sheet". 2b. An example of the axes for the calibration graph is in the Training Guide. 	<p>IX. Sheet 1 (p. 33)</p> <p>IX. Sheet 2 (p. 34)</p>
J. Checking the Calibration Curve	<ol style="list-style-type: none"> 1. On the calibration curve, locate the absorbance value recorded for the check standard (0.50 mg/liter) in H, Step 8. 2. Draw a dotted line over to the calibration curve. 	<ol style="list-style-type: none"> 1a. This information should be available from the "Example Laboratory Data Sheet" in the column next to "Boron, mg/l in Standards". 1b. The curve was developed in Procedure I, Making a Calibration Curve. 	<p>IX. Sheet 1 (p. 33)</p> <p>IX. Sheet 2 (p. 34)</p>

EFFLUENT MONITORING PROCEDURE: Determination of Boron, Curcumin Method

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>J. Checking the Calibration Curve (continued)</p>	<p>3. From that point on the calibration curve, draw a perpendicular line down to the concentration line.</p> <p>4. Record the concentration of boron, mg/liter at this point for the check standard.</p> <p>5. Compare this observed concentration of boron, mg/liter to the expected concentration of 0.50 mg/liter.</p> <p>6. If the observed concentration of boron, mg/liter is within the acceptable range, the calibration curve can be used for K., Finding the Concentration of Boron, mg/liter for the Sample(s).</p>	<p>4a. Record the value on the "Example Laboratory Data Sheet" in the column next to the absorbance value for the 0.50 mg/liter standard.</p> <p>5a. The observed concentration should be within $\pm 2\%$ of the expected concentration of 0.50 mg/liter. 1) 2% of 0.50 is 0.01. 2) Thus the acceptable range of the observed value is 0.49 - 0.51 mg/liter.</p> <p>6a. If the observed concentration is not 0.49 - 0.51, you can re-run a 0.50 mg/liter check standard and repeat the comparison in step 5 above.</p> <p>6b. If the observed concentration still is not 0.49 - 0.51, discard the calibration curve and establish a new one by running standards and constructing a curve. Begin at E.1.</p>	<p>IX.J.4. (p. 33)</p>
<p>K. Finding the Concentration of Boron, mg/liter for the Sample(s)</p>	<p>1. On the calibration curve, locate the absorbance value recorded for the sample in H. Step 9.</p> <p>2. Draw a dotted line over to the calibration curve.</p> <p>3. From that point on the calibration curve, draw a perpendicular line down to the concentration line.</p>	<p>1a. This information should be available from the "Example Laboratory Data Sheet" in the column next to the "Sample Identification Code."</p> <p>2a. The curve was developed in Procedure I, Making a Calibration Curve.</p>	<p>IX. Sheet 1 (p. 33)</p> <p>IX. Sheet 2 (p. 34)</p>



OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
K. Finding the Concentration of Boron, mg/liter for the Sample(s) (continued)	4. Record the concentration of boron, mg/liter at this point for the sample. 5. Repeat steps 1 through 4 for each sample. 6. Sign the "Example Laboratory Data Sheet".	4a. Record the value on the "Example Laboratory Data Sheet" in the column next to the absorbance value for that sample. 6a. On the line marked "Analyst" in the top section.	IX.K.4. (p. 33) IX.K.6. (p. 33)

EFFLUENT MONITORING PROCEDURE: Determination of Boron, Curcumin Method

TRAINING GUIDE

<u>SECTION</u>	<u>TOPIC</u>
I*	Introduction
II	Educational Concepts-Mathematics
III	Educational Concepts-Science
IV	Educational Concepts-Communications
V	Field & Laboratory Equipment
VI	Field & Laboratory Reagents
VII	Field & Laboratory Analysis
VII*	Safety
IX*	Records & Reports

Training guide materials are presented here under the heading marked. These standardized headings are used throughout this series of procedures.

INTRODUCTION

TRAINING GUIDE NOTE

REFERENCES/RESOURCES

Boron may occur naturally in some waters or may be added to water in certain cleaning compounds or as an industrial waste effluent.

The determination of boron in waters, industrial wastes, and sewage effluents is particularly important to agriculture. Boron in small quantities is an essential element for plant growth. However, boron in excess of 2.0 mg/liter in irrigation water is harmful to most plants; some are affected by concentrations as low as 0.75 mg/liter. Drinking waters generally contain less than 0.1 mg/liter. Rarely do they contain more than 1.0 mg/liter which is still considered innocuous for human consumption.

The test described in this instruction can be found in Standard Methods on page 287, entitled 405A. Curcumin Method. This is the only procedure for Boron acceptable for NPDES purposes.

1. Water Quality Criteria, 1972. U.S. EPA Government Printing Office, No. 5501-00520, Washington, DC, p. 341.
2. Standard Methods for the Examination of Water and Wastewater, 14th ed., 1976, APHA, Washington, DC, p. 287.

EFFLUENT MONITORING PROCEDURE: Determination of Boron, Curcumin Method

RECORDS AND REPORTS

SECTION IX

SHEET 1
EXAMPLE LABORATORY DATA SHEET

Name of Plant _____

STEP		SAMPLE	SAMPLE	SAMPLE	SAMPLE
A.5	Identification Code				
A.5	Type (grab or composite)				
A.5	Date and Time Collected				
A.5	Sample Collector				
A.6	Date and Time Analysis Began				
F.15	Time Evaporation Ended				
H.7	Time Absorbances Read				
K.6	Analyst				

E.2 E.14	Concentration of Boron, mg/l in Standards	H.8	Absorbance	J.4	Concentration of Boron, mg/l in Check Standard
	0.00				
	0.20				
	0.50				
	0.80				
	1.00				

F.8	Sample Identification Code	H.9	Absorbance	K.4	Concentration of Boron, mg/l in Sample

EFFLUENT MONITORING PROCEDURE: Determination of Boron, Curcumin Method

RECORDS AND REPORTS

SECTION IX

SHEET 2

DETERMINATION OF BORON, CURCUMIN METHOD

CALIBRATION GRAPH

SIGNATURE OF PREPARER: _____

DATE GRAPH WAS PREPARED: _____

ABSORBANCE

0.60

0.50

0.40

0.30

0.20

0.10

0.00

0.00 0.10 0.20 0.30 0.40 0.50 0.60 0.70 0.80 0.90 1.00

CONCENTRATION OF BORON mg/l

61

62

A PROTOTYPE FOR DEVELOPMENT OF
ROUTINE OPERATIONAL PROCEDURES

for the

DETERMINATION OF TOTAL CALCIUM
(VOLUMETRIC METHOD)

as applied in

WASTEWATER TREATMENT FACILITIES
and in the
MONITORING OF EFFLUENT WASTEWATERS

Developed by the

National Training and Operational Technology Center
Municipal Operations and Training Division
Office of Water Program Operations
U.S. Environmental Protection Agency

EFFLUENT MONITORING PROCEDURE: Volumetric Determination of Total Calcium

This Operational Procedure was developed by:

NAME William T. Engel

ADDRESS Charles County Community College
P. O. Box 910
La Plata, Maryland 20646

POSITION Associate Professor of Chemistry

EDUCATION AND TECHNICAL BACKGROUND

BS - Saint Francis College, Loretto, Pennsylvania

MS - Xavier University, Cincinnati, Ohio

4 years Instructor (Chemistry)

2 years Assistant Director Pollution Abatement
Technology Department

EFFLUENT MONITORING PROCEDURE: Volumetric Determination of Total Calcium

1. Objective:

To determine the total calcium content of a wastewater effluent sample.

2. Description of Analysis:

The procedure involves a pretreatment of the sample with nitric acid which eliminates interferences from suspended material normally present in wastewater. The sample is subsequently analyzed by the EDTA titrimetric method. The results are reported as Total Calcium in mg/l.

Source of Procedure: Standard Methods for the Examination of Water and Wastewater, 13th Edition, APHA, AWWA, WPCF, 1971. pages 84, 85, and 416.

Methods for Chemical Analysis of Water and Wastes, 1974 Edition, Methods Development and Quality Assurance Research Laboratory, National Environmental Research Center, Cincinnati, Ohio 45268. pages 82-83.

EFFLUENT MONITORING PROCEDURE: Volumetric Determination of Total Calcium

Equipment and Supply Requirements

A. Capital Equipment

1. Balance - analytical 0.1 mg accuracy
2. Balance - triple beam 0.1 g accuracy
3. pH meter
4. Hot plates must heat to above 100°C

B. Reusable Supplies

1. Beakers - 50 ml, 250 ml, and 600 ml
2. Bottles - storage B/S 8 and 32 oz.
3. Buret - 50 ml
4. Cylinders - graduated, 25 ml, 50 ml and 500 ml
5. Flask - Erlenmeyer, 125 ml, and 500 ml
6. Pipets - Mohr, 2 ml, 5 ml, and 10 ml
7. Pipets - Volumetric, 10 ml and 25 ml
8. Ring stand
9. Flask - Volumetric, 100 ml, 500 ml, and 1000 ml
10. Clamp - Buret
11. Funnel - 60°
12. Plastic wash bottle
13. Measuring spoon - 0.2 g

C. Consumable Supplies

1. EDTA (Ethylenedinitriloacetic acid disodium salt)
($\text{Na}_2\text{H}_2\text{C}_{10}\text{H}_{12}\text{O}_8\text{N}_2 \cdot 2\text{H}_2\text{O}$)
2. Calcium carbonate (CaCO_3) anhydrous powder
3. Hydrochloric acid - concentrated (HCl)
4. Ammonium hydroxide - concentrated (NH_4OH)
5. Methyl red indicator
6. Nitric acid - concentrated (HNO_3)
7. Filter paper (#42)
8. Sodium hydroxide (NaOH)
9. Sodium chloride (NaCl)
10. Eriochrome blue black R indicator (Calcon)

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
DETERMINATION OF TOTAL CALCIUM (VOLUMETRIC METHOD), mg/l			
<p>A. Equipment Preparation</p> <p>1. Glassware Wash-up</p> <p>2. Balance Inspection</p> <p>3. pH Meter Inspection</p>	<p>1. Clean all glassware in a suitable detergent.</p> <p>1. Clean all balances - analytical balance and triple beam balance.</p> <p>1. Use EMP on pH to calibrate the pH meters.</p>	<p>1a. Distilled water drains without leaving any droplets.</p> <p>1a. Free of dust and dirt.</p> <p>1a. See EPA-430/1-74-015, Outline #5.</p>	
<p>B. Reagent Preparation</p> <p>1. Distilled Water</p> <p>2. Hydrochloric Acid (HCl 1:1)</p>	<p>1. Deionized distilled water should be prepared by passing distilled water through a mixed bed of cation and anion exchange resins.</p> <p>1. If metal impurities are present, follow the same distillation procedure as for the nitric acid.</p> <p>2. Pour 400 ml of distilled water in a 1 liter volumetric flask.</p> <p>3. Add 500 ml of concentrated hydrochloric acid to the flask and mix.</p>	<p>1a. All reagents and calibration standards should be prepared with this water.</p> <p>3a. Transfer the acid in fume hood.</p>	<p>68</p>

EFFLUENT MONITORING PROCEDURE: Volumetric Determination of Total Calcium

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>B. Reagent Preparation (Continued)</p>	<p>4. Mix thoroughly. 5. Dilute to 1 liter. 6. Transfer to a 1 liter storage bottle and label as 1:1 hydrochloric acid (HCl).</p>		
<p>3. EDTA Titrant (0.01M) (Ethylenedini- triolo tetra- acetic acid disodium salt) $\text{Na}_2\text{H}_2\text{C}_{10}\text{H}_{12}\text{O}_8\text{N}_2 \cdot 2\text{H}_2\text{O}$</p>	<p>1. Weigh 3.723 grams of EDTA in a weighing boat. 2. Dissolve the EDTA in approximately 800 ml of distilled water in a 1 liter volumetric flask. 3. Dilute to 1 liter. 4. Transfer to a storage bottle and label.</p>	<p>1a. One milliliter of 0.01 M EDTA titrant is equivalent to 400.8 μg calcium.</p>	
<p>4. Methyl Red Indicator</p>	<p>1. Weigh 20 mg of methyl red indicator in a weighing boat. 2. Dissolve in a mixture of 60 ml of ethyl alcohol and 40 ml of distilled water. 3. Transfer the indicator solution to a 100 ml storage bottle and label appropriately.</p>		

OPERATING PROCEDURES	STEP SEQUENCE	
<p>B. Reagent Preparation (Continued)</p> <p>5. Ammonium Hydroxide Solution (NH_4OH 3N)</p>	<p>1. Measure 203 ml of reagent grade ammonium hydroxide (NH_4OH) in a graduated cylinder.</p> <p>2. Transfer to 500 ml of distilled water in a 1 liter volumetric flask.</p>	
<p>6. Standard Calcium Solution</p>	<p>3. Mix the solution and dilute to 1 liter.</p> <p>4. Transfer to a storage bottle and label.</p> <p>1. Weigh 1.000 gram anhydrous calcium carbonate (CaCO_3) in a weighing boat.</p> <p>2. Transfer to a 500 ml Erlenmeyer flask.</p> <p>3. Add 1:1 HCl to the flask until all the calcium carbonate has dissolved.</p> <p>4. Add 200 ml distilled water to the flask.</p> <p>5. Boil the flask for a few minutes.</p> <p>6. Cool the solution and add 4 drops of methyl red indicator.</p>	5a.

INFORMATION/OPERATING GOALS/SPECIFICATIONS

T
GU

agent
oxide
ed

f

lask.

dilute

e

a

lask

ved.

water

few

5a. This will expel excess carbon dioxide (CO₂).

add

72

TRAINING
GUIDE NOTES

72

EFFLUENT MONITORING PROCEDURE: Volumetric Determination of Total Calcium

OPERATING PROCEDURES.	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS.	TRAINING GUIDE NOTES
<p>B. Reagent Preparation (Continued)</p>	<p>7. The color must now be adjusted to an intermediate orange color by adding 3N NH₄OH or 1:1 HCl as required.</p>		
<p>7. Sodium Hydroxide Solution (NaOH 1N)</p>	<p>8. Transfer the solution to a 1 liter volumetric flask and dilute to 1 liter with distilled water.</p> <p>9. Transfer to a storage bottle and label.</p> <p>1. Weigh out 40 grams of sodium hydroxide.</p> <p>2. Dissolve in 600 ml of distilled water in a 1 liter volumetric flask.</p> <p>3. Dilute to 1 liter.</p> <p>4. Transfer to a storage bottle and label.</p>	<p>9a. This solution is equivalent to 1.00 mg CaCO₃ per 1.00 ml.</p>	
<p>8. Eriochrome Blue Black R Indicator</p>	<p>1. Grind together 200 mg of Eriochrome Blue Black R dye with 100 grams of solid sodium chloride (NaCl).</p> <p>2. Store in tightly stoppered bottle.</p>	<p>1a. Use a mortar and pestle.</p> <p>1b. 0.2 grams will be used for each titration, it is therefore advantageous to have a 0.2 gram transfer spoon available.</p>	

73

74

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>C. Standardization of EDTA Titrant</p>	<p>1. Measure 25.0 ml of the standard calcium solution into a 125 ml Erlenmeyer flask.</p>		
	<p>2. Add 2.0 ml NaOH solution to the flask.</p> <p>3. Transfer 0.2 grams of the Eriochrome Blue Black R indicator to the flask.</p> <p>4. Add the EDTA titrant from the buret slowly and with swirling until the color changes from red through purple to bluish purple to a pure blue.</p> <p>5. Record the ml of EDTA on the data sheet.</p>	<p>2a. This solution should now be in the pH range of 12-13.</p> <p>2b. A pH meter may be used to check the solution.</p> <p>5a. If the concentration of EDTA was exactly 0.01M, then the titration should require 25 ml and the value of B in the formula:</p> $\text{mg/l Ca} = \frac{A \times B \times 400.8}{\text{ml sample}}$ <p>would be equal to (1).</p> <p>5b. B is the mg of CaCO₃ equivalent to 1.00 ml of EDTA titrant.</p> <p>(Continued)</p>	



EFFLUENT MONITORING PROCEDURE: Volumetric Determination of Total Calcium

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
C. Standardization of EDTA Titrant (Continued)		5c. Therefore: $B = \frac{\text{ml of EDTA titrant required in your standardization}}{25 \text{ ml}}$	
		If ml of titrant = 25 ml then $B = 25/25 = 1$ If ml of titrant = 24 then $B = \frac{24}{25} = 0.96$	
D. Analysis of Calcium in the sample. 1. Pretreatment of Sample	1. Acidify the sample at the time of sample collection by the addition of conc. Nitric acid to a pH of 2. 2. Transfer 50 ml of well mixed sample to a 150 ml beaker. 3. Add 5 ml of 1:1 hydrochloric acid. 4. Place the beaker on a hot plate and heat at 95°C for 15 minutes.	1a. The nitric acid should be free from Ca or re-distilled. 1b. Use a pH meter to measure. 2a. Use a graduated cylinder. 3a. Use a graduated pipet. 3b. The hydrochloric acid should be free from calcium or redistilled. 4a. The sample should not boil during this time.	

77

78

OPERATING PROCEDURES	STFP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>D. Analysis of Calcium in the sample (continued)</p> <p>2. Titration</p>	<p>5. Cool to room temperature.</p> <p>6. Filter the sample.</p> <p>7. Adjust the sample volume to 50 ml with distilled water.</p> <p>8. Transfer the sample to a 125 ml Erlenmeyer flask.</p> <p>1. Using the digested sample, add 2.0 ml of NaOH (sodium hydroxide) solution to the beaker.</p> <p>2. Transfer 0.2 grams of the Eriochrome Blue Black R indicator to the beaker.</p> <p>3. Add the EDTA titrant from the buret slowly, and with thorough mixing, until the color changes from red through purple to bluish purple to a pure blue.</p>	<p>6a. Use filter paper.</p> <p>6b. Filter into a 50 ml graduated cylinder</p> <p>6c. Wash down the walls of the beaker and add to the flask.</p> <p>7a. Dilute to the mark on the volumetric flask.</p> <p>7b. The sample is now ready for analysis.</p> <p>8a. For titration.</p> <p>1a. This solution should now be in the pH range of 12-13.</p> <p>1b. A pH meter may be used to check the solution.</p> <p>2a. Use a 0.2 gram spoon.</p>	

EFFLUENT MONITORING PROCEDURE: Volumetric Determination of Total Calcium

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>D. Analysis of Calcium in the sample (continued)</p>	<p>4. Record the ml of EDTA on the data sheet.</p>	<p>4a. Calculation of Total Calcium (mg/l)</p> $\text{mg/l Total Ca} = \frac{A \times B \times 400.8}{\text{ml of sample}}$ <p>A = ml of EDTA used in titration B = mg CaCO₃ equivalent to 1.00 ml EDTA (See C.5 for the value)</p> <p>Example: A = 5.2 ml B = 0.96 ml of sample = 50</p> $\text{mg/l Total Calcium (Ca)} = \frac{5.2 \times 0.96 \times 400.8}{50} = 40 \text{ mg/l}$	

A PROTOTYPE FOR DEVELOPMENT OF
ROUTINE OPERATIONAL PROCEDURES

for the

DETERMINATION OF COPPER (Cu^{++}), MAGNESIUM (Mg^{++})
MANGANESE (Mn^{++}), AND ZINC (Zn^{++})

as applied in

WASTEWATER TREATMENT FACILITIES
and in the
MONITORING OF EFFLUENT WASTEWATERS

Developed by the

National Training and Operational Technology Center
Municipal Operations and Training Division
Office of Water Program Operations
U.S. Environmental Protection Agency

Ch.MET.aa.EMP.1b.7.77

EFFLUENT MONITORING PROCEDURE: Determination of Cu^{++} , Mg^{++} , Mn^{++} , and Zn^{++}

This instructional sequence was developed by:

NAME Paul F. Hallbach

ADDRESS EPA, OWPO, NTOTC, Cincinnati, Ohio 45268

POSITION Chemist-Instructor

EDUCATION AND TECHNICAL BACKGROUND

B.S. Chemistry

14 years Industrial Chemist

16 years HEW-FWPCA-EPA-Chemist

EFFLUENT MONITORING PROCEDURE * Determination of Cu^{++} , Mg^{++} , Mn^{++} , and Zn^{++}

1. Analysis Objective:

To determine the copper, magnesium, manganese and zinc concentration of an effluent.

2. Brief Description of Analysis:

The sample is digested with concentrated nitric acid and evaporated to dryness. The residue is treated with hydrochloric acid, silicates and other insoluble material are removed by filtration and the sample is analyzed for the total metals of interest by atomic absorption spectrophotometry.

Source of Procedure: Methods for Chemical Analysis of Water and Wastes, 1974, Environmental Protection Agency, Methods Development and Quality Assurance Research Laboratory, Cincinnati, Ohio, p. 82

EFFLUENT MONITORING PROCEDURE: Determination of Cu^{++} , Mg^{++} , Mn^{++} , and Zn^{++}

Operating Procedures:

- A. Sample Digestion
- B. Reagent Preparation
- C. Instrument Calibration
- D. Instrumental Analysis
- E. Calculations

EFFLUENT MONITORING PROCEDURE: Determination of Cu^{++} , Mg^{++} , Mn^{++} , and Zn^{++}

General Description of Equipment and Supplies Used in the Process

A. Capital Equipment

1. Balance, analytical - sensitivity 0.1 milligram
2. Atomic absorption spectrophotometer
3. pH meter
4. Hot plate, 110 V

B. Reusable Supplies

1. Flasks, volumetric 100 ml, 1000 ml pipets
2. Pipets, volumetric, 50 ml, 3 ml, 1 ml
3. Reagent bottles, glass with glass stopper
4. Anion and cation exchange resin cartridges
5. Beakers, 100 ml
6. pH paper
7. Watch glass
8. Funnel, 80 mm diameter
9. Ring stand and 3 inch ring
10. Graduates 50 ml, 10 ml

C. Consumable Supplies

1. Reagents

Copper metal (analytical reagent grade)
Magnesium oxide, analytical reagent grade
Manganese metal
Zinc metal

EFFLUENT MONITORING PROCEDURE: Determination of Cu^{++} , Mg^{++} , Mn^{++} , and Zn^{++}

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
A. Sample Digestion	<ol style="list-style-type: none"> 1. Transfer 50 ml of sample into a clean 100 ml beaker. 2. Check the pH using pyridron paper. 3. Add 3.0 ml of concentrated nitric acid. 4. Place the beaker on a hot plate. 5. Evaporate to dryness cautiously. 6. Remove the beaker from the hot plate. Allow it to cool to room temperature and add another 3 ml aliquot of concentrated nitric acid. 7. Cover the beaker with a watch glass and place it on the hot plate. 8. Increase the temperature of the hot plate so that a gentle reflux action occurs. 9. Add 5 ml of 1:1 HCl and again warm the beaker to dissolve the residue. 	<ol style="list-style-type: none"> 1a. Use a 50 ml pipet. 2a. The pH should be 2.0. If the sample was not acidified upon collection, add 1:1 nitric acid dropwise until the pH is adjusted to 2.0. 3a. Use a 3 ml pipet. Use a rubber bulb on the pipet. 4a. Adjust the hot plate for medium heat. 5a. Make certain that the sample does not boil. 6a. Use a 3.0 ml pipet. 8a. Continue heating adding additional acid as necessary, until the digestion is complete (generally indicated by a light colored residue). 	

88

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
A. Sample Digestion (Cont'd.)	10. Wash down the beaker walls and watch glass with distilled water. 11. Filter the sample through Whatman #42 filter paper into a clean 100 ml volumetric flask. 12. Dilute the flask to the mark with distilled water.	10a. Use a plastic wash bottle.	
B. Reagent Preparation			
1. Deionized Distilled Water	1. Prepare by passing distilled water through a mixed bed of cation and anion exchange resins.	1a. Use deionized distilled water for the preparation of all reagents, calibration standards and as dilution water.	
2. Nitric Acid Concentrated (HNO ₃)	1. Commercially available reagent grade.		
3. Hydrochloric Acid (HCl) 1:1	1. Prepare a 1:1 solution of reagent grade hydrochloric acid by adding 25 ml of commercially available reagent grade hydrochloric acid to 25 ml of deionized water.	1a. Use a 50 ml graduate	
4. Copper Stock Standard Solution	1. Carefully weigh 1.00 grams of electrolyte copper (analytical reagent grade) on an analytical balance.	1a. Use a plastic weighing dish.	

EFFLUENT MONITORING PROCEDURE: Determination of Cu^{++} , Mg^{++} , Mn^{++} , and Zn^{++}

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/ SPECIFICATIONS	TRAINING GUIDE NOTES
B. Reagent Preparation (Cont'd.)	2. Transfer the copper into a clean 100 ml beaker. Add 15 ml distilled water and 5 ml HNO_3 and dissolve.		
	3. Transfer the solution to a 1000 ml volumetric flask and dilute to the mark with deionized distilled water.	3a. Use a plastic wash bottle to rinse the beaker into the volumetric flask. 3b. One ml equals 1 mg copper (1000 mg/l):	
5. Magnesium Stock Standard Solution	1. Weigh 0.829 grams of magnesium oxide on an analytical balance.	1a. Use a plastic weighing dish.	
	2. Transfer the reagent into a 100 ml beaker and add 15 ml distilled water and 10 ml of concentrated HNO_3		
	3. After dissolution transfer into a clean 1000 ml volumetric flask and dilute to the mark with deionized water.	3a. Use a plastic wash bottle to rinse the beaker into the volumetric flask. 3b. One ml equals 0.50 mg Mg (500 mg/l).	
6. Manganese Stock Standard Solution	1. Weigh 1.000 grams of manganese metal on an analytical balance.	1a. Use a plastic weighing dish.	
	2. Transfer into a 100 ml beaker. Add 15 ml distilled water.		
	3. Add 10 ml HNO_3 and dissolve.	3a. Warm if necessary.	
	4. Transfer the solution into a 1000 ml volumetric flask.		

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>B. Reagent Preparation (Cont'd.)</p> <p>7. Zinc Stock Standard Solution</p> <p>8. Fuel and Oxidant</p>	<p>5. Dilute to the mark with 1 percent (v/v) HCl.</p> <p>1. Weigh 1.00 gram of zinc metal on an analytical balance.</p> <p>2. Transfer into a 100 ml beaker. Add 15 ml distilled water and 10 ml of concentrated HNO₃ and dissolve.</p> <p>3. Transfer the solution into a 1000 ml volumetric flask and dilute to the mark with deionized distilled water.</p> <p>1. Commercial grade acetylene is generally acceptable.</p> <p>2. Air may be supplied from a compressed air line, a laboratory compressor, or from a cylinder of commercial air.</p>	<p>5a. One percent (v/v) HCl can be made by adding 10 ml HCl to 990 ml of deionized water.</p> <p>5b. One ml equals 1 mg Mn (1000 mg/l)</p> <p>1a. Use a plastic weighing dish.</p> <p>2a. Use a 10 ml graduate.</p> <p>3a. Use a plastic wash bottle to rinse the beaker during the transfer.</p> <p>3b. One ml equals 1 mg Zn (1000 mg/l)</p> <p>2a. Caution: Air supply must be free from oil or other contaminants.</p>	
<p>C. Instrument Calibration</p>	<p>1. Turn on air supply.</p> <p>2. Turn on acetylene supply.</p> <p>3. Turn on instrument and ignite flame.</p>	<p>3a. See instruction manual for your particular instrument.</p>	

EFFLUENT MONITORING PROCEDURE: Determination of Cu⁺⁺, Mg⁺⁺, Mn⁺⁺, and Zn⁺⁺

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>C. Instrument Calibration (Cont'd.)</p>	<p>4. Turn on power to hollow cathode lamp.</p> <p>5. Select wave length for appropriate metal.</p> <p>6. Prepare a series of standard solutions for each metal as follows:</p> <p><u>Copper</u> Transfer 10.0 ml of stock copper solution into a 100 ml volumetric flask and dilute to the mark with deionized distilled water and shake well.</p> <p>Transfer 0.2, 0.4, 0.6, 0.8, and 1.0 ml of the diluted standard solution into each of five 100 ml volumetric flasks respectively. Dilute to the mark with deionized distilled water and shake well. The concentration of these solutions will be 0.2, 0.4, 0.6, 0.8, and 1.0 mg/l respectively.</p> <p><u>Magnesium</u> Transfer 1.0 ml of stock magnesium standard into a 100 ml volumetric flask and dilute to the mark with distilled water.</p>	<p>4a Select lamp for proper metal analysis.</p> <p>5a. Copper (324.7 nm), magnesium (285.2 nm), manganese (279.5 nm) and zinc (213.9 nm)</p> <p>Use a 1 ml micro pipet graduated in 0.1 ml.</p>	<p>96</p> <p>97</p>

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>C. Instrument Calibration (Cont'd.)</p>	<p>Transfer 0.2, 0.4, 0.6, 0.8, and 1.0 ml of the diluted standard solution into each of five 100 ml volumetric flasks respectively. Dilute to the mark with deionized distilled water and shake well. The concentration of these solutions will be 0.02, 0.04, 0.06, 0.08 and 0.10 mg/l respectively.</p> <p><u>Manganese</u> Transfer 10.0 ml of stock manganese solution into a 100 ml volumetric flask. Dilute to the mark with deionized distilled water and shake well.</p> <p>Transfer 0.2, 0.4, 0.6, 0.8, and 1.0 ml of the diluted standard into each five 100 ml volumetric flasks respectively. Dilute to the mark with deionized distilled water and shake well. The concentration of these solutions will be 0.2, 0.4, 0.6, 0.8, and 1.0 mg/l respectively.</p> <p><u>Zinc</u> Transfer 10.0 ml of stock zinc solution into a 100 ml</p>		



EFFLUENT MONITORING PROCEDURE: Determination of Cu⁺⁺, Mg⁺⁺, Mn⁺⁺, and Zn⁺⁺

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>C. Instrument Calibration (Cont'd.)</p>	<p>volumetric flask. Dilute to the mark with deionized distilled water and shake well.</p> <p>Transfer 0.2, 0.4, 0.5, 0.8, and 1.0 ml of the diluted standard into each of five 100 ml volumetric flasks respectively. Dilute the mark with deionized distilled water and shake well. The concentration of these solutions will be 0.2, 0.4, 0.6, 0.8, and 1.0 mg/l respectively.</p> <p>7. Ignite flame and aspirate standard solutions into the flame.</p> <p>8. Prepare a calibration curve by plotting the concentration of the respective metals against the response for each concentration.</p>	<p>8a. Record the response on a recorder or use the readout provided on the instrument.</p>	
<p>D. Instrumental Analysis</p>	<p>1. Aspirate the unknown solution into the instrument immediately following the aspiration of the standards.</p> <p>2. Record the response.</p>	<p>1a. Flame characteristics and instrumental settings should be the same for standards and unknowns.</p>	

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
E. Calculations	1. Determine the concentration of the metal in the sample by substituting the observed instrumental response on the appropriate calibration curve.		



A PROTOTYPE FOR DEVELOPMENT OF
ROUTINE OPERATIONAL PROCEDURES

for the

DETERMINATION OF LEAD BY ATOMIC ABSORPTION
USING THE EXTRACTION PROCEDURE

as applied in

WASTEWATER TREATMENT FACILITIES
and in the
MONITORING OF EFFLUENT WASTEWATERS

Developed by the

National Training and Operational Technology Center
Municipal Operations and Training Division
Office of Water Program Operations
U.S. Environmental Protection Agency

104

7 EFFLUENT MONITORING PROCEDURE: Determination of Lead by Atomic Absorption
Using the Extraction Procedure

This operational procedure was developed by:

NAME John D. Pfaff

- ADDRESS EPA, ONPO, NTOTC, Cincinnati, Ohio 45268

POSITION Chemist-Instructor

EDUCATION AND TECHNICAL BACKGROUND

B.S. Chemistry

3 years - Research Chemist

15 years - Training Instructor

**EFFLUENT MONITORING PROCEDURE: Determination of Lead by Atomic Absorption
Using the Extraction Procedure**

1. Analysis Objectives:

The learner will determine the lead content of a sample by extracting the sample and determining the results on an Atomic Absorption Spectrophotometer.

2. Brief Description of Analysis:

The sample is concentrated through chelation and extraction with organic solvents. The method produces a solution which can be run under standard conditions for atomic absorption. The solution is then aspirated into any atomic absorption instrument of sufficient sensitivity. This procedure contains directions for the use of an Instrumentation Laboratory Inc.* Model 153 Atomic Absorption Spectrophotometer.

3. Applicability of this Procedure:

a. Range of Concentration:

The extraction procedure is recommended for levels of lead below 100 micrograms per liter.

b. Pretreatment of Samples:

This EMP includes the digestion procedure specified in the Federal Register Guidelines to determine total lead concentrations.

c. Treatment of Interferences in Samples:

Since this EMP procedure involves chelation and extraction of lead from a water sample into organic solvents, the lead is separated from dissolved materials that might interfere in a direct aspiration. The EMP includes procedures or information about glassware preparation, reagent purity and the exceptional sensitivity of this analysis to turbulence and absorption bands in the flame. The choice of using a 217.0 or 283.3 nm wavelength is also discussed. A section about interferences to atomic absorption spectrophotometry (chemical, dissolved solids, ionization and spectral) can be found in the Source of Procedure.*

* Source of Procedure: Methods for Chemical Analysis of Water and Wastes, 1974, Environmental Protection Agency, Methods Development and Quality Assurance Research Laboratory, Cincinnati, Ohio 45268, Page 89 and 112.

*Mention of particular brand name does not constitute endorsement by the U.S. Environmental Protection Agency.

**EFFLUENT MONITORING PROCEDURE: Determination of Lead by Atomic Absorption
Using the Extraction Procedure**

- A. Preparing to Test the Solution
- B. Reagent Preparation
- C. Sample Preservation and Handling
- D. Instrument Set-Up
- E. Solubilization of Sample
- F. Preparation of Standard Dilutions
- G. Extraction Procedure
- H. Instrument Operation
- I. Calibration
- J. Instrument Shut-down
- K. Maintenance

SEE FIGURE 1 ON PAGE 31 IN THE TRAINING GUIDE FOR A "FLOW DIAGRAM FOR THE
EXTRACTION METHOD".

EFFLUENT MONITORING PROCEDURE: Détermination of Lead by Atomic Absorption
Using the Extraction Procedure

General Description of Equipment Used in the Process

A. Capital Equipment

1. Instrumentation Laboratories* Model 153 Atomic Absorption or equivalent
2. Hollow cathode lamp - Lead
3. Balance, analytical with a 0.1 milligram sensitivity
4. Pressure regulator valves:
 - a. Two stage regulator designed to deliver acetylene with an Inlet CGA 510 connector
 - b. Two stage regulator designed to deliver air from cylinder with an Inlet CGA 1340 connector
5. Balance, with a 0.1 or 0.01 gram sensitivity
6. Magnetic stirrer - hot plate and magnet retriever
7. Still, borosilicate glass distillation apparatus
8. Steam bath for 100 ml beakers, 6 plus X samples
9. pH meter (optional) to adjust pH for test

B. Reusable Supplies

1. Six plus one/sample. Beaker 100 ml capacity
2. One for each sample plus one more. Beaker 250 ml capacity
3. Three bottles, dropper, brown glass, 100 ml capacity
4. One bottle reagent, brown glass, 1000 ml capacity
5. Two bottles reagent, clear glass, 500 ml capacity
6. One cylinder, graduated, 250 ml capacity
7. Two cylinders, graduated, 500 ml capacity
8. Two flasks, Erlenmeyer, graduated, 500 ml capacity.
9. Two flasks, volumetric, 1000 ml capacity
10. Four flasks, volumetric, 100 ml capacity
11. Six plus one/sample. Flask, volumetric, wide base, 10 ml capacity
12. One for each sample, very small funnel to filter 3-5 ml
13. Six plus one/sample. Funnel separatory, 250 ml capacity
14. One pipet volumetric, 1 ml capacity
15. One pipet volumetric, 2 ml capacity
16. One pipet volumetric, 3 ml capacity
17. Three pipets volumetric, 5 ml capacity
18. Five pipets volumetric, 10 ml
19. One pipet volumetric, 20 ml
20. Three pipets graduated 1/10, 10 ml
21. One pipet graduated 1/10, 1 ml
22. Two pipets, graduated 1/10, 25 ml
23. One pipet, volumetric, 15 ml
24. One for each sample. Watch glasses, 3.5 inch diameter
25. Instrument manufacturer's manual on the atomic absorption instrument
26. Safety glasses
27. One separatory funnel rack
28. One wash bottle, plastic

* Mention of a particular brand name does not constitute endorsement by the U.S. Environmental Protection Agency.

EFFLUENT MONITORING PROCEDURE: Determination of Lead by Atomic Absorption
Using the Extraction Procedure

C. Consumable Supplies

1. Deionizing column - mixed bed type, such as Barnstead-Ultrapur-red cap (D0809) or Fisher #9-034-3
2. Gases:
 - Fuel - acetylene (C_2H_2) - for use with the atomic absorption instrument, "purified" grade, cylinder size (DOT-8AL) 380 cuf, CGA size 510
 - Oxidant-air - for use with the atomic absorption instrument, "dry" grade, cylinder size (DOT-3AA2015) 2200 cuf, CGA size 1340
3. Labels for reagent bottles - at least 7
4. Laboratory notebooks
5. Pencil, wax-marking
6. Reagents
 - a. Ammonium hydroxide (NH_4OH) - a test reagent, purchase reagent grade - 1 pt (473 ml) minimum
 - b. Bromphenol blue* - ($C_{19}H_{10}Br_4O_5S$) - a test reagent, purchase ACS grade - 5 gram minimum, may not be needed if pH meter is used to adjust pH
 - c. Carbon disulfide - (CS_2) - used to prepare a reagent, purchase reagent grade - 1 pint (473 ml) minimum
 - d. Chloroform ($CHCl_3$) - used to prepare a reagent, purchase reagent grade, 1 gal (3800 ml) minimum
 - e. Ethyl alcohol* - (CH_3CH_2OH) - used to prepare a reagent, purchase denatured - 1 pt (473 ml) minimum (May not be needed if pH meter is used to adjust pH)
 - f. Hydrochloric acid (HCl) - used as a reagent, purchase ACS grade, 1 pt (473 ml) minimum
 - g. Lead nitrate, anhydrous ($Pb(NO_3)_2$) - used to prepare standards, purchase ACS grade 1/4 pound or 100 grams minimum
 - h. Nitric acid (HNO_3) - used as a reagent, purchase ACS grade, 1 pt (473 ml) minimum
 - i. Pyrrolidine (C_4H_9N) - a test reagent, purchase reagent grade, 1/4 lb or 100 g minimum can be purchased from Aldrich Chemical Co.
940 W. St. Paul Ave.
Milwaukee, WI 53233
7. Filter paper, Whatman #42 to fit very small funnel

* If Bromphenol blue solution is purchased, these reagents need not be purchased.

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>A. Preparing to Test the Solution.</p>	<ol style="list-style-type: none"> 1. Clean all glassware to remove all metals. 2. Assemble equipment to be used. 3. Prepare all reagents for the test. 4. Optimize the atomic absorption instrument. 5. Solubilize the sample(s) and extract the sample(s) and standard dilutions. 6. Aspirate all into the AA. 7. Calculate results and record data. 	<ol style="list-style-type: none"> 1a. All glassware including sample bottles should be washed. 2a. See the equipment lists contained in this EMP. 3a. See procedures, Section B. 4a. See procedures, Sections D and H and manufacturer's manual. 5a. See procedures, Sections E, F, and G. 6a. See procedures, Section I. 7a. Record sample information in a lab notebook along with date and time of analysis. 7b. After analysis, record the analytical information beside sample information. 	<p>V.A.1. (p. 29)</p> <p>I.A.2. (p. 27)</p>
<p>B. Reagent Preparation</p> <ol style="list-style-type: none"> 1. Deionized distilled water 2. Nitric acid concentrated 	<ol style="list-style-type: none"> 1. Prepare by passing distilled water through a mixed bed of cation and anion exchange resins. 1. No preparation necessary. 	<ol style="list-style-type: none"> 1a. Use deionized distilled water for the preparation of all reagents, standards, blanks, and as dilution water. 1b. Use bed such as "Barnstead-Ultrapure" red cap (D0809) or Fisher #9-034-3. 1a. Use only reagent grade of analyzed purity. 	

EFFLUENT MONITORING PROCEDURE:

Determination of Lead by Atomic Absorption Using the Extraction Procedure

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>B. Reagent Preparation (continued)</p>			
<p>3. Nitric acid (1:1 dilution) HNO_3</p>	<p>1. Prepare by adding an equal volume of acid to an equal volume of water (i.e., 250 ml acid to 250 ml water).</p>	<p>1a. Caution: acid should always be added to water so as to avoid spattering. 1b. A graduated Erlenmeyer flask can be used for this preparation.</p>	
<p>4. Hydrochloric acid (0.60 N)</p>	<p>1. Add about 50 ml water to a 100 ml volumetric flask.</p> <p>2. Pipet 5 ml concentrated HCl to the flask.</p> <p>3. Cool and dilute to 100 ml with water.</p>	<p>1a. Use a reagent grade analyzed acid to avoid contamination. 1b. Caution: protective equipment should be used.</p> <p>2a. Use a 5 ml volumetric pipet.</p> <p>3a. Store in brown dropper-bottle (about 100 ml vol.) before use.</p>	
<p>5. Stock lead solution</p>	<p>1. Dissolve 1.599 g of analytical reagent grade lead nitrate ($Pb(NO_3)_2$) in water.</p> <p>2. Add 10 ml concentrated nitric acid (HNO_3)</p> <p>3. Dilute to 1 liter with deionized distilled water.</p>	<p>1a. Weighing should be done on an analytical balance. 1b. All reagents should be labeled with reagent name, concentration, date of preparation. 1c. Concentration of stock is 1 mg Pb/ml. 1d. Use a 1000 ml volumetric flask.</p> <p>2a. Use a graduate.</p>	
<p>6. Working lead standard solution.</p>	<p>1. Add 500 ml water to a volumetric flask.</p> <p>2. Add 10 ml concentrated Nitric Acid (HNO_3).</p>	<p>1a. Use a 1000 ml volumetric flask. 1b. Should be prepared fresh each day an analysis is to be done.</p> <p>2a. Use a graduate.</p>	

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>B. Reagent Preparation (continued)</p> <p>7. Pyrrolidine dithiocarbamic acid (PDCA)</p>	<p>3. Add 10.0 ml stock lead solution (Reagent 5) to the flask.</p> <p>4. Dilute to the mark with water.</p> <p>1. Add 500 ml chloroform to a liter flask.</p>	<p>3a. Use a volumetric pipet.</p> <p>4a. Concentration is 0.01 mg Pb/ml.</p> <p>1a. This reagent should be prepared in a well ventilated area (or hood)</p> <p>1b. Measure 500 ml with graduated cylinder.</p>	
	<p>2. Add 18 ml of analytical grade pyrrolidine.</p> <p>3. Add 15 ml of carbon disulfide (CS₂) in small portions with swirling.</p> <p>4. Dilute to 1 liter with chloroform.</p>	<p>2a. Pipet with a graduated pipet.</p> <p>2b. Generates heat -- cool before proceeding.</p> <p>2c. For supplier, see chemical list.</p> <p>2d. CAUTION -- reagent is flammable, toxic and corrosive.</p> <p>3a. Carbon disulfide is very odorous. Prepare in hood or well ventilated area.</p> <p>3b. Use a measuring pipet.</p> <p>3c. CAUTION - Heat generated - cool before proceeding.</p> <p>4a. This solution can be stored for several months if stored in a brown bottle in a refrigerator.</p>	
<p>8. Ammonium hydroxide (NH₄OH) concentrated</p>	<p>1. Pour the concentrated NH₄OH into a glass dropper bottle.</p>	<p>1a. Use a hood to prevent inhalation of fumes. Avoid contact with skin. Wear protective equipment.</p> <p>1b. Only some drops of this are needed for the pH adjustment of acidified samples.</p> <p>1c. A brown glass dropper bottle conserves the stability of this reagent.</p>	

EFFLUENT MONITORING PROCEDURE:

Determination of Lead by Atomic Absorption Using the Extraction Procedure

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>B. Reagent Preparation (continued)</p> <p>9. Ammonium hydroxide (NH₄OH) 2N</p> <p>10. Bromphenol blue indicator</p> <p>11. Hydrochloric Acid (1:1 dilution) HCl</p>	<p>1. Dilute 3 ml of concentrated NH₄OH to 100 ml with water.</p> <p>1. Dissolve 0.1 g of the solid in 100 ml of 50% ethyl alcohol.</p> <p>1. Prepare by adding an equal volume of acid to an equal volume of water (i.e., 250 ml acid to 250 ml water).</p>	<p>1a. This reagent should be prepared in the hood to prevent inhalation of fumes. Avoid contact with skin. Wear protective equipment.</p> <p>1b. Only several drops are needed per sample.</p> <p>1c. Use a brown dropper bottle to store.</p> <p>1a. A platform balance can be used for weighing.</p> <p>1b. In order to prepare this solution 95% ethyl alcohol should be diluted in half (i.e., 50 ml alcohol to 50 ml water).</p> <p>1c. This solution is stable indefinitely so long as it is kept tightly stoppered to prevent evaporation.</p> <p>1a. CAUTION: Acid should always be added to water so as to avoid spattering.</p> <p>1b. CAUTION: Heat is generated.</p> <p>1c. Use safety glasses.</p>	
<p>C. Sample Preservation and Handling</p>	<p>1. Collect a representative sample.</p> <p>2. Preserve the sample by adjusting to pH 2. Use 3 ml of 1:1 redistilled nitric acid per liter.</p>	<p>1a. Collect about 1 liter or 1 quart to assure ample amount for analysis.</p> <p>2a. Usually this amount is sufficient. However, some highly buffered waters might require more.</p> <p>2b. Check the pH of the sample with litmus paper to assure that the pH is 2 or below.</p>	

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>D. Instrument Set-Up</p> <p>1. Pre warm-up</p> <p>2. Lamp installation</p>	<p>1. Prepare the instrument for initial operation.</p> <p>1. Install lead hollow cathode lamp.</p>	<p>Ta. Reference is made to the manufacturer's manual of instruction.</p> <p>Tb. Check power requirements and availability.</p> <p>Tc. Provide adequate ventilation, including vent over instrument burner.</p> <p>Td. Provide adequate space for instrument and work area.</p> <p>Te. Provide drain facility.</p> <p>1a. The trimpot for the position into which the lamp is installed should be adjusted until the lamp draws its optimum lamp current. Set the lead lamp to 5 mA.</p> <p>1b. Do not exceed the maximum current rating for the lamp as this can seriously affect its life and stability.</p> <p>1c. Refer to the operation manual for proper installation procedure.</p>	<p>I.D.2. (p. 28)</p>
<p>3. Burner optimization</p>	<p>1. Attach the necessary tanks of gases to the instrument.</p> <p>2. Follow manufacturer's instructions and optimize the burner.</p>	<p>1a. For the lead procedure, acetylene and air are required. Use purified grades of gases.</p> <p>1b. Attach pressure regulator to each cylinder. Care should be taken to match C.G.A. numbers.</p> <p>1c. Connect cylinders through the regulators to the rear of the instrument at the marked positions.</p> <p>1d. All cylinders should be securely chained to prevent them from tipping over.</p> <p>2a. The analysis of lead is exceptionally sensitive to turbulence and absorption bands in the flame. Therefore, some care should be taken to position the light beam in the most stable, center portion of the flame. To do this, first adjust the burner to maximize the absorption reading with a lead standard. Then aspirate a water blank and make minute adjustments in the burner alignment to minimize the signal.</p>	<p>V.D.3.1a. (p. 29)</p>

EFFLUENT MONITORING PROCEDURE: Determination of Lead by Atomic Absorption Using the Extraction Procedure

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>E. Solubilization of Sample</p>	<ol style="list-style-type: none"> 1. Acidify the sample with 1:1 nitric acid to a pH of 2 at the time of sample collection. 2. Transfer 200 ml of well mixed sample to a 250 ml beaker. 3. Add 3 ml of concentrated nitric acid to the sample. 4. Place the beaker on a hot plate and cautiously evaporate to dryness. 5. Cool the beaker and add another 3 ml portion of the concentrated nitric acid. 6. Cover the beaker with a watch glass and return to the hot plate. 7. Increase the temperature of the hot plate so that a gentle reflux action occurs. 8. Continue heating until a light colored residue forms. 9. Add sufficient 1:1 hydrochloric acid (HCl) solution and again warm the beaker to dissolve the residue. 	<ol style="list-style-type: none"> 1a. For most samples, 3 ml of 1:1 HNO₃ per liter of sample lowers the pH sufficiently. 1b. Highly buffered samples require more acid. If in doubt, check the pH. 2a. 200 ml is the usual sample volume for concentrations containing less than 100 µg/liter. 2b. Choose a volume appropriate to the expected level of metals. 3a. Use a 10 ml graduated pipet. 4a. The sample should not boil during this time. 5a. Use a 10 ml graduated pipet. 7a. This will be indicated by droplets forming on the underside of the watch glass. 8a. This will indicate that the digestion is complete. 8b. Additional acid may be required to complete this digestion. 9a. Use a 1 ml graduated pipet. 9b. Use the watch glass again to cover the beaker. 	<p>VII.E.1. (p. 30)</p>

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES																																								
<p>E. Solubilization of Sample (continued)</p>	<p>10. Wash down the beaker walls and watch glass with distilled water.</p> <p>11. Filter the sample into a 250 ml graduated beaker.</p> <p>12. Adjust the volume to 200 ml (or the starting volume size).</p> <p>13. The sample is now ready for the Extraction Procedure (G).</p>	<p>11a. Use a filter paper such as Whatman #42.</p> <p>11b. To remove silicates and other insoluble material that could clog the atomizer of the atomic absorption instrument.</p> <p>12a. The sample is now ready for extraction.</p> <p>12b. The graduations on the beaker can be used to measure the volume.</p> <p>12c. Results shall be reported as "total".</p>																																									
<p>F. Preparation of Standard Dilutions</p>	<p>1. Prepare a blank and series of standard dilutions in 250 ml separatory funnels.</p>	<p>1a. Using the working lead standard (Reagent 6) prepare the blank and standards using volumetric pipets (1, 5, 10, 20 ml) for measuring the working lead standard and a graduated cylinder for the water. Prepare as follows, using Columns A and B.</p> <table border="1" data-bbox="1099 994 1854 1350"> <thead> <tr> <th>A</th> <th>B</th> <th>C</th> <th>D</th> <th>E</th> </tr> <tr> <th>ml's of Std. in 200 ml.</th> <th>ml's of water</th> <th>mg Pb/l conc. in 200 ml</th> <th>mg Pb/l conc. in final 10 ml</th> <th>Instrument Reading</th> </tr> </thead> <tbody> <tr> <td>0.0</td> <td>200</td> <td>0.0</td> <td>0.0</td> <td>_____</td> </tr> <tr> <td>1.0</td> <td>199</td> <td>0.050</td> <td>1.0</td> <td>_____</td> </tr> <tr> <td>5.0</td> <td>195</td> <td>0.250</td> <td>5.0</td> <td>_____</td> </tr> <tr> <td>10.0</td> <td>190</td> <td>0.500</td> <td>10.0</td> <td>_____</td> </tr> <tr> <td>15.0</td> <td>185</td> <td>0.750</td> <td>15.0</td> <td>_____</td> </tr> <tr> <td>20.0</td> <td>180</td> <td>1.000</td> <td>20.0</td> <td>_____</td> </tr> </tbody> </table>	A	B	C	D	E	ml's of Std. in 200 ml.	ml's of water	mg Pb/l conc. in 200 ml	mg Pb/l conc. in final 10 ml	Instrument Reading	0.0	200	0.0	0.0	_____	1.0	199	0.050	1.0	_____	5.0	195	0.250	5.0	_____	10.0	190	0.500	10.0	_____	15.0	185	0.750	15.0	_____	20.0	180	1.000	20.0	_____	<p>VII.F.1 (p. 30)</p> <p>123</p>
A	B	C	D	E																																							
ml's of Std. in 200 ml.	ml's of water	mg Pb/l conc. in 200 ml	mg Pb/l conc. in final 10 ml	Instrument Reading																																							
0.0	200	0.0	0.0	_____																																							
1.0	199	0.050	1.0	_____																																							
5.0	195	0.250	5.0	_____																																							
10.0	190	0.500	10.0	_____																																							
15.0	185	0.750	15.0	_____																																							
20.0	180	1.000	20.0	_____																																							

122

EFFLUENT MONITORING PROCEDURE:

Determination of Lead by Atomic Absorption Using the Extraction Procedure

OPERATING PROCEDURE	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
F. Preparation of Standard Dilutions (continued)	2. Carry these standards through the extraction procedure, beginning at G.1, pH adjustment.	1b. Make sure stopcock on separatory funnel is closed before any addition. 1c. Mark the mls of standard used (A) on the funnels. 1d. These standard dilutions should be prepared fresh for each analysis.	
G. Extraction Procedure	1. Place a 250 ml separatory funnel into the ring on the stand. 2. Add 200 ml of sample to the separatory funnel.	1a. Check the stop-cock on the separatory funnel to assure that it is closed. NOTE: If a pH meter with two electrodes is to be used to adjust the pH in the next steps, do those steps before adding sample to the separatory funnel. If using a single combination electrode, this can be done in the separatory funnel. 2a. If you have a solubilized sample from Procedure E, pour the 200 ml from the beaker into the funnel, and rinse the beaker into the funnel. 2b. In other cases, use a 500 ml graduated cylinder to measure the necessary volume. Add rinse. 2c. Record sample volume used for analysis. 2d. If use of 200 ml of sample gives results too high to be on scale, a dilution can be made. However, the dilution should be made to provide a 200 ml final volume. 2e. If use of a 200 ml sample gives too low results, more than one sample aliquot can be treated and the extracts combined.	

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>G. Extraction Procedure (continued)</p> <p>1. pH adjustment</p> <p>2. Chelation and Extraction</p>	<p>1. Add 2 drops bromphenol blue indicator to the sample(s) and to each standard.</p> <p>2. Mix well.</p> <p>3. Add ammonium hydroxide drop-wise until a very pale blue color persists.</p> <p>4. Add 0.6 N hydrochloric acid (Reagent 4) drop-wise until the blue color just disappears.</p> <p>5. Add 2.0 ml 0.6 N hydrochloric acid (Reagent 4). Stopper and shake.</p> <p>1. Add 5.0 ml pyrrolidine dithiocarbamic acid (PDCA) (Reagent 7)</p> <p>2. Shake vigorously for 2 min.</p>	<p>1a. Should a pH meter be used to adjust the pH; omit steps 1 through 4 and begin at step 5.</p> <p>1b. Apply each step to each sample and standard.</p> <p>3a. Use concentrated NH_4OH for acidified samples and standards made with 10 or more ml of acidified working lead standard. Use 2 N NH_4OH for more dilute standards.</p> <p>3b. The reagents should be in glass dropper bottles for this addition. Use a hood.</p> <p>4a. Use a glass dropper bottle for this addition. The bottle should be labeled HCl-.6N.</p> <p>4b. Pale yellow color may appear.</p> <p>5a. Use a 2.0 ml volumetric pipet.</p> <p>5b. If a pH meter is used, add acid until the pH is 2.3. Then, the next procedure will bring the pH to the optimum of 2.8.</p> <p>1a. Use a 5.0 ml volumetric pipet for this step.</p> <p>1b. This reagent should be allowed to come to room temperature before pipetting, since it will be stored in a refrigerator.</p> <p>1c. The bottle should be restoppered immediately after use and returned to the refrigerator to prolong usefulness.</p> <p>2a. CAUTION: Use proper technique with the separatory funnel. The reagent contains volatile solvents and pressure is formed.</p>	

EFFLUENT MONITORING PROCEDURE:

Determination of Lead by Atomic Absorption Using the Extraction Procedure

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
G. Extraction Procedure (continued)	3. Allow the reagent to settle to the bottom of the separatory funnel.	3a. Enough time should be allowed for complete separation of the two phases. 3b. It may take up to 3 minutes.	
	4. Open stopcock and slowly drain off the reagent phase into a 100 ml beaker.	4a. Mark the beaker with the number of ml used to prepare the <u>standard</u> or with the <u>sample</u> identification code.	
	5. Add a second 5.0 ml of PDCA reagent to the separatory funnel.	5a. The same volumetric pipet can be used for all additions of PDCA to all samples and standards provided caution is used to prevent contamination.	
	6. Shake vigorously for two minutes.		
	7. Allow reagent to settle and separate.	7a. This should take about 2-4 minutes.	
	8. Open stopcock and slowly drain off reagent phase.	8a. The reagent should be drained into the same beaker used in step 4. This will combine both the extraction volumes. 8b. Pale pink color may show in extracts.	
3. Recovery of Complex	1. Evaporate to dryness on a steam bath in hood.	1a. Residue is light color with possible pale green or blue tinges. 1b. Do not "bake" the residue. 1c. Should take about 10-15 minutes.	
	2. Remove and cool 2 minutes.		
4. Digestion of Complex	1. Add 2 ml concentrated nitric acid (HNO ₃)	1a. The concentrated HNO ₃ must be a good grade as any lead in the acid will be concentrated along with the sample. 1b. Best carried out in a hood. This is a violent reaction; bubbling, dark brown fumes given off.	

(Continued)

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>G. Extraction Procedure (continued)</p>	<p>2. Place the beaker on a low temperature hot plate and evaporate just up to dryness.</p> <p>3. Remove from hot plate and cool for 2 minutes.</p>	<p>1c. Use a measuring pipet and add acid down walls, dropwise at first until most residue dissolves.</p> <p>1d. Hold beaker at a 45 degree angle and slowly add the acid. Rotate the beaker while adding to effect thorough contact of the acid with the residue.</p> <p>2a. Care should be taken to remove the beaker when only a very small amount of deep brown liquid remains in the beaker.</p> <p>2b. The evaporation should take about 8 minutes.</p>	
<p>5. Dissolving the Residue</p>	<p>1. Add 2 ml of 1:1 nitric acid (HNO_3) - Reagent 3.</p> <p>2. Return sample to the low temperature hot plate and heat for 1 minute.</p> <p>3. Cool and quantitatively transfer the solution to a 10 ml volumetric flask -- bring to final volume mark.</p> <p>4. The sample is now ready for aspiration into the atomic absorption instrument.</p>	<p>1a. Use a measuring pipet.</p> <p>1b. Down inside walls at first.</p> <p>2a. Both standards and samples should be at this point together and carried through this step as it could affect the final concentration of acid.</p> <p>3a. A stirring rod and a plastic wash bottle containing deionized distilled water should be used to wash the beaker for transfer.</p> <p>3b. A wide base 10 ml volumetric flask is suggested or place the volumetric flask in a beaker to prevent tipping it over. A 10 ml stoppered graduated cylinder can be used instead of a volumetric flask.</p> <p>3c. Mark the flask or cylinder with the number of ml used to prepare the standard. This is also the mg Pb/l concentration in this final 10 ml. For samples, mark the identification code.</p>	

EFFLUENT MONITORING PROCEDURE:

Determination of Lead by Atomic Absorption Using the Extraction Procedure

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>H. Instrument Operation</p>	<ol style="list-style-type: none"> 1. Rotate the lamp turret until the lead lamp is in the light path. 2. Depress the power switch to the on position. 3. Depress the hollow cathode power switch to the on position. 4. Set mode switch at "A". 5. Set absorption/emission switch for channel "A" at absorbance. 6. Adjust channel "A" hollow cathode lamp current. 7. Adjust burner position. 8. Set monochromator to proper wavelength. 	<ol style="list-style-type: none"> 1a. See instrument manual for location and explanation of each control or switch. 1b. Figures 2 through 5 in the Training Guide show locations of controls. 2a. Properly connected to the power line, this switch controls all power except that of the hollow cathode lamps. 2b. Figure 3 - location of controls. 3a. In the on position, connects the hollow cathode lamps to their power supply. 3b. Figure 3. 4a. Selects the single channel direct readout mode of operation. 4b. See Figure 5. 5a. Selects the negative logarithmic absorbance for channel "A". 5b. See Figure 5. 6a. For the lead lamp, this should be 5.0 mA. 6b. See Figure 5. 7a. For the lead analysis a Rolling Burner should be used. This is the standard burner head supplied with the IL-153 instrument. 7b. The burner height should be adjusted to 8 mm for lead analysis. 7c. See instrument manual for procedures on burner adjustment. 8a. For the lead analysis the wavelength should be set for 283.3 nm. <p style="text-align: right;">(Continued)</p>	<p>VII.H.1: (pp. 32-35)</p>

132

133



OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>H. Instrument Operation (continued)</p>	<p>9. Set slit width.</p> <p>10. Peak the wavelength.</p> <p>11. Set I and I₀ indicators in the green balance zones.</p> <p>12. Ignite the flame.</p>	<p>8b. The 283.3 nm lead line is about half as sensitive as that of the 217.0 nm line. However, the 283.3 line has less background absorbance and is correspondingly less noisy. Consequently, the operator should experiment to decide which line to use.</p> <p>8c. See Figure 4 for location of controls.</p> <p>9a. The slit setting for the lead procedure should be 80 or position 4 on the instrument.</p> <p>9b. See Figure 4.</p> <p>10a. Turn the wavelength selector and watch the channel "A" Incident Intensity Indicator. Adjust the wavelength selector knob past the indicated wavelength in both directions and note where the maximum upward deflection on the indicator occurs. Place the wavelength adjustment knob at the position where the maximum upward deflection is obtained.</p> <p>11a. Adjust the channel "A" incident intensity control to a position so that the indicator needle on the channel "A" indicator rests approximately in the center of the green balance zone at the top of the meter. Adjust the transmitted intensity control so that the indicator needle on the indicator is parallel to the needle on the channel "A" incident intensity indicator.</p> <p>11b. See Figure 5 for location.</p> <p>12a. The section in the manufacturer's manual describing burner operation should be read thoroughly before attempting to ignite the burner.</p>	

EFFLUENT MONITORING PROCEDURE:

Determination of Lead by Atomic Absorption Using
the Extraction Procedure

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
H. Instrument Operation (continued)	13. Adjust fuel and oxidant pressures.	13a. For the lead procedures it is recommended that the acetylene pressure be set at 4.5 psig and the air pressure be set at 7.0 psig. This will produce a stoichiometric flame that will produce best results. 13b. The analysis of lead is exceptionally sensitive to turbulence and absorption bands in the flame, particularly for the 283.3 nm band. Therefore, some care must be taken to position the light beam in the most stable, center portion of the flame. To do this, first adjust the burner to maximize the absorbance reading with a lead standard. Then, aspirate a water blank and make minute adjustments in the burner alignment to minimize the signal. 13c. See Figure 3 for location of controls.	
	14. Set scale expansion at 2.5.	14a. Graduated in relative absorbance units, this control allows selection of the instrumental sensitivity. 14b. See Figure 5 for location.	
	15. Set calibration control fully clockwise.	15a. This control is in parallel with the channel "A" Digital display, used with the automatic calibration button to set the appropriate absorbance or concentration level on the display. 15b. See Figure 4 for location.	
	16. Set dampening selector switch at M position.	16a. Selects the settling time of both the digital displays or selects the integration sampling period. The "M" position is equivalent to a ten second time constant in the real-time damping of the readout. 16b. See Figure 5 for location.	

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
H. Instrument Operation (continued)	17. Aspirate blank and depress auto-zero. 18. Set read/hold switch to read. 19. Set 0000 with auto-zero zero control.	17a. Activates the channel "A" zeroing servos, electronically setting I equal to 10 when the blank is aspirated. 17b. See Figure 4 for location. 18a. Activates the digital display. 18b. See Figure 5 for location. 19a. The zero control sets the channel "A" digital display to zero. 19b. See Figure 4 for location.	
I. Calibration	1. Calibrate the instrument by aspirating the 10 mg/l standard. 2. Depress the auto-cal button while the standard is being aspirated. 3. Set the digital display to read 1000. 4. Release the auto-cal button. 5. Final adjustment to the concentration read-out should be done with the calibration control without depressing the auto-cal button.	1a. The IL Model 153 atomic absorption instrument can be calibrated either in absorbance or concentration. The procedure given is for concentration. 1b. It is assumed the steps in sections F and G have been carried out. 2a. See Figure 4 for location. 2b. This activates the calibration mechanism to establish the slope of the absorbance concentration. 3a. Use the calibration control to set the digital display if it does not read 1000. 3b. See Figure 4 for location of controls. 5a. Should different sensitivities be required, the scale expansion switch can provide it.	

EFFLUENT MONITORING PROCEDURE:

Determination of Lead by Atomic Absorption Using the Extraction Procedure

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>I. Calibration (continued)</p>	<p>6. Aspirate the rest of the standards.</p> <p>7. Multiply the resultant concentration by the appropriate dilution factor if the sample has been diluted.</p> <p>8. Aspirate the samples.</p> <p>9. Occasionally re-aspirate the 10.0 mg/l standard.</p>	<p>6a. Record the instrument reading for each standard in Column E on the Table in Procedure F.1a.</p> <p>6b. A standard curve, plotting the meter display vs. the known concentration (column E vs. column C, Table I) can be made to prove if a straight line relationship is being obtained.</p> <p>6c. So long as 200 ml portions are used, the calibration as described here will give correct values for the samples in mg/l. Since the standard used to calibrate the instrument was set to read its concentration in mg/l, the dilution factors needed because of using 200 ml and concentrating the sample to 10 ml are incorporated into the instrument calibration. Consequently, no change in the sample size can be done unless the value obtained from the instrument read-out is multiplied by a dilution factor.</p> <p>7a. Concentration of sample before dilution = $\frac{\text{Concentration found from instrument} \times \text{total mls. after dilution}}{\text{mls. of sample taken for dilution}}$</p> <p>8a. Leave the parameters as established by the aspiration of the standards.</p> <p>9a. This will indicate whether the instrument calibration has remained constant.</p> <p>9b. On re-checking the same calibrating standard, depress the auto-cal button to re-establish the slope of the absorption-concentration relationship.</p>	

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
I. Calibration (continued)	10. Record all analytical data in a laboratory notebook.	10a. Report results to the nearest whole microgram per liter.	
J. Instrument Shut-down	1. If a flame is burning aspirate water for about 15 seconds. 2. Close the acetylene cylinder valve. 3. Close the air cylinder valve. 4. Depress H.C. switch to off. 5. Depress power switch to off.	1a. This will prevent build-up of solids in the capillary. 2a. The flame will automatically extinguish itself, leaving about 9 psig in the acetylene supply line. 5a. Caution: Exercise care in touching the burner head and vent area. These will be hot enough to cause serious burns.	
K. Maintenance	1. Clean the instrument regularly. 2. Insure the drain cup is filled each day prior to ignition.	1a. A regular program of care and maintenance will prolong the life-time and maximize its utility. Such items as filters in gas lines, air intakes, burner compartment, burner, and nebulizer should be cleaned. 2a. See the instrument manufacturer's manual for exact procedures.	143

EFFLUENT MONITORING PROCEDURE:

**Determination of Lead by Atomic Absorption Using
the Extraction Procedure**

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
K. Maintenance (continued)	3. Change the acetylene tank when its pressure falls to 75 psig. 4. Change O rings in burner body when aspirating organic solvents.	3a. This will insure a minimum carry over of any acetone. 4a. The manual will provide instructions.	

TRAINING GUIDE

<u>SECTION</u>	<u>TOPIC</u>
I*	Introduction
II	Educational Concepts - Mathematics
III	Educational Concepts - Science
IV	Educational Concepts - Communications
V*	Field and Laboratory Equipment
VI	Field and Laboratory Reagents
VII*	Field and Laboratory Analysis
VIII	Safety
IX	Records and Reports

*Training guide materials are presented here under the headings marked *.

INTRODUCTION	Section I	
	TRAINING GUIDE NOTE	REFERENCES/RESOURCES
A.2.	<p>The atomic absorption determination of lead should be considered as part of the overall determination of the heavy metals. The procedure includes a standard procedure for the metal which utilizes direct aspiration into the atomic absorption instrument. This procedure, however, is limited as to the lower limit which can be measured. The detection limit for the standard procedure is given as 0.5 mg Pb/l; however, the maximum level permitted in potable water is 0.05 mg/l. Consequently, an additional procedure must be carried out--that is, some type of concentration step.</p> <p>When the concentration of the metal is not sufficiently high to determine directly or when considerable dissolved solids are present in the sample, certain of the metals may be chelated and extracted with organic solvents.</p> <p>There is a general procedure available which utilizes a chelating and extraction procedure at distinct pH values to remove all heavy metals from one sample. The procedure here is a modification of this procedure and is specific for the determination of lead.</p> <p>The easiest concentration procedure would involve an evaporation at a low pH of a large volume of sample and then direct aspiration. This procedure is recommended for waters where little interference is present and having a low concentration of solids. The extraction procedure contained in this EMP should be used if the two previous conditions cannot be met in a sample. By using the extraction technique both solids and interfering ions are removed. To just filter such a sample would remove the suspended portion and not give total results.</p> <p>Atomic absorption spectroscopy is similar to flame emission photometry in that a sample is atomized and aspirated into a flame. In atomic absorption a light beam is directed through the flame into a monochromator and onto a detector that measures the amount of light absorbed by the sample. Absorption is more sensitive in most cases because it depends upon the presence of free unexcited atoms which usually have a higher ratio of existence than excited atoms. Since the wavelength of the light beam is characteristic of only the metal being</p>	

EFFLUENT MONITORING PROCEDURE: Determination of Lead by Atomic Absorption
Using the Extraction Procedure

INTRODUCTION

Section I (Cont'd.)

TRAINING GUIDE NOTE

REFERENCES/RESOURCES

determined the light energy absorbed by the sample in the flame is a measure of the concentration of that metal in the sample. This principle is the basis of atomic absorption spectroscopy.

D.2.

The spectral source of monochromatic light for atomic absorption analyses is provided by a hollow cathode lamp. Each lamp emits only the spectra of the element it is designed for. A different lamp is used for each element with few exceptions. The lamps are enclosed in a glass envelope filled with an inert gas. The lead lamp uses neon gas at a low pressure (1 to 10 mm Hg). Once sufficient voltage is applied across the electrodes within the lamp, the inert gas ionizes and current begins to flow. When this happens, positive gas ions bombard the cathode and heating occurs. As the inner surface of the cathode heats, it sputters and the metal from which it is made vaporizes and fills the cathode volume. Charged gas particles collide with the metal atoms raising their valence electrons to higher energy states. When these excited electrons return to their ground state, they emit light. The spectrum thus emitted contains the same wavelength of light required for absorption of that metal (lead) atom in the flame.

FIELD AND LABORATORY EQUIPMENT

Section V

	TRAINING GUIDE NOTE	REFERENCES/RESOURCES
<p>A.1.</p>	<p>The procedure for washing glassware should include the following steps:</p> <ul style="list-style-type: none"> a) Wash with detergent b) Rinse with tap water c) Wash with 1:1 nitric acid (HNO_3) Prepare by adding equal volume of concentrated nitric acid to water--that is, add 500 ml of acid to 500 ml of distilled water. d) Rinse with tap water e) Wash with 1:1 hydrochloric acid (HCl) Prepare by adding equal volume of concentrated hydrochloric acid (HCl) to water--that is, add 500 ml of acid to 500 ml of distilled water. f) Rinse with tap water g) Rinse with deionized distilled water <p>Chromic acid may be useful to remove organic deposits from glassware; however, the analyst should be cautioned that the glassware must be thoroughly rinsed with water to remove the last traces of chromium and any other metals. Chromic acid should not be used with plastic bottles.</p>	<p>Methods for Chemical Analysis of Water and Wastes, 1974. EPA, EMSL. Page 81.</p> <p>EPA Methods Manual, page 81.</p>
<p>D.3</p>	<p>As acetylene (CH_2CH) is packed dissolved in acetone (CH_3COCH_3), cylinders should be stored only in an upright position. The acetone content of the gas typically depends on the cylinder temperature and pressure. Avoid introducing acetone into the instrument. Should this occur the normal flame obtained will have a slight pink tinge and yield an abnormally high background signal. To reduce acetone carry-over, it is desirable to allow acetylene cylinders to stand undisturbed for at least twenty-four (24) hours before use.</p>	<p>Instrumentation Laboratory, Inc., page 3.</p> <p>Instrumentation Handbook 113 Hartwell Ave. Lexington, MA 02173</p>

FIELD AND LABORATORY ANALYSIS

Section VII

	TRAINING GUIDE NOTE	REFERENCES/RESOURCES
E.1	<p>The "E" section must be carried out to produce a result for the determination of the total metal. If Section "E" is not carried out, the sample should be filtered through a 0.045 micron filter, extracted and the value reported as dissolved metal.</p>	
F.1	<p>In an effort to reduce the number of separatory funnels necessary for this determination, it is possible to prepare the blank and standard 200 ml volumetric flasks. Then allow them remain in the flasks while the sample or samples are carried through steps solubilization and extraction. The standards could then be extracted.</p> <p>Then all, standards, blank, and samples, could be carried through the remaining steps, and brought together at step G.5.2.</p>	

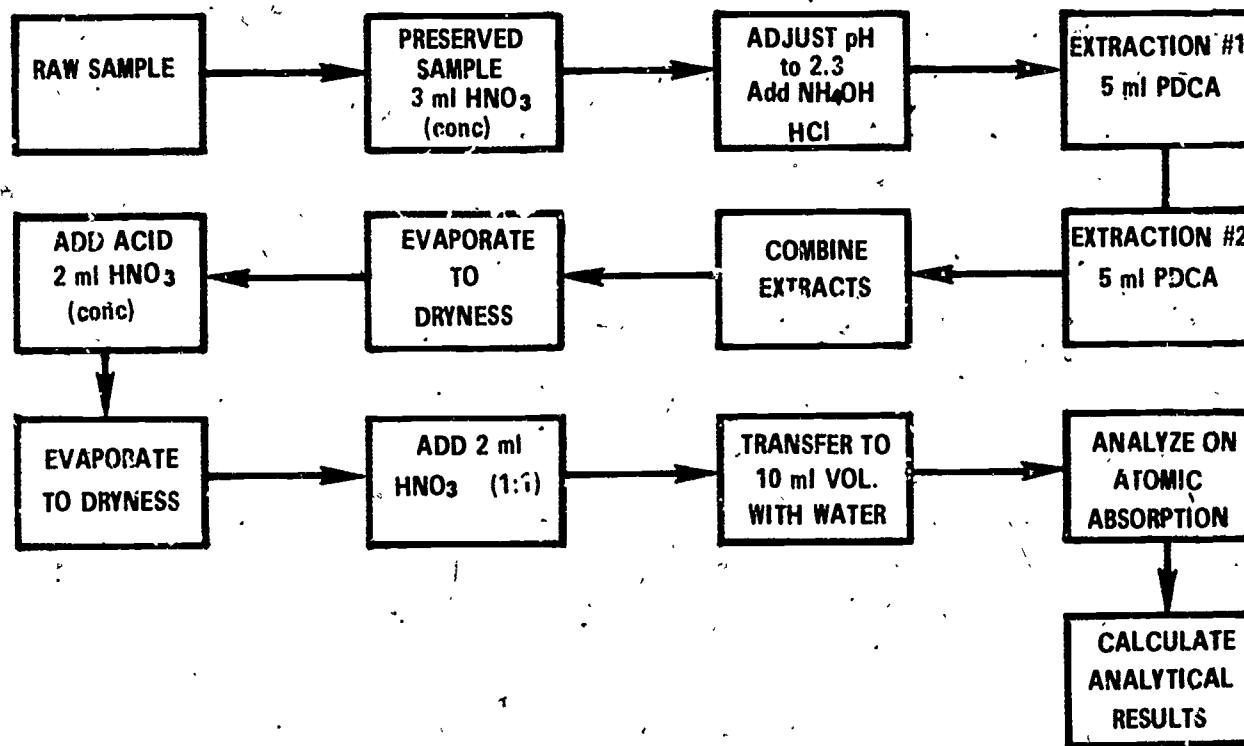


Figure 1. FLOW DIAGRAM FOR THE EXTRACTION METHOD

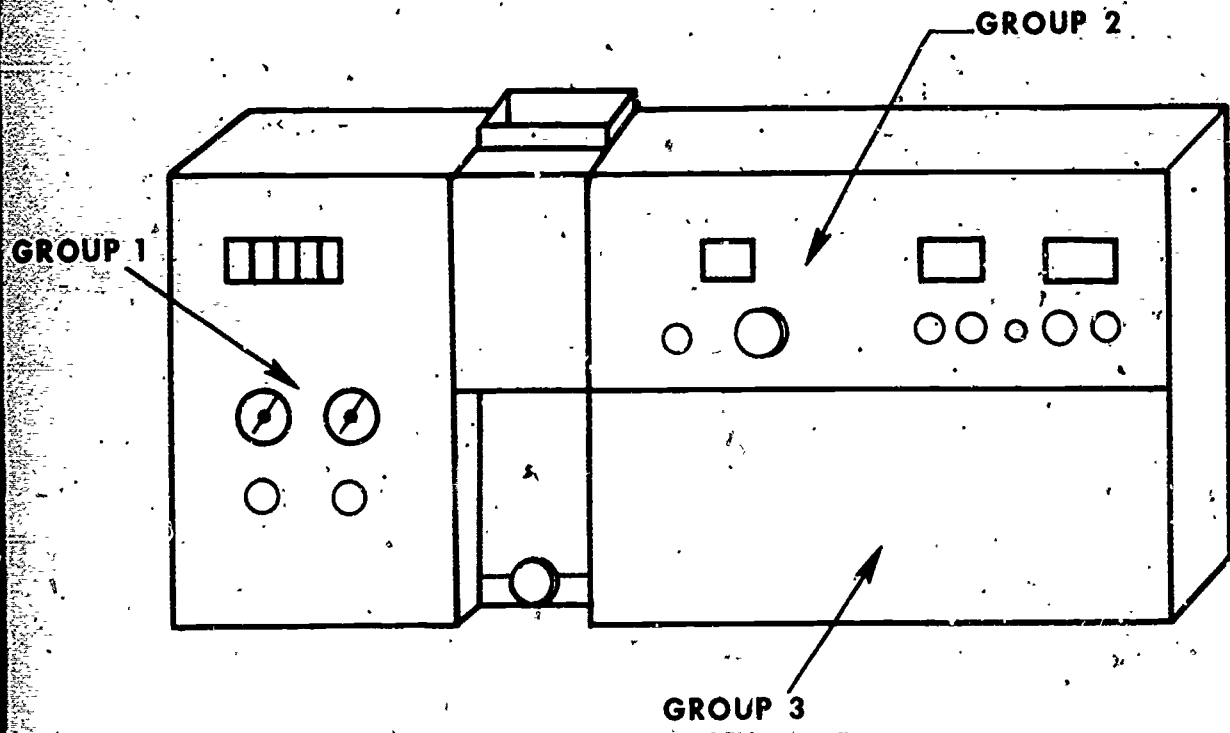


Figure 2. OPERATING CONTROL GROUPS

152

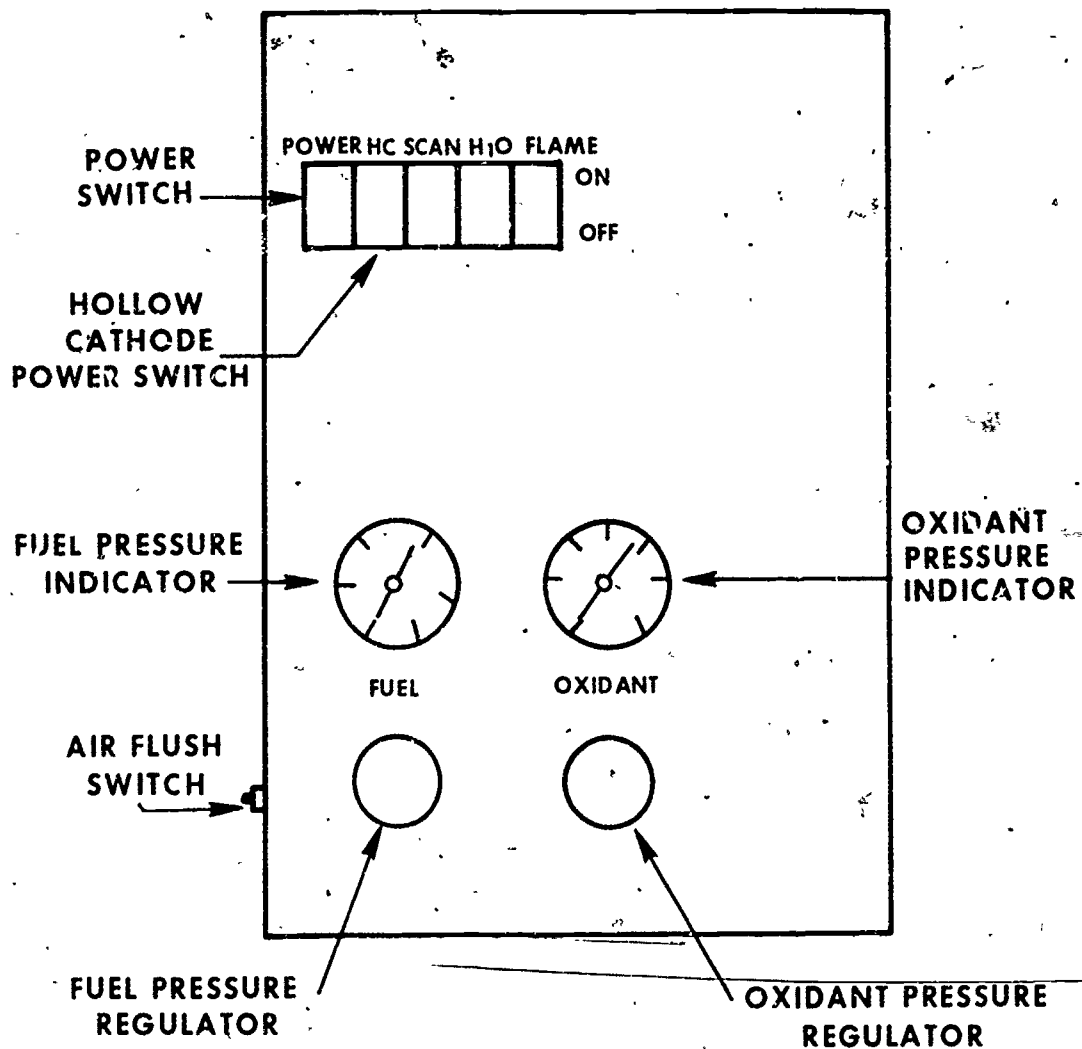


Figure 3. LOCATION OF GROI #1 CONTROLS

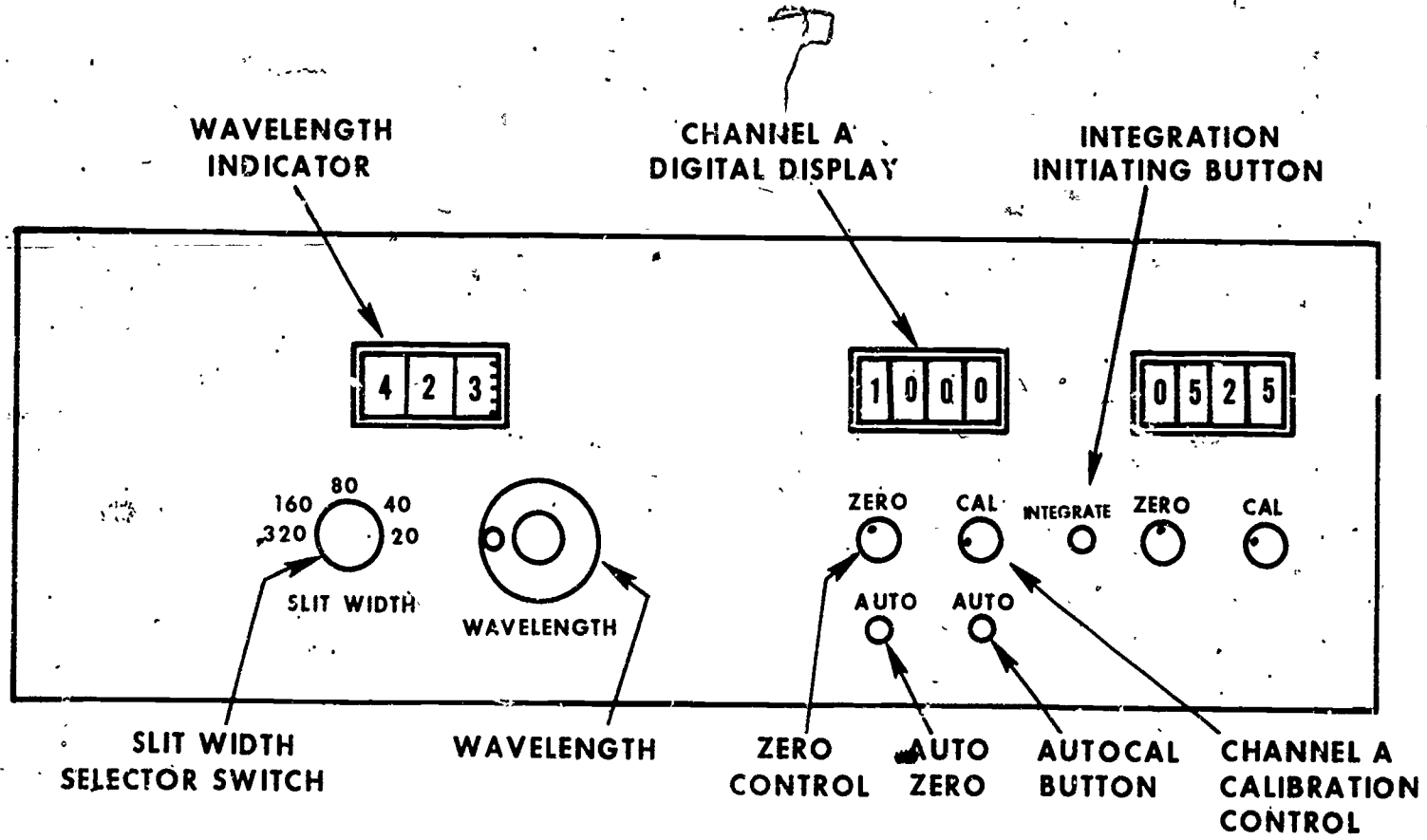


Figure 4. LOCATION OF GROUP 2 CONTROLS

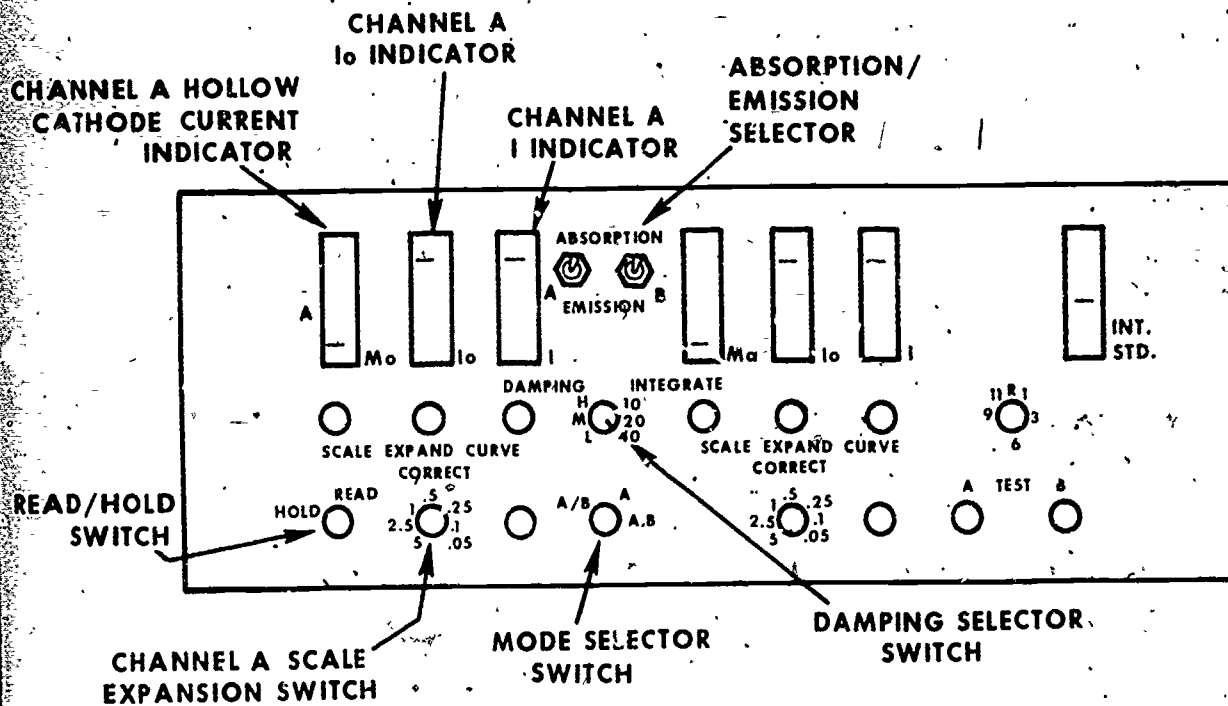


Figure 5. LOCATION OF GROUP 3 CONTROLS

A PROTOTYPE FOR DEVELOPMENT OF
ROUTINE OPERATIONAL PROCEDURES

for the

DETERMINATION OF MERKURY USING THE FLAMELESS
ATOMIC ABSORPTION (COLD VAPOR) TECHNIQUE

as applied in

WASTEWATER TREATMENT FACILITIES
and in the
MONITORING OF EFFLUENT WASTEWATERS

Developed by the

National Training and Operational Technology Center
Municipal Operations and Training Division
Office of Water Program Operations
U.S. Environmental Protection Agency

EFFLUENT MONITORING PROCEDURE: Determination of Mercury Using the Flameless Atomic Absorption (Cold Vapor) Technique

This operational procedure was developed by:

NAME John D. Pfaff

ADDRESS EPA, OWPO, NTOTC, Cincinnati, Ohio 45268

POSITION Chemist-Instructor

EDUCATION AND TECHNICAL BACKGROUND

B.S. Chemistry.

3 years - Research Chemist

15 years - Training Instructor

EFFLUENT MONITORING PROCEDURE: Determination of Mercury Using the Flameless Atomic Absorption (Cold Vapor) Technique

1. Analysis Objectives:

The learner will use the attached EMP to place the Coleman Model MAS-50 Mercury Analyzer System into operation including calibration, reagent and sample preparation, and use of the instrument to determine the concentration of mercury in a sample.

2. Brief Description of Analysis:

This procedure is a three step procedure which 1) chemically vaporizes the sample, 2) introduces the mercury and 3) determines the mercury by flameless atomic absorption techniques.

EFFLUENT MONITORING PROCEDURE: Determination of Mercury Using the Flameless Atomic Absorption (Cold Vapor) Technique

Operating Procedures:

- A. Equipment Preparation
- B. Instrument Set-up
- C. Reagent Preparation
- D. Sample Handling and Preservation
- E. Solubilization of Sample
- F. Calibration
- G. Sample Determination
- H. Calculation

EFFLUENT MONITORING PROCEDURE: Determination of Mercury Using the Flameless Atomic Absorption (Cold Vapor) Technique

General Description of Equipment Used in the Process

A. Capital Equipment

1. Flameless Mercury Analyzer System - Coleman MAS-50*
The Coleman MAS-50 is a self contained unit designed to analyze mercury in water and other environments. It is line operated and complete manufacturer's specifications are as follows:

Sensitivity: 0.01 μg
Saturation: 9 μg
Weight: 17 lbs.

Reproducibility: less than 3%
Power Requirements: 110/220 volts

2. Analytical balance, 200 gram capacity
3. Trip balance, 500 gram capacity
4. Water bath, capable of maintaining 95°C temperature
5. Recorder (optional) - any multi-range, variable speed recorder that is compatible with the system
6. pH meter and electrodes

B. Reusable Supplies

1. Twelve BOD bottles (one bottle is needed per sample)
2. Volumetric flasks
Two 1000 ml
Four 100 ml
One 250 ml
3. Pipets
Three 10 ml volumetric
Three 10 ml graduated
Three 5 ml graduated
Two 1 ml graduated
4. One 100 ml graduated cylinder; one 25 ml graduated cylinder
5. One laboratory apron or coat
6. One pair safety glasses
7. One spatula
8. One pipet bulb
9. One wash bottle for distilled water
10. One glass stirring rod (about 6 inches long)
11. One powder funnel
12. Rubber stoppers - two size #2 (for drying tube)
13. Fifteen ft. Tygon tubing
14. One glass tubing - 6 inches x 3/4 inch diameter
15. One Rotometer (any unit capable of measuring air flow of 1 liter/min.)
16. One set cork hole borers
17. One brush (for cleaning balance)
18. One watch glass for each sample and standard
19. One 150 ml beaker for each sample

*Mention of a specific brand name does not constitute endorsement by the U.S. Environmental Protection Agency.

EFFLUENT MONITORING PROCEDURE: Determination of Mercury Using the Flameless Atomic Absorption (Cold Vapor) Technique

The following equipment is needed depending on which method is chosen to trap the mercury.

a. Liquid Trap

1. Straight glass frit, coarse porosity, such as Corning #404260
2. Filtering flask, such as Corning #40058
3. Rubber stopper, one hole to accept frit
4. Reagents, KMnO_4 and H_2SO_4

b. Solid Trap

1. Activated carbon such as Barnebey and Cheney #580-13 or #580-22
from: Barnebey and Cheney
E. 8th Avenue & Cassidy St.
Columbus, OH 43219

or

Coleman Instruments
42 Madison St.
Maywood, IL 60153
Item #50-160

2. Glassware: can be assembled similar to the drying tube (B-3) or it can be purchased as catalog no. 50-807 from Coleman Instrument Co. (will include adsorbent)

c. Closed System

The following equipment is needed when using the closed system with a trap.

1. Two position valve, or stopcock, such as Corning #442838
2. Glass "Y" shaped tubing connector
3. Pinch clamp, type used for stopping flow in tubing

C. Consumable

1. Sulfuric acid (H_2SO_4) concentrated
2. Nitric acid (HNO_3) concentrated
3. Potassium Permanganate, KMnO_4
4. Potassium Persulfate, $\text{K}_2\text{S}_2\text{O}_8$
5. Sodium Chloride, NaCl

EFFLUENT MONITORING PROCEDURE: Determination of Mercury Using the Flameless Atomic Absorption (Cold Vapor) Technique

C. Consumable (Cont'd.)

6. Hydroxylamine Sulfate $(\text{HONH}_2)_2 \cdot \text{H}_2\text{SO}_4$ or Hydroxylamine Hydrochloride $\text{NH}_2\text{OH} \cdot \text{HCl}$
7. Stannous Sulfate, SnSO_4 or Stannous Chloride, SnCl_2
8. Mercuric Chloride, HgCl_2
9. Potassium dichromate, $\text{K}_2\text{Cr}_2\text{O}_7$ (for cleaning glassware)
10. Magnesium Perchlorate, $\text{Mg}(\text{ClO}_4)_2$ for drying tube, 20 g.
11. Distilled water
12. Sponges (for cleaning laboratory table tops)
13. Notebook for recording weights and readings
14. Graph paper, arithmetic (for plotting standard curve)
15. Two pieces of glass tubing (5mm diameter, about two inches long) for the drying tube
16. Glass wool (for drying tube)
17. Plastic weighing boats
18. Pen or pencil

EFFLUENT MONITORING PROCEDURE: Determination of Mercury Using the Flameless Atomic Absorption (Cold Vapor) Technique

Sample

Solubilization

Chemical Sample Preparation.

- a. Oxidation of all mercury to mercuric form
- b. Reduction of all mercuric mercury to metallic mercury

Aeration

The metallic mercury is circulated as a vapor through the system

Flameless Atomic Absorption

Absorption of energy at 253.7nm from a hollow cathode lamp measurement by a photodetector

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>A. Equipment Preparation</p> <p>1. Cleaning of Glassware</p> <p>2. Balance Preparation</p>	<p>1. Wash with detergent</p> <p>2. Rinse with tap water</p> <p>3. Rinse with 1:1 Nitric Acid</p> <p>4. Rinse with tap water</p> <p>5. Rinse with 1:1 Hydrochloric acid</p> <p>6. Rinse with tap water</p> <p>7. Rinse with distilled water</p> <p>1. Check all balances for cleanliness and proper operation.</p>	<p>1a. Cleaning should be carried out in this order.</p> <p>1b. Care should be taken to insure clean glassware as mercury is a common contaminant. All glassware should be kept covered after cleaning.</p> <p>3a. Add 500 ml conc. nitric acid (HNO_3) to 500 ml distilled water.</p> <p>5a. Add 500 ml conc. hydrochloric acid (HCl) to 500 ml distilled water.</p>	<p>V.A.1.1</p>
<p>B. Instrumental Set-up</p> <p>1. Mercury Trap - Liquid Type</p>	<p>1. Before operation of the instrument, four additions to the system should be considered (Fig. 1).</p> <p>2. One of the following mercury traps should be included in the system.</p>	<p>1a. There are two ways the flow system can be set up. It can be operated as a closed or open system. In the closed system the mercury vapor continuously passes through the system until wasted in the mercury trap by the operator. In the open system the vapor passes through the absorption tube only once and goes directly to the trap. Which system is chosen will dictate what equipment is necessary. Figure 1 shows the choices and the equipment necessary for each.</p> <p>2a. Because of the toxic nature of mercury vapor precaution must be taken to avoid contamination. The vapor will be held in the trap after it has been measured.</p>	

EFFLUENT MONITORING PROCEDURE: Determination of Mercury Using the Flameless Atomic Absorption (Cold Vapor) Technique

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>1. Mercury Trap - Liquid Type (Cont'd)</p>	<p>3. For a liquid type trap use a 250 ml side arm filtering flask.</p> <p>4. Assemble as shown in Figure 2.</p> <p>5. Insert straight gas dispersion tube or frit through the hole so that the bottom or fritted end is about one inch above bottom of the flask.</p> <p>6. Insert into filtering flask.</p> <p>7. Connect tygon tubing to top end of frit and a second piece of tygon tubing to the side arm of filtering flask.</p> <p>8. Add 200 ml of 1:1 potassium permanganate ($KMnO_4$) -- sulfuric acid (H_2SO_4) Reagent (Reagent #16)</p>	<p>3a. Use a filtering flask such as Corning #400580 or its equivalent.</p> <p>4a. Use a #3 cork hole borer to make the hole.</p> <p>5a. Frit should have a coarse porosity such as Corning #404260 or equivalent. The frit should always fall below liquid level in the flask. Should the level become low add more liquid (Reagent #10). The nonfritted end should be lubricated and care taken when the frit is inserted through the stopper so as not to break the frit and injure the worker.</p> <p>7a. Care should be taken so that the liquid level does not come close to the opening of the side arm of the flask. This could flood the instrument if allowed to do so. If flooding should occur, disassemble the absorption tube and clean it and the tubing immediately.</p> <p>8a. A solution of 0.25% iodine in a 3% potassium iodide (KI) solution may also be used.</p> <p>8b. Filling the flask can be postponed until all of the apparatus is assembled.</p>	<p>VIII.</p>
<p>2. Mercury Trap - Solid Type</p>	<p>1. The apparatus can be prepared similar to the drying tube (B-3) but packed with 2-3 grams of activated carbon.</p>	<p>1a. Locate after 2 position valve in closed system, figure 1 (system two) or after the analyzer in an open system, figure 1 (system three).</p>	

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>2. Mercury Trap - Solid Type (Cont'd.)</p>	<p>2. The equipment can be purchased with adsorbent as an option from the analyzer manufacturer.</p>	<p>2a. Position as above.</p>	
<p>3. Drying Tube</p>	<p>1. Construct as shown in Figure 3.</p> <p>2. Bore a hole through a number 2 stopper with a number 2 cork hole borer. Repeat with a second stopper.</p> <p>3. Insert a 2 inch long piece of glass tubing (5 mm diameter) through each stopper allowing about 1/2 inch protruding from each end.</p> <p>4. Fill a 6 inch piece of 3/4 inch diameter tubing with 20 grams of magnesium perchlorate ($Mg(ClO_4)_2$).</p> <p>5. Use a small piece of glass wool in each end of the tube to prevent loss of granules.</p> <p>6. Insert stopper prepared above in each end of tube.</p> <p>7. Replace drying agent when needed.</p>	<p>1a. Place between sample container and instrument.</p> <p>3a. Care should be taken when inserting glass tubing.</p> <p>4a. Other drying agents such as calcium chloride ($CaCl_2$) may be used.</p> <p>5a. The tube should not be packed so tight as to restrict gas flow.</p> <p>7a. Replace magnesium perchlorate or any drying agent regularly. These materials tend to cake and form a plug when their limit of saturation is approached. The length of time the material will</p>	



EFFLUENT MONITORING PROCEDURE: Determination of Mercury Using the Flameless Atomic Absorption (Cold Vapor) Technique

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>3. Drying Tube (Cont'd.)</p> <p>4. Rotometer</p>	<p>1. Must be capable of measuring a gas flow of 1 liter per minute.</p> <p>2. Connect one length of tubing between the sample container and the drying tube, then out of the drying tube through the rotometer to the fitting on the side of the instrument marked "in".</p> <p>3. A second length of tubing should begin at the fitting of the instrument marked "out" and proceed to the next piece of equipment.</p>	<p>last will vary with use and samples. Experience will dictate a routine.</p> <p>1a. Place between water trap and instrument. See Figure 1 for location.</p> <p>1b. The rotometer may be removed from the circuit after the instrument pump rate is checked.</p> <p>1c. The flow rate should be checked periodically to insure flow rate has not changed.</p> <p>2a. The connection must be made to the sample container by side arm. Reverse tubing connections may flood the instrument with liquid.</p> <p>3a. See Figure 1 for gas flow path.</p>	
<p>5. Two Position Valve</p>	<p>1. A two position valve is necessary when using a closed system and a trap. Use stopcock, Corning No. 442838 or equivalent for the two way valve or stopcock.</p> <p>2. One position of the valve should go through the trap to the sample container. The other position should by-pass the trap.</p>	<p>2a. It is important to maintain a specific air volume in the system. Once the system is calibrated this volume cannot be changed unless the system is recalibrated.</p>	

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>5. Two Position Valve (Cont'd.)</p> <p>5. MAS-50</p>	<p>and be connected to the frit of the sample container.</p> <ol style="list-style-type: none"> 1. Place the instrument on a flat surface where it will be isolated from shock and undue temperature variation. 2. Before turning on the instrument a check should be made of the power requirements. 3. Remove back panel and note the position of the 110-220 volt switch. Place the switch in the position desired. Then close the rear panel. 4. Meter mechanical zero adjustment <ul style="list-style-type: none"> ● with power switch off ● Check to see if meter pointer indicates exactly 0 micrograms. ● Using the mechanical zero adjust screw (Figure 4). Adjust meter until it reads 0 micrograms. 5. Place the On-Off switch in the On position. Allow 15 minute warm-up. 	<ol style="list-style-type: none"> 1a. Do not block or cover the ventilation slots at the rear and base of the instrument. 2a. The instrument requires a grounded power connection. Preferably a three prong receptacle. If not then a grounded two prong set-up. 3a. Should 220 V. position be used an appropriate 220 V plug should be installed in place of the plug supplied with the analyzer. 4a. If meter pointer does indicate zero, no further action is necessary. 	

EFFLUENT MONITORING PROCEDURE: Determination of Mercury Using the Flameless Atomic Absorption (Cold Vapor) Technique

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>6. MAS-50 (Cont'd.)</p> <p>7. Recorder (optional)</p>	<p>6. During warm-up check the connections in the external set-up of the gas flow path as in Figure 1.</p> <p>1. An external recorder may be connected to the instrument to give indications coinciding with the instruments meter.</p> <p>2. Adjust recorder by placing instrument controls as follows:</p> <ul style="list-style-type: none"> ● Meter switch - % T ● Memory - OFF ● Recorder - plugged in ● Shutter - open <p>3. Use recorder zero to align recorder with 100% T.</p> <p>4. Use analyzer's recorder range control to achieve full scale indication on 0% T.</p>	<p>6a. Any leaks in the system will cause low readings and allow hazardous mercury vapors to escape.</p> <p>1a. Any recorder having 50 K ohm impedance or more - the output of the instrument is variable from 0 to 300 mV full scale. (Such as a Coleman Model 56 or 165 chart recorder)</p> <p>4a. Shutter closed for this adjustment.</p>	
<p>C. Reagent Preparation</p> <p>1. Sulfuric Acid 0.5 N</p>	<p>1. Add 14.0 ml conc. sulfuric acid (H_2SO_4) to approximately 500 ml water and mix. Then dilute with water to 1 liter volume.</p>	<p>1a. The concentrated H_2SO_4 should be of low mercury concentration.</p> <p>1b. Unless specified the term water means distilled water.</p>	

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
2. Potassium Permanganate Soln. 5% solution w/v	1. Prepare 100 ml of solution containing 5.0 grams potassium permanganate (KMnO ₄)	1a. Should a larger amount of reagent solution be needed the same ratio should be maintained. For example: prepare 1000 ml of solution containing 50 grams KMnO ₄ .	
3. Potassium Persulfate Soln. 5% solution w/v	1. Dissolve 5.0 g potassium persulfate (K ₂ S ₂ O ₈) in water and dilute to 100 ml.		
4. Sodium Chloride - Hydroxylamine Sulfate Soln. Solution 12% NaCl 12% ((HONH ₂) ₂ · H ₂ SO ₄) w/v	1. Dissolve 12.0 g of sodium chloride (NaCl) and 12.0 g of hydroxylamine sulfate ((HONH ₂) ₂ · H ₂ SO ₄) in water and dilute to 100 ml.	1a. Hydroxylamine hydrochloride (NH ₂ OH · HCl) may also be used. It should be prepared in the same manner.	
5. Stannous Sulfate soln. 10% solution w/v	1. Dissolve 25.0 g stannous sulfate (SnSO ₄) in 0.5 N sulfuric acid and dilute with 0.5 sulfuric acid to 250 ml.	1a. Stannous chloride (SnCl ₂ · 2H ₂ O) may be used and be prepared in the same manner. 1b. The acid is reagent no. 1. 1c. This a suspension and should be stirred continuously during use.	
6. Sulfuric Acid Concentrated (H ₂ SO ₄)	1. No preparation necessary.	1a. This should be reagent grade and low in mercury concentration. 1b. Caution: this is corrosive.	
7. Nitric Acid Concentrated (HNO ₃)	1. No preparation necessary.	1a. This should be reagent grade and low in mercury content. 1b. Caution: this is corrosive. 1c. If a high reagent blank is obtained, it may be necessary to distill the nitric acid.	
8. Stock Mercury Soln. (HgCl ₂) (Cont'd.)	1. Dissolve 0.1354 g of mercuric chloride (HgCl ₂) in water.	1a. Caution: solution will increase in temperature.	



EFFLUENT MONITORING PROCEDURE: Determination of Mercury Using the Flameless Atomic Absorption (Cold Vapor) technique

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
8. Stock Mercury Soln. (HgCl ₂) (Cont'd.)	2. Add 10 ml conc. nitric acid (HNO ₃) 3. Cool to room temperature. 4. Dilute to 100 ml with water.	2a. Concentration of stock solution is now 1 ml = 1 mg Hg	
9. Intermediate Mercury Solution - (HgCl ₂) Dilution of Solution 8	1. This solution is a dilution of the stock solution to adjust the conc. of Hg to 10.0 µg/ml. Proceed as follows. 2. Add about 700 ml water to a 1000 ml vol. flask. 3. Add 0.5 ml conc. HNO ₃ 4. Add 10 ml stock Hg soln. 5. Dilute to 1000 ml mark. This solution contains 10 µg/ml.	1a. Prepare fresh before use. 3a. The nitric acid conc. of the dilutions including the working solution should be maintained at 0.15%. This acid should be added to the flask before addition of the aliquot.	
10. Working Mercury Solution - (HgCl ₂) Dilution of Solution 9	1. Add about 700 ml water to a 1000 ml vol. flask. 2. Add 1.5 ml conc. HNO ₃ . 3. Add 10 ml of intermediate solution (10 µg/ml). 4. Dilute to 1000 ml mark. This is working solution and contains 0.1 µg/ml.	1a. Prepare fresh before use.	

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS.	TRAINING GUIDE NOTES
11. Potassium Permanganate 0.1 N (KMnO_4) and Sulfuric Acid 10% Solution (For Mercury Trap)	<ol style="list-style-type: none"> 1. Dissolve .316 g potassium permanganate (KMnO_4) in 100 ml water. 2. Add 10 ml conc. sulfuric acid (H_2SO_4) to about 80 ml water. Dilute to 100 ml with water. 3. Mix equal volumes of each solution; KMnO_4 (1) and H_2SO_4 (2). 	<ol style="list-style-type: none"> 1a. Let stand until following solution is prepared. 2a. Caution: heat generated. 2b. Should be at room temperature before volume adjustment. 	
12. Nitric Acid (1 + 1 dilution) (HNO_3)	<ol style="list-style-type: none"> 1. Prepare by adding an equal volume of acid to an equal volume of water (ie. 50 ml HNO_3 to 50 ml water). 	<ol style="list-style-type: none"> 1a. Caution: acid should always be added to water to avoid spattering. 1b. Caution: Heat is generated. 1c. Use safety glasses 	
13. Hydrochloric Acid (1 + 1 dilution HCl)	<ol style="list-style-type: none"> 1. Prepare by adding an equal volume of acid to an equal volume of water (ie. 50 ml HCl to 50 ml water). 	<ol style="list-style-type: none"> 1a. See cautions above. 1b. Use safety glasses. 	
D. Sample Handling and Preservation	<ol style="list-style-type: none"> 1. Upon collection the sample pH should be lowered to 2 or lower by the addition of concentrated nitric acid. 	<ol style="list-style-type: none"> 1a. If only dissolved mercury is to be determined the sample should be filtered before addition of the acid. For total mercury the filtration is omitted. 	

EFFLUENT MONITORING PROCEDURE:

Determination of Mercury Using The Flameless Atomic Absorption (Cold Vapor) Technique

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
E. Solubilization of Sample	<ol style="list-style-type: none">1. Acidify the sample with 1:1 nitric acid to a pH of 2 at the time of sample collection.2. Transfer 100 ml of well mixed sample to a 150 ml beaker.3. Add 3 ml of concentrated nitric acid to the sample.4. Place the beaker on a hot plate and cautiously evaporate to dryness.5. Cool the beaker and add another 3 ml portion of the concentrated nitric acid.6. Cover the beaker with a watch glass and return to the hot plate.7. Increase the temperature of the hot plate so that a gentle reflux action occurs.8. Continue heating until a light colored residue forms.	<ol style="list-style-type: none">2a. 100 ml is the usual sample volume. If more metals are to be determined, a larger sample volume must be used.3a. Use a 10 ml graduated pipet.4a. The sample should not boil during this time.5a. Use a 10 ml graduated pipet.7a. This will be indicated by droplets forming on the underside of the watch glass.8a. This will indicate that the digestion is complete.	

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
E. Solubilization of Sample (Continued)	9. Add sufficient [1:1 hydrochloric acid (HCl) solution] and again warm the beaker to dissolve the residue. 10. Wash down the beaker walls and watch glass with distilled water. 11. Adjust the volume to 100 ml (or the starting volume size).	9a. Use a 1 ml graduated pipet. 11a. The sample is now ready for analysis.	
F. Calibration	1. Turn on power and pump switches. 2. Set range switch on %T. 3. Memory switch off. 4. Shutter closed. 5. Adjust to 0%T with 0%T adjustment knob.	1a. Refer to figure 4 and 5. 1b. Allow 15 minutes warm-up time. 1c. Calibration is necessary initially and thereafter only if mercury lamp or optical filter is replaced or if readings with standard samples constantly deviate more than 5% of full-scale in one direction. 2a. Refer to Figure 5 for location of various switches.	

185

186

EFFLUENT MONITORING PROCEDURE: Determination of Mercury Using the Flameless Atomic Absorption (Cold Vapor) Technique

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
F. Calibration (Cont'd.)	<p>6. Open shutter.</p> <p>7. Adjust 100% T with 100% T adjustment knob.</p> <p>8. Repeat steps 4 through 7 until 0% T adjustment does not change full-scale (100% T) setting.</p> <p>9. Prepare four samples of 100 ml volume containing 1.0 µg of mercury.</p> <p>10. Place the meter switch on % T and open shutter, memory switch off, check 0% T and 100% T settings.</p> <p>11. Determine as if these four standards were samples. Record the % T value for each standard sample.</p> <p>12. Adjust meter switch to X1.</p> <p>13. Adjust the 100% knob so that meter indicates value of mercury standard.</p> <p>14. Set meter switch to % T.</p>	<p>9a. Use the working solution (Reagent #10) (0.1 µg/ml) and dilute 10 ml of that solution to 100 ml with distilled water. This 100 ml volume will now contain 10 µg Hg.</p> <p>10a. This step should be performed just after placing the aerator in the sample BOD bottle.</p> <p>11a. Use steps 18-32 of Section F - Sample Determination.</p> <p>11b. When the sample contains only distilled water and an inorganic mercury (i.e., mercuric chloride) the heating step (steps 23 and 24 of sample determination) can be omitted.</p> <p>12a. This step should only be done after the needle of the meter has returned to 100% T.</p> <p>13a. The meter should be adjusted on the X1 setting to read 1 mg on the X1 scale.</p>	

187

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>F. Calibration (Cont'd.)</p>	<p>15. Using the calibration adjust % T to average value found in step 11.</p> <p>16. Switch meter switch between X1 and % T settings.</p> <p>17. Reset 0% T and 100% T with meter switch at % T and memory off.</p> <p>18. Transfer 0, 0.5, 1.0, 2.0, 5.0, and 10 ml portions of the working mercury solution (sol. 10) to a series of 300 ml BOD bottles.</p> <p>19. Add enough water to each bottle to make a total volume of 100 ml.</p> <p>20. Treat as samples using the procedure under sample determination.</p> <p>21. Construct a standard curve.</p>	<p>15a. Located on the back of the analyzer. This is a screwdriver adjustment.</p> <p>16a. The two scales should agree on the location of the needle. The X1 scale showing 1 μg and the % T scale showing 55% T (the corresponding % T value just below 1 μg on the X1 scale).</p> <p>18a. Working solution contains 0.1 μg per ml. The series will contain from 0 to 1.0 μg of mercury.</p> <p>19a. The 0 μg or blank is a check on the purity of the reagents.</p> <p>21a. Use arithmetic graph paper and plot maximum meter reading versus micrograms of mercury in standard.</p>	<p>II.E.21a</p>
<p>G. Sample Determination Operation</p>	<p>1. Plug the line cord into power receptacle.</p>	<p>1a. This procedure provides start-to-shutdown instructions. It is assumed that the analyzer is properly installed (Instrument Set-Up) adjusted, and calibrated (Calibration).</p>	

EFFLUENT MONITORING PROCEDURE: Determination of Mercury Using the Flameless Atomic Absorption (Cold Vapor) Technique

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>G. Sample Determination Operation (Cont'd.)</p>	<ol style="list-style-type: none"> 2. Allow at least 15 minute warm-up. 3. Memory switch OFF. 4. Meter switch to % T. 5. Close shutter. 6. Turn 0% T knob to adjust meter pointer to approximately 0% T. 7. Open shutter. 8. Turn 100% T knob to adjust meter pointer to exactly 100% T. 9. Close shutter. 0. Readjust 0% T. 1. Open shutter. 2. Meter switch to X5. 3. Adjust 100% T. 4. Depress reset button. 5. Set memory switch as desired. 	<p>14a. This step is necessary only on older model instruments.</p> <p>15a. With memory switch on, the meter will hold the maximum value. At the OFF position the analyst will have to watch for the highest value. The memory switch must be turned off after each run to cancel the meter reading and allow instrument</p>	

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>G. Sample Determination Operation (Cont'd.)</p>	<p>16. Select desired scale - X1, X5, % T.</p> <p>17. Pour 100 ml of sample into BOD bottle.</p> <p>18. Add 5 ml conc. sulfuric acid (H_2SO_4). Mix.</p> <p>19. Add 2.5 ml conc. nitric acid (HNO_3). Mix.</p> <p>20. Add 15 ml potassium permanganate ($KMnO_4$) - reagent no. 2.</p> <p>21. Allow to stand 15 minutes.</p> <p>22. Add 8 ml potassium persulfate ($K_2S_2O_8$) - reagent no. 3.</p> <p>23. Heat in a water bath at 95°C for 2 hours.</p>	<p>to return to zero. In all cases the memory function does not affect the recorder.</p> <p>16a. Scale X1 - Direct reading 0 to 9 μg of Hg. X5 - Direct reading 0 to 0.28 μg of Hg % T - Energy transmission scale 100 to 0% T.</p> <p>17a. If the sample has more than 9 μg of mercury a smaller portion of sample should be used and diluted to 100 ml. The resulting value should be multiplied by the dilution factor.</p> <p>17b. Use a separate BOD bottle for each sample or standard. The reagents can be added to the bottle and bottles can be heated.</p> <p>18a. Use caution when adding concentrated acids.</p> <p>20a. Shake and add additional portions of potassium permanganate solution, if necessary, until the purple color persists for at least 15 minutes.</p> <p>20b. A small (25 ml) graduated cylinder can be used.</p> <p>21a. This is a standing time and the sample should be allowed to actually stand for the 15 minutes.</p> <p>23a. The heat step is required for methyl mercuric chloride when present in or spiked to a natural system. For standards prepared with distilled water and mercuric chloride, the heating step is not necessary.</p>	<p>VII 17.</p>

EFFLUENT MONITORING PROCEDURE: Determination of Mercury Using the Flameless Atomic Absorption (Cold Vapor) Technique

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>G. Sample Determination Operation (Cont'd.)</p>	24. Cool to room temperature.	24a. Tap water can be used to adjust this temperature.	
	25. Add 6 ml of sodium chloride hydroxylamine sulfate, $(\text{NaCl} \cdot (\text{HO} \cdot \text{NH})_2 \cdot \text{H}_2\text{SO}_4)$ reagent no. 4.	25a. This is done to reduce the excess potassium permanganate.	
	26. Allow to stand at least 30 seconds.	26a. Up to this point all samples to be run could be treated together as a group. From this point each must be done individually as the mercury is generated and must be measured immediately.	
	27. Add 5 ml of stannous sulfate (SnSO_4) solution. reagent no. 5.		
	28. Immediately insert bubbler into the bottle.	28a. The shutter should be open and the valve in the position to by-pass the trap.	
	29. Record the highest value indicated by the meter pointer.	<p>29a. The meter will increase and reach a maximum within 1 minute.</p> <p>29b. With the memory switch on, the meter will automatically retain the highest value reached.</p> <p>29c. With the memory switch off, either a recorder must be used or the meter watched continuously.</p>	
	30. In the closed system, turn one stopcock valve to the trap position and allow pump to continue operating.	30a. Once this has been done the mercury will be absorbed and the sample cannot be reused.	
	31. Position memory to OFF if it was used.	<p>31a. On new instruments this will clear memory. On older instruments the reset button should be depressed.</p> <p>31b. Before the next sample can be started the meter must be returned to the 0 μg position.</p>	

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>G. Sample Determination Operation (Cont'd.)</p>	<p>32. Turn stopcock valve to the sample position.</p> <p>33. Remove frit from sample container.</p> <p>34. Determine all other samples by repeating steps 7 through 31.</p> <p>35. When last sample has run, shut down analyzer.</p>	<p>33a. Pump should be allowed to continue operating.</p> <p>35a. To shut down, keep valve to trap until meter reads 0%, and then remove aerator from BOD bottle. It is recommended that the analyzer be left ON during the day if routine sampling is to continue.</p>	
<p>H. Calculations</p>	<p>1. Use values determined from the standards.</p> <p>2. Plot the meter reading (μg) versus the known concentration of the standards (μg).</p> <p>3. Calculate the mercury concentration in the sample</p>	<p>1a. Step 20 under Calibration.</p> <p>2a. This will show if the meter is in calibration. If all readings agree, i.e., if for example, a known standard of 1.0 μg concentration reads 1.0 μg on the X1 scale, the unknown sample concentrations can be read directly in μg. Otherwise the standard curve should be used to obtain the correct value.</p> <p>3a. Use the equation:</p> $\mu\text{g Hg}^*/\text{l} = \left(\frac{\mu\text{g reading}}{\text{from meter}} \right) \times \left(\frac{1000}{\text{ml's of sample used}} \right)$ <p>*Hg = Mercury</p>	

EFFLUENT MONITORING PROCEDURE: Determination of Mercury Using the Flameless Atomic Absorption (Cold Vapor) Technique

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
H. Calculations (Cont'd.)	4. Report mercury conc. as follows: Below 0.2 µg/l as <0.2 µg/l Between 1.0 and 10.0 µg/l using one place after decimal Above 10.0 µg/l using only whole numbers	4a. <= less than.	

EFFLUENT MONITORING PROCEDURE: Determination of Mercury Using the Flameless Atomic Absorption (Cold Vapor) Technique

Section I

TRAINING GUIDE, NOTE

REFERENCES/RESOURCES

Theoretical Concepts

The method is based on the method developed by Hatch and Ott. The Mercury in the sample is oxidized to the mercuric ion with potassium permanganate in a nitric-sulfuric acid medium. Hydroxylamine sulfate is then added to remove the excess permanganate. Stannous sulfate, is then added to reduce the mercury to metallic form. Then the mercury is vaporized and circulated by the bubbler system. This consists of a circulating pump and the bubbler. Measurement is made with a flameless atomic absorption spectrophotometer. The energy at the 253.7 nm mercury line emitted by a mercury lamp is absorbed by the mercury vapor in the flow-through absorption cell. The change in transmittance is detected by a phototube.

Analytical Chemistry,
Volume 40, pg. 2085
December 1968

Instruction Manual
Coleman Model MAS-50
Mercury Analyzer
Coleman Instruments
Maywood, IL

EFFLUENT MONITORING PROCEDURE: Determination of Mercury Using the Flameless Atomic Absorption (Cold Vapor) Technique

Section II

TRAINING GUIDE NOTE	REFERENCES/RESOURCES
<p>E.21.A</p> <p>The standard curve is a reference to a fundamental law of absorption chemistry known as the Beer-Lambert law. Simply, this law states that the amount of energy absorbed by a solution is proportional to the concentration of the absorbing material in the solution. Applied to this outline the amount of energy absorbed at the wavelength of 253.7 nm is proportional to the amount of mercury present in a solution.</p> <p>If a series of known solutions are prepared as in step 18 under Calibration and the meter reading, i.e. the maximum reading on either the % T or the XI scale are plotted, a straight line should result.</p> <p>When an unknown sample value is obtained its mercury content can be determined from the straight line or standard curve.</p> <p>The two meter setting, i.e., % T and XI will give the same results. Percent T will require a standard curve to convert the % T reading to concentration while the XI scale will convert and express the reading directly into micrograms of mercury.</p> <p>To use the attached graph paper prepare the standards as in step 18 and 19 of the Calibration section and run them as samples. The known concentrations are plotted along the bottom of the graph. Then plot the values obtained from the meter, XI is plotted on the left of the graph and % T is plotted on the right.</p> <p>Where the known concentration line intersects with the appropriate meter reading a mark is made. After all six standards are plotted, draw a line through the marks.</p>	

202

EFFLUENT MONITORING PROCEDURE: Determination of Mercury Using the Flameless Atomic Absorption (Cold Vapor) Technique

Section V

TRAINING GUIDE NOTE

REFERENCES/RESOURCES

A.1.1

The use of chromic acid is a recommended procedure.

1. Pour 35 ml of distilled water in a 250 ml beaker.
2. Add about 1/8 teaspoon (simply estimate this quantity) of sodium dichromate, $\text{Na}_2\text{Cr}_2\text{O}_7$, to the water.
3. Swirl the beaker until the sodium dichromate has dissolved.
4. Keep repeating steps 2 and 3 until no more sodium dichromate will dissolve.
5. Pour the solution into a 2 liter beaker.
6. Slowly pour 1 liter of concentrated sulfuric acid, H_2SO_4 , into the 2 liter beaker.
Caution: Use eyeglasses and protective clothing.
7. Stir the mixture thoroughly.
8. Store it in a glass stoppered bottle.
9. The cleaning solution should be at a temperature of about 50°C when it is used.
10. It may therefore be necessary to warm the cleaning solution.
11. When using the warm cleaning solution, fill the piece of glassware with the solution.
12. Allow it to soak for 2-3 minutes (or longer).
13. Pour the cleaning solution back into the storage bottle.
14. Rinse the piece of glassware ten times with tap water.
15. The cleaning solution may be reused until it turns green.
16. It should then be discarded.

13th Standard Methods,
p. 135, section 2.c.2

EFFLUENT MONITORING PROCEDURE: Determination of Mercury Using the Flameless Atomic Absorption (Cold Vapor) Technique

Section VIII

TRAINING GUIDE NOTE


REFERENCES/RESOURCES

Maintenance Practices

Coleman MAS-50 Operation Manual


1. Immediately clean all spilled materials from the instrument and wipe dry.
2. Do not leave vapors in absorption cell. Remove bubbler from the sample flask and operate pump after passing air through trap for 2 minutes.
3. Whenever the instrument is not in relatively continuous use, turn it off and cover it with the plastic cover. If it is not to be used for several days, the line cord may be disconnected from the power source.
4. Circulating air may eventually cause dust to accumulate on exposed optical parts. Parts, such as the lamp filter and cell windows may be wiped with clean, lintless, non-scratching tissue.
5. If a liquid sample should be drawn into the Analyzer disconnect the tubing at both ends of the absorption tube. Remove the tube, disassemble, rinse with distilled water, dry and reassemble with new windows. Flush the system with distilled water introduced at the cell end of both inlet and outlet tubing. Then gently blow out both lines with air. Re-install the absorption cell.
6. If problems arise in attaining either 0% T or 100% T with the appropriate adjustment controls, the windows in the absorption tube should be checked particularly the one opposite the mercury lamp. If the cell window has become cloudy it should be changed. Care should be taken to keep finger prints off the windows. After exposure to the ultra-violet light the prints cannot be removed. To replace follow the manufacturer's directions.

X1 Readings

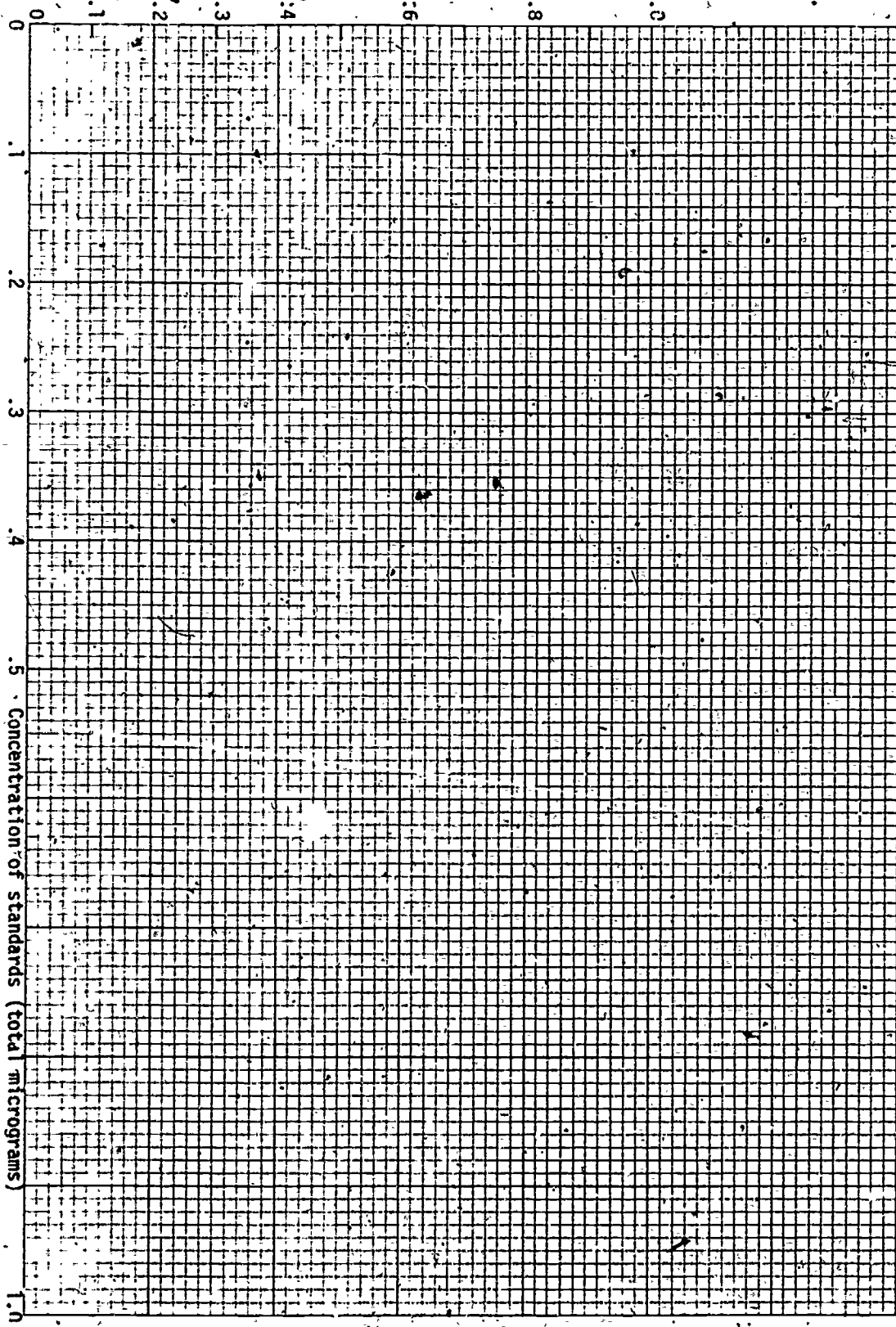


* T Readings


205

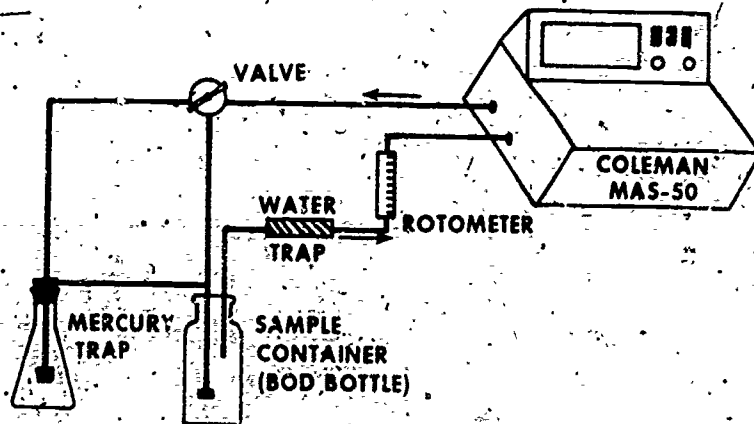

SYSTEM ONE: LIQUID MERCURY TRAP CLOSED SYSTEM

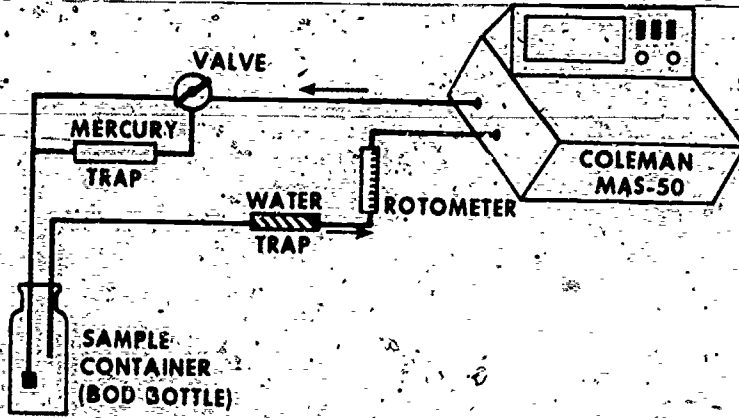

SYSTEM TWO: SOLID MERCURY TRAP CLOSED SYSTEM

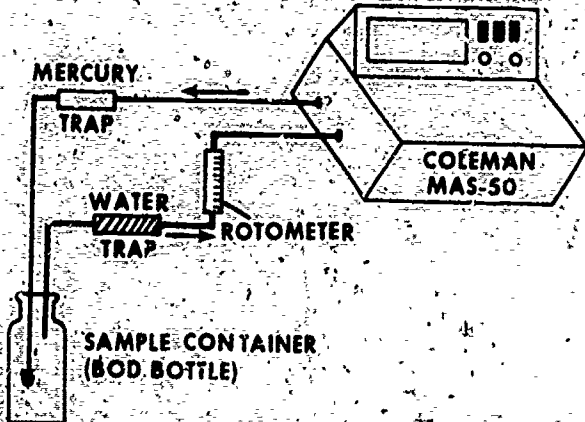

SYSTEM THREE: SOLID MERCURY TRAP OPEN SYSTEM

Figure 1. FLOW SYSTEMS FOR THE COLD VAPOR TECHNIQUE FOR MERCURY

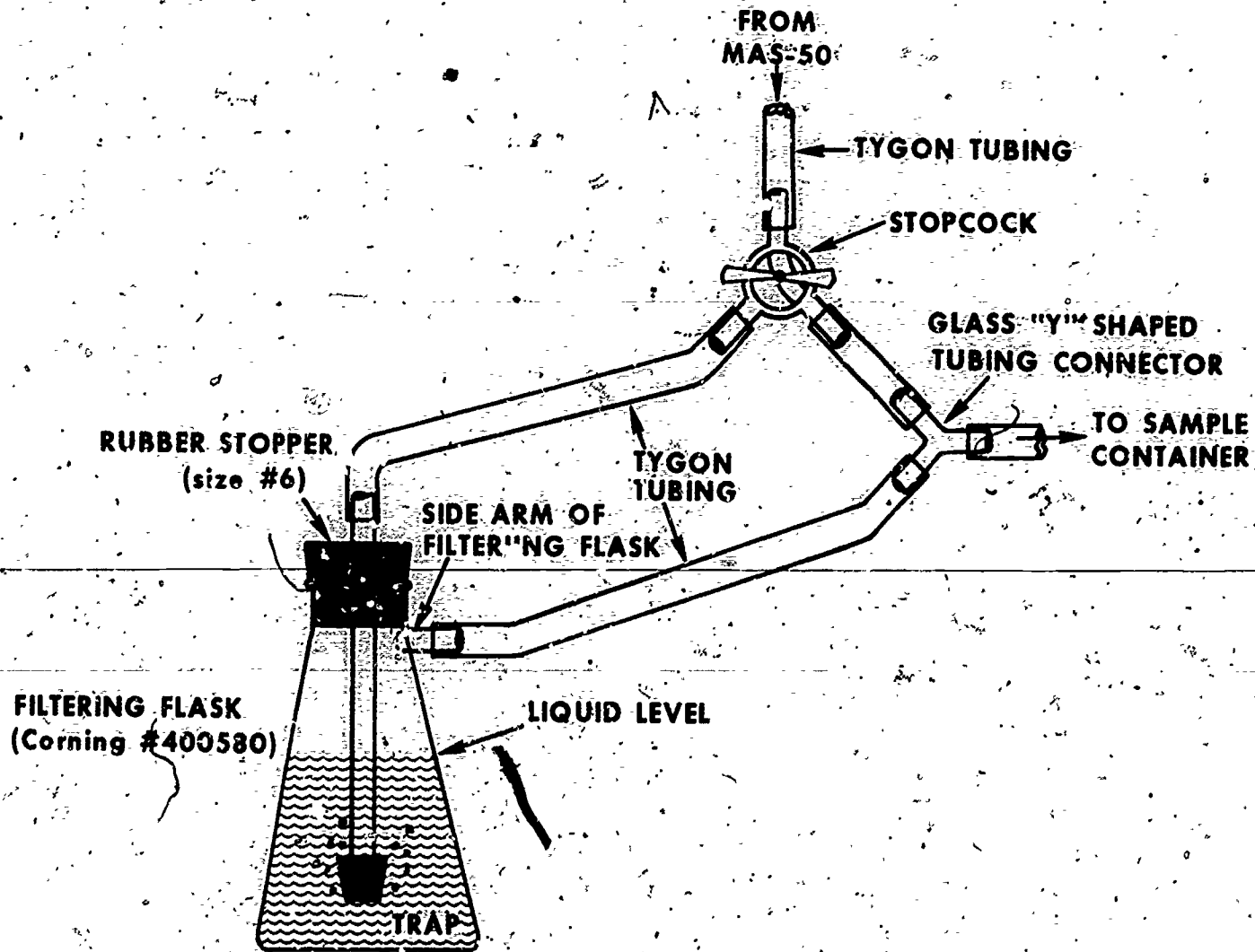


Figure 2. ARRANGEMENT OF TWO-POSITION STOPCOCK AND MERCURY TRAP

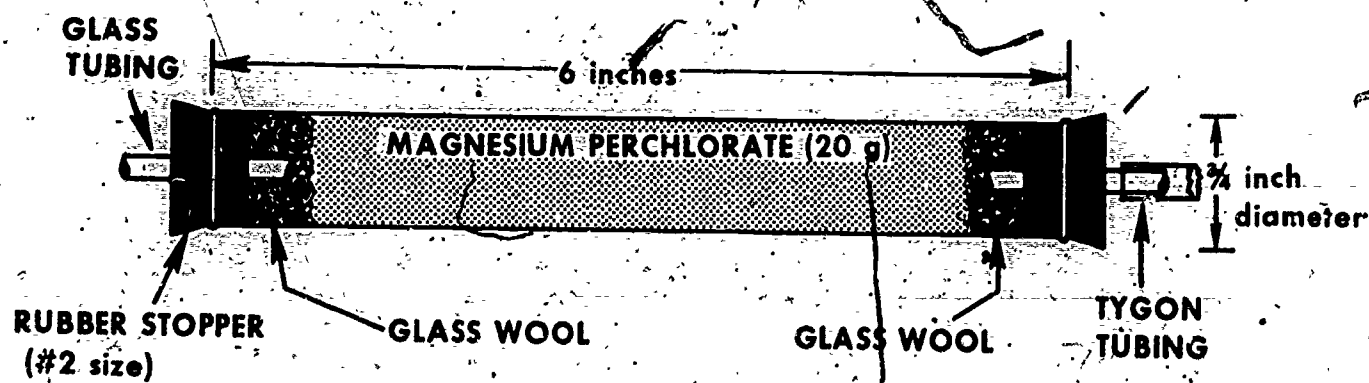


Figure 3. DRYING TUBE

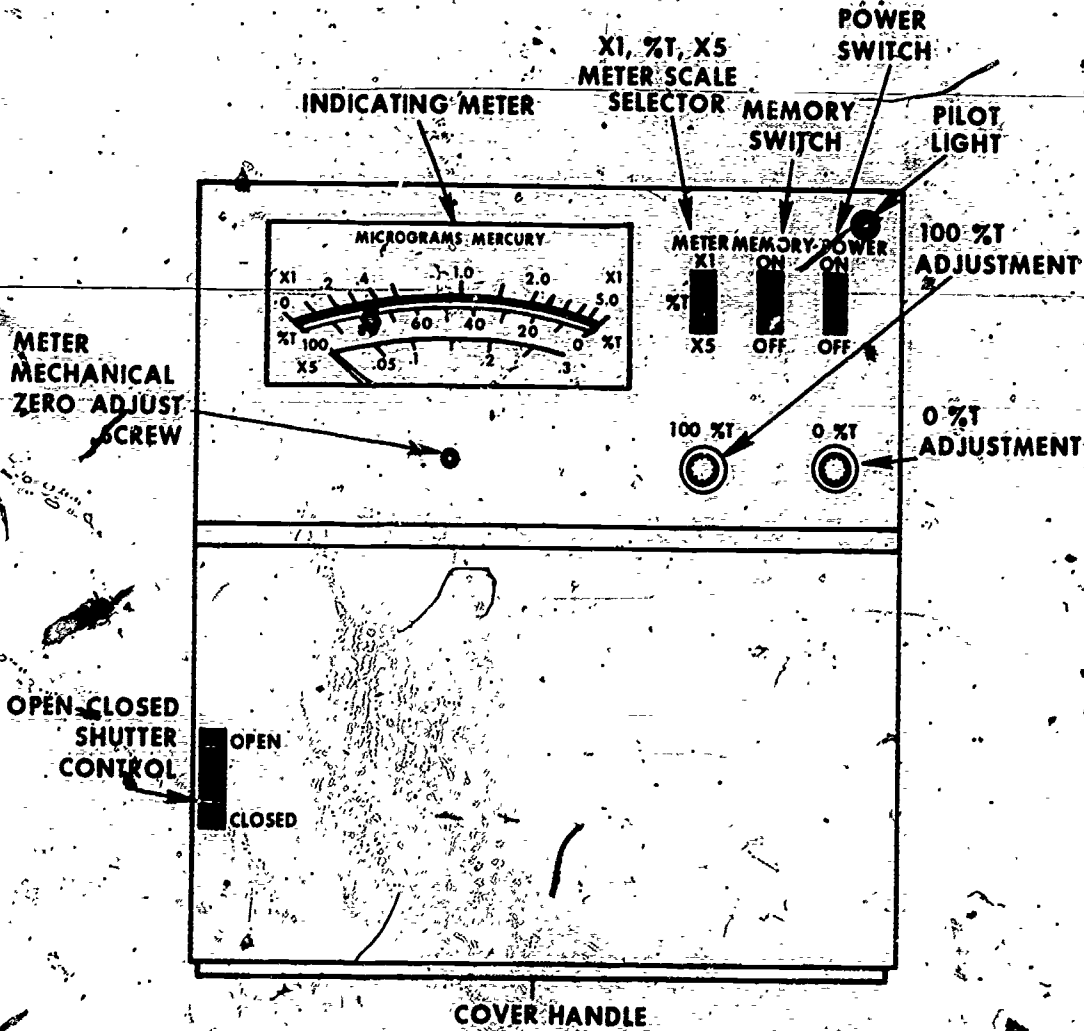


Figure 4. FRONT CONTROLS AND OPERATIONAL COMPONENTS

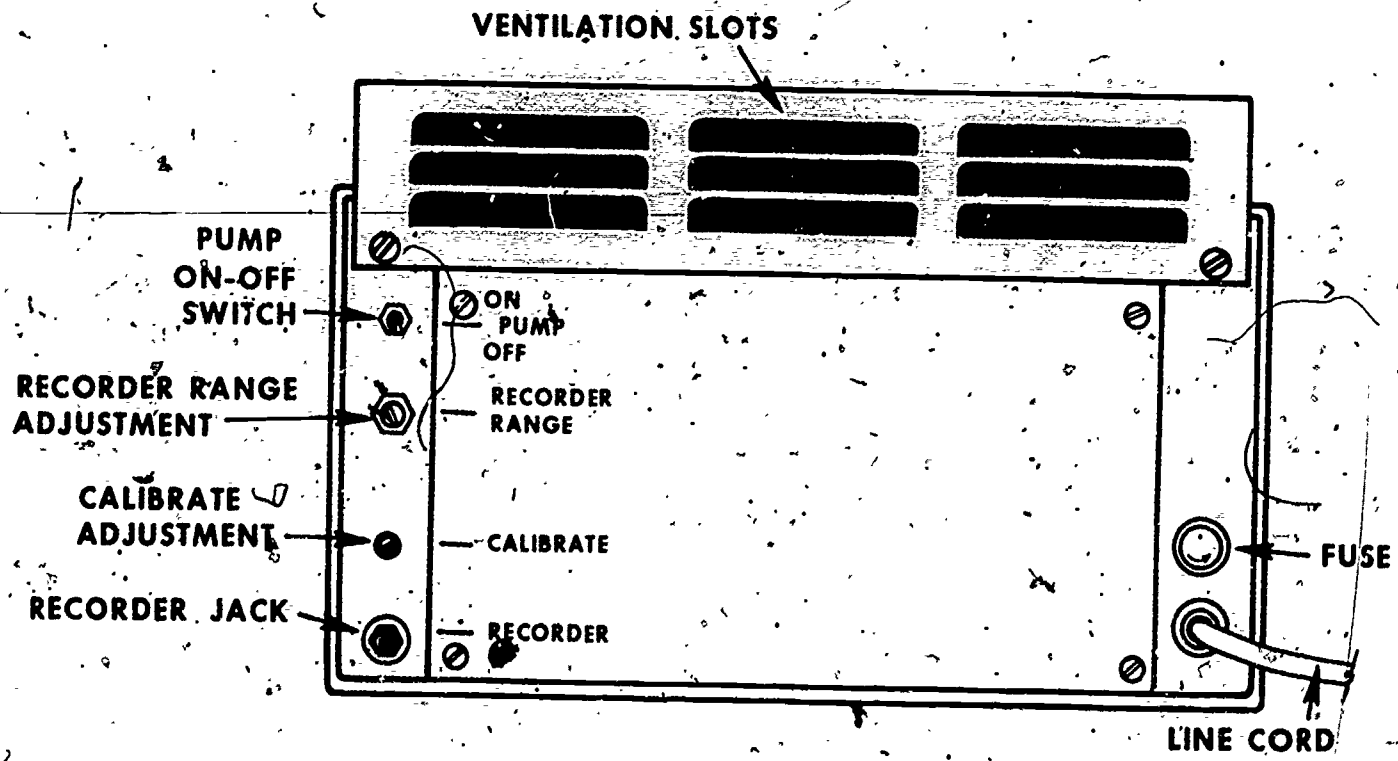


Figure 5. REAR PANEL

A PROTOTYPE FOR DEVELOPMENT OF
ROUTINE OPERATIONAL PROCEDURES

for the

DETERMINATION OF POTASSIUM USING FLAME PHOTOMETRY

as applied in

WASTEWATER TREATMENT FACILITIES
and in the
MONITORING OF EFFLUENT WASTEWATERS

Developed by the

National Training and Operational Technology Center
Municipal Operations and Training Division
Office of Water Program Operations
U.S. Environmental Protection Agency

214

EFFLUENT MONITORING PROCEDURE: Determination of Potassium Using Flame Photometry

This operational procedure was developed by:

NAME Charles R. Feldmann

ADDRESS EPA, OHPO, NTOTC, Cincinnati, Ohio 45268

POSITION Chemist-Instructor

EDUCATION AND TECHNICAL BACKGROUND

B.S. - Chemistry

M.S. - Chemistry

1-1/2 years Industrial Chemist

4 years additional Graduate School

4 years College Chemistry Instructor

1-1/2 years DHEW - Air Pollution Program, Chemist

10 years DI - EPA, Chemist-Instructor

EFFLUENT MONITORING PROCEDURE: Determination of Potassium Using Flame Photometry

1. Analysis Objectives:

The learner will determine the potassium content of a sewage effluent.

2. Brief Description of Analysis:

The learner will prepare a calibration graph using the percent transmittance and the concentration of a series of standard potassium solutions. The percent transmittance of a sample of sewage effluent will next be determined, and the potassium concentration found by using the calibration graph.

The work will be done using a Beckman Model B spectrophotometer with flame photometry attachments. Although this procedure has been prepared for use with a specific instrument, other equivalent commercially available instruments may of course be used. Mention of a particular brand name does not constitute endorsement by the U.S. Environmental Protection Agency.

3. Applicability of this Procedure:

a. Range of Concentration:

Concentration levels as low as 0.1 mg/liter can be detected in a good flame photometer. Although the cited reference (below) does not specifically state the highest potassium concentration which may be determined, a limit of 10 mg/liter may be inferred. Therefore, sample dilution may be required.

b. Pretreatment of Samples:

The digestion procedure referenced in the Federal Register, Wednesday, December 1, 1976, Part II, table I, footnote 15 (Methods for Chemical Analysis of Water & Wastes, 1974, page 83, par. 4.1.4.) is given in this effluent monitoring procedure.

c. Treatment of Interferences in the Sample:

When the following ratios are exceeded, an interference in the determination of potassium is present, and the cited reference must be consulted.

<u>Ratio</u>	<u>Value of the Ratio</u>
sodium/potassium	5/1
calcium/potassium	10/1
magnesium/potassium	100/1

This procedure was taken from the Standard Methods, 14th ed., pg. 234, Method 317 A, and from the Beckman Instrument Manual.

EFFLUENT MONITORING PROCEDURE: Determination of Potassium Using Flame Photometry

Equipment and Supply Requirements

A. Capital Equipment:

1. Beckman Model B. spectrophotometer with flame photometer attachments and instrument manual
2. Source of distilled water
3. Source of deionized water
4. Oven, for use at 110°C and at 103°C (the oven used in the sodium procedure CH.MET.es.EMP.1b.7.77, Determination of Sodium Using Flame Photometry, may be used here)
5. Analytical balance, 100 g capacity

B. Reusable Supplies:

1. Trip balance, 100 g capacity
2. Pencil or pen
3. Twelve inch ruler
4. Graduated cylinder, 50 ml (for glassware cleaning)
5. Beaker, 150 ml
6. Beaker, 250 ml (for glassware cleaning)
7. Beaker, 2 liter (for glassware cleaning)
8. Laboratory apron
9. Safety glasses
10. Glass stoppered bottle, 1 liter (for storage of cleaning solution)
11. Three polyethylene bottles, 200 ml
12. One or two polyethylene bottles, 1 liter. (See page 6-13, steps 9 and 10.)
13. Graduated cylinder, 250 ml
14. Erlenmeyer flask, 125 ml
15. Erlenmeyer flask, 500 ml
16. Graduated cylinder, 100 ml
17. Graduated cylinder, 10 ml
18. Plastic weighing boat, about 2 inches square
19. Small spatula
20. One or two volumetric flasks, 1 liter (See page 6-13, 9a.)
21. Pipet, 100 ml
22. Two plastic squeeze bottles
23. Funnel, about 60 mm diameter
24. One piece Whatman number 40 (or equivalent) filter paper (to fit the funnel)
25. One ring (to support the funnel)
26. Six, or twelve, volumetric flasks, 100 ml (See page 6-17, 3a.)
27. One pipet bulb
28. One or two graduated pipets, 10 ml (See page 6-17, 3a.)
29. Eight, or sixteen, small beakers (depending on whether one or both sets of standards are prepared); either glass (supplied with the flame photometer) or disposable plastic, and used to hold solution for aspiration.
30. Two pressure regulators (one for hydrogen, one for oxygen)
31. Wrenches for use in attaching pressure regulators to gas cylinders
32. Wrench for use in making hose connections
33. Two clamps (for safely anchoring gas cylinders)
34. Screwdriver, medium size
35. One ring stand
36. Hot plate

EFFLUENT MONITORING PROCEDURE: Determination of Potassium Using Flame Photometry

B. Reusable Supplies (Continued)

37. Thermometer, 100°C
38. One piece of rubber or plastic tubing to fit onto the tube extending from the bottom of the atomizer burner, 1 ft.
39. Asbestos gloves or crucible tongs
40. One piece of pH sensitive paper (to measure a pH of 2)

C. Consumable Supplies:

1. Potassium chloride, KCl, 3 g.
2. One piece graph paper; divided into squares of equal size
3. Sodium dichromate, Na₂Cr₂O₇ (for cleaning glassware)
4. Concentrated sulfuric acid, H₂SO₄ (for cleaning glassware)
5. Soap (for cleaning glassware)
6. Brush (for cleaning glassware)
7. Brush (for cleaning balance)
8. Paper towels
9. Matches
10. Concentrated nitric acid, HNO₃
11. Cylinder of hydrogen gas (for use with the flame photometer)
12. Cylinder of oxygen gas (for use with the flame photometer)
13. Concentrated hydrochloric acid, HCl
14. Distilled concentrated nitric acid, HNO₃*

All reagents should be of high quality. Different chemical manufacturers may have different ways of indicating a high quality reagent. While no endorsement of one chemical manufacturer over another is intended, the following are some designations used in four chemical catalogs to indicate high quality reagents.

<u>Catalog</u>	<u>Designations</u>
Thomas	Reagent, ACS, Chemically Pure (CP)
Matheson; Coleman & Bell	Reagent, ACS
Curtin Matheson Scientific, Inc.	Primary Standard, ACS, AR
Fisher	Certified, ACS

* This reagent is used for acidifying the sample at the time of collection. It must be of higher purity than Reagent, ACS, etc. J. T. Baker "Ultrex" nitric acid is an example of a higher purity acid.

Beckman Model B spectrophotometer
with flame photometry attachments

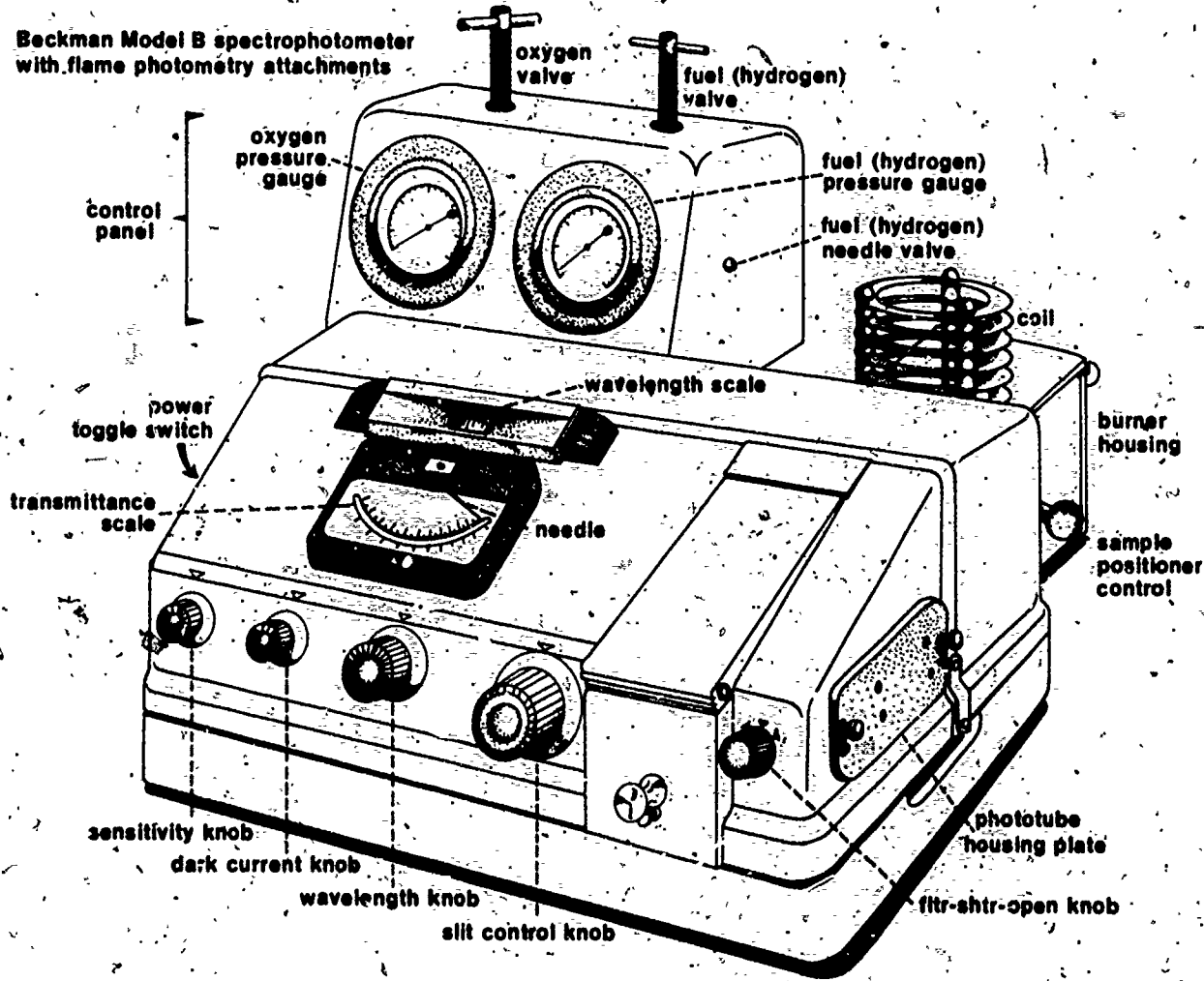


Figure 1

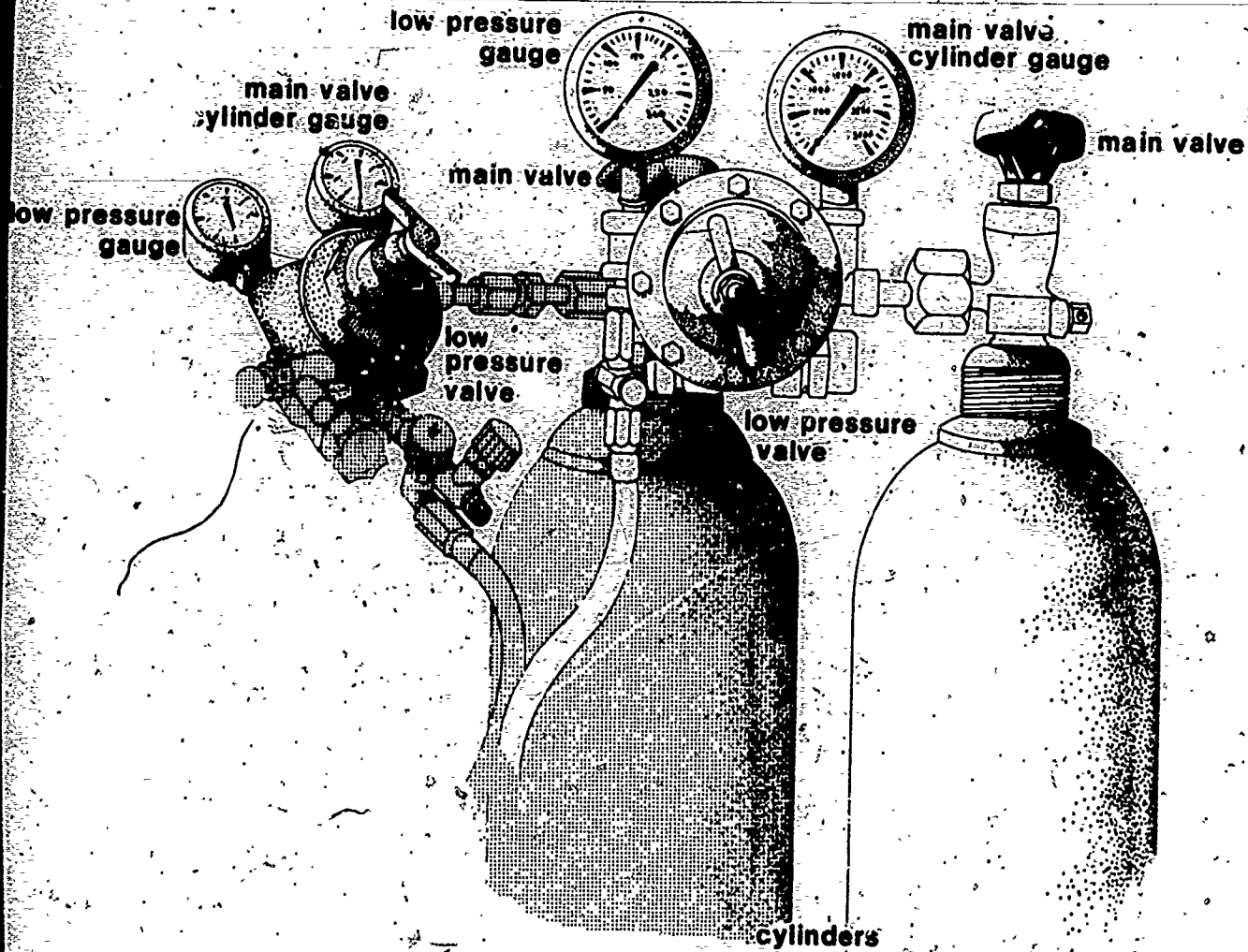


Figure 2

220

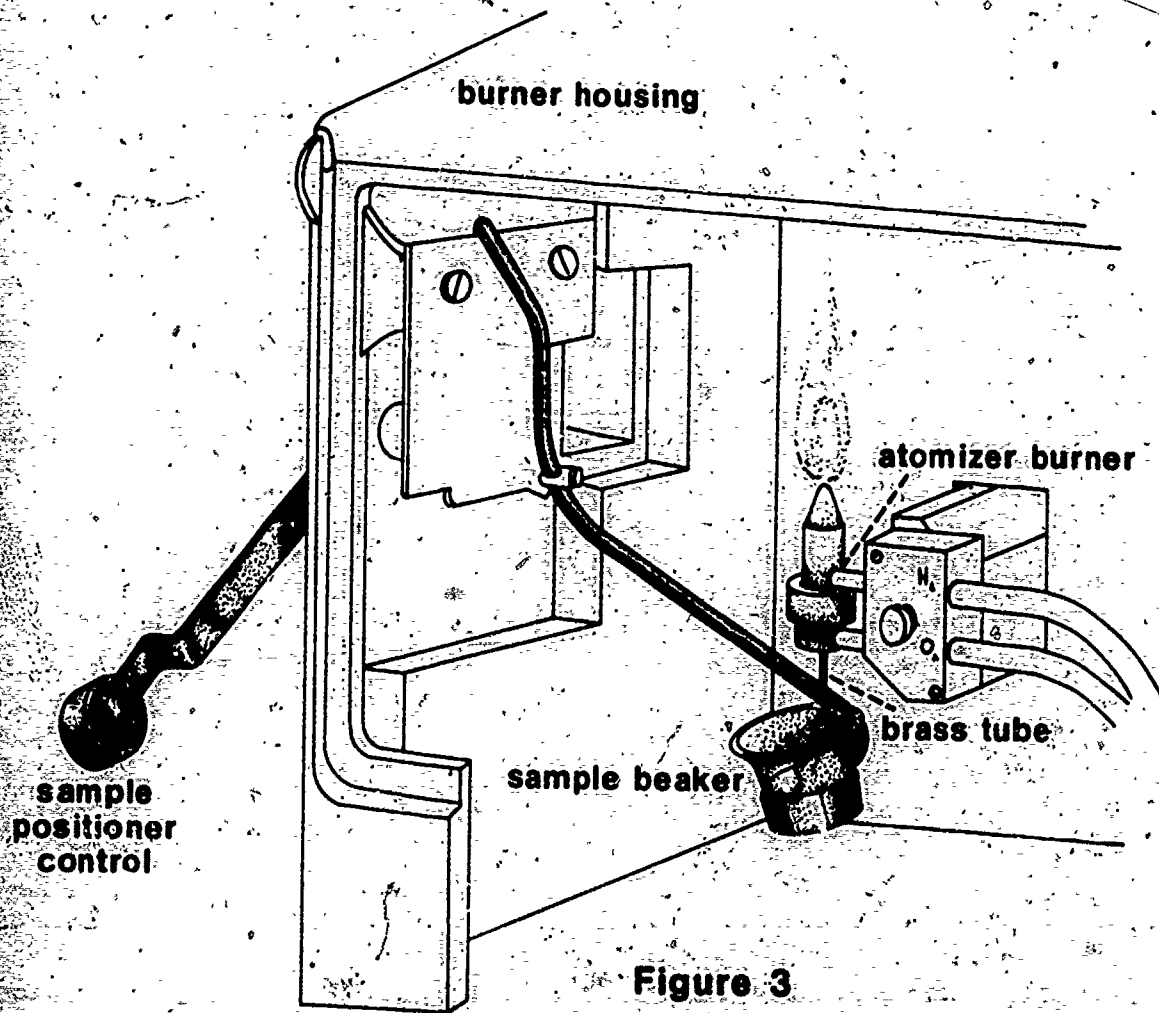


Figure 3

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
A. Equipment Preparation			
1. Cleaning of Glassware	<ol style="list-style-type: none"> 1. Clean all glassware and polyethylene bottles. 2. Rinse with nitric acid solution. 3. Rinse with tap water. 4. Rinse with deionized water. 	<ol style="list-style-type: none"> 2a. Preparation of the acid is described in B.1. 3a. Ten times. 4a. Five times. 	V.A.1.1
2. Balance Preparation	<ol style="list-style-type: none"> 1. Check the analytical balance for cleanliness and proper operation. 	<ol style="list-style-type: none"> 1a. Consult the manufacturer's manual if the balance does not operate properly. 	
3. Spectrophotometer	<ol style="list-style-type: none"> 1. Attach the oxygen and hydrogen cylinders to the Beckman Model B spectrophotometer with flame photometry attachments. 2. Turn the wavelength knob to a reading of 760 nanometers. 3. Turn the slit control knob to a reading of 0.4 mm. 4. Turn the sensitivity knob to the standby position. 5. Turn the fltr-shtr-open knob to the shtr position. 	<ol style="list-style-type: none"> 1a. Consult the manufacturer's manual for specific instructions. 1b. For the remainder of this procedure, the words instrument, or flame photometer, will be used to mean this particular instrument. 5a. This is the closed position. 	

222

223

EFFLUENT MONITORING PROCEDURE: Determination of Potassium Using Flame Photometry

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>A. Equipment Preparation (Continued)</p> <p>4. Phototube</p>	<p>1. Remove the phototube housing plate (see Figure 1).</p> <p>2. Check the condition of the desiccant in the phototube compartment.</p>	<p>1a. The red sensitive phototube should be used for potassium determination.</p> <p>1b. Refer to figures 1, 2, & 3 when other parts of the instrument are mentioned in this procedure.</p> <p>2a. It may need drying at 103-105°C for 1 hour, followed by cooling in a desiccator for 30 min.</p> <p>2b. The spectrophotometer will be turned on in E.</p>	
<p>B. Reagent Preparation</p> <p>1. Nitric Acid, HNO₃ Solution, 1 + 15</p>	<p>1. Measure 150 ml of distilled water.</p> <p>2. Pour it into a 500 ml Erlenmeyer flask.</p> <p>3. Measure 10 ml of concentrated nitric acid, HNO₃.</p> <p>4. Pour the nitric acid slowly into the distilled water.</p> <p>5. Swirl the flask.</p> <p>6. Store the solution in a polyethylene bottle.</p>	<p>1a. Use a 250 ml graduated cylinder.</p> <p>3a. Use a 10 ml graduated cylinder.</p> <p>5a. To mix the acid and water.</p> <p>6a. This solution is used for rinsing glassware and polyethylene bottles.</p> <p>6b. Larger amounts may be prepared using the same proportion of acid and water.</p>	

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
B. Reagent Preparation (continued)			
2. Nitric Acid, HNO ₃ , Distilled	1. See 1a.	1a. This reagent is used undiluted for acidifying the sample at the time of collection.	
3. Hydrochloric Acid, HCl, Distilled, 1 + 1	1. Measure 2.5 ml of distilled hydrochloric acid, HCl.	1a. Use a 10 ml graduated cylinder. 1b. Five ml of the 1 + 1 mixture are needed for each sample. Therefore, larger volumes may be prepared as necessary.	
	2. Pour it into a 125 ml Erlenmeyer flask.	2a. Use a larger flask if larger amounts of the mixture are being prepared.	
	3. Measure 2.5 ml of deionized water.	3a. Use the 10 ml graduated cylinder from 1a.	
	4. Pour it into the same flask.		
	5. Swirl the flask.	5a. To mix the acid and water.	
	6. Store the solution in a polyethylene bottle.		
4. Stock Potassium Solution	1. Weigh 3 g. of potassium chloride, KCl.	1a. Use a trip balance. 1b. This is slightly more than is needed.	
	2. Dry it at 110°C.	2a. For 1 hour.	
	3. Cool it in a desiccator.	3a. For 30 minutes.	
	4. Weigh 1.907 g. of the potassium chloride, KCl.	4a. Use an analytical balance.	
	5. Transfer it to a 1 liter volumetric flask.	5a. Use deionized water in the plastic squeeze bottle.	

226

227

EFFLUENT MONITORING PROCEDURE: Determination of Potassium Using Flame Photometry

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>B. Reagent Preparation (continued)</p>	<p>6. Dissolve the salt in de-ionized water in the flask.</p> <p>7. Dilute to the 1 liter mark.</p> <p>8. Mix thoroughly.</p> <p>9. If the expected potassium concentration in the sample is between 1.0 and 10.0 mg/liter, prepare the intermediate potassium solution (B.5 below).</p> <p>10. If the expected potassium concentration is between 0.1 and 1.0 mg/liter, prepare the standard potassium solution (B.6 below).</p>	<p>7a. With deionized water.</p> <p>8a. The concentration of this solution is 1.0 mg of potassium per 1.0 ml of solution.</p> <p>9a. If there is doubt about the expected potassium concentration, both the intermediate and standard potassium solutions should be prepared, along with two calibration graphs.</p>	
<p>5. Intermediate Potassium Solution</p>	<p>1. Pipet 100.0 ml of the stock potassium solution (B.4.) into a 1 liter volumetric flask.</p> <p>2. Dilute to the 1 liter mark.</p> <p>3. Mix thoroughly.</p>	<p>2a. With deionized water.</p> <p>3a. The concentration of this solution is 0.1 mg of potassium per 1.0 ml of solution.</p>	

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>B. Reagent Preparation (continued)</p> <p>6. Standard Potassium Solution</p>	<ol style="list-style-type: none"> 1. Pipet 100.0 ml of the intermediate potassium solution (B.5.) into a 1 liter volumetric flask. 2. Dilute to the 1 liter mark. 3. Mix thoroughly. 4. Transfer the potassium solutions you have prepared to polyethylene bottles. 	<ol style="list-style-type: none"> 2a. With deionized water. 3a. The concentration of this solution is 0.01 mg of potassium per 1.0 ml of solution. 	
<p>C. Sample Collection</p>	<ol style="list-style-type: none"> 1. Collect the sample in a polyethylene bottle. 2. Acidify the sample with distilled nitric acid, HNO_3. 3. Gently swirl the bottle. 4. Check the pH of the sample. 	<ol style="list-style-type: none"> 1a. Although only a few ml of sample are actually needed, a larger volume, such as 200 ml, is more convenient to work with. 2a. Use 5 ml of acid/liter of sample; 1.0 ml for 200 ml of sample. 2b. Measure the acid in a 10 ml graduated cylinder. 3a. To mix the acid and sample. 4a. Use pH sensitive paper. 4b. It must be less than 2. If it is not, add a few more drops of acid, swirl the bottle, and recheck the pH. Repeat the procedure until the pH is less than 2. 4c. Do not store the sample for more than 6 months before completing the analysis. 	

230

231

EFFLUENT MONITORING PROCEDURE: Determination of Potassium Using Flame Photometry

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>D. Sample Digestion</p>	<ol style="list-style-type: none"> 1. Shake the sample container. 2. Measure 100 ml of the acidified sample. 3. Pour it into a 250 ml Erlenmeyer flask. 4. Measure 5 ml of 1 + 1 distilled hydrochloric acid, HCl. 5. Pour it into the flask. 6. Swirl the flask. 7. Put the flask on a hot-plate. 8. Turn on the hotplate and adjust the temperature setting so the sample comes to a temperature of 95°C. 9. After the sample has come to 95°C, continue the heating for 15 min. 10. Assemble a filtration apparatus. 	<ol style="list-style-type: none"> 1a. To ensure that it is homogeneous. 2a. Use a 100 ml graduated cylinder. 3a. If any solids remain in the cylinder, rinse them into the flask with a few ml of deionized water. 4a. Use a 10 ml graduated cylinder. 6a. To mix the acid and sample. 8a. Check the temperature of the sample with a thermometer. 8b. Do not allow the temperature of the sample to go above 95°C. 9a. During the 15 minute period, proceed with Step 10. 10a. A ring stand, a funnel (approximately 60 mm diameter), a ring to hold the funnel, and a piece of medium porosity filter paper (such as Whatman number 40). 10b. Use a clean 100 ml graduated cylinder as receiver. 	

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>D. Sample Digestion (continued)</p>	<p>11. After the 15 min. heating period, remove the flask from the hotplate.</p> <p>12. Allow it to cool at room temperature for about 15 min.</p> <p>13. Filter the sample through the filter paper into the cylinder.</p> <p>14. Wash the filter paper with small portions (a few ml) of deionized water, and catch the washings in the cylinder.</p> <p>15. Cover the cylinder with an inverted 150 ml beaker.</p>	<p>11a. Be cautious of the hot flask. Use a glove or tongs.</p> <p>13a. Solids in the sample would clog the burner on the flame photometer.</p> <p>14a. When the volume in the cylinder is 100 ml, remove the cylinder from under the funnel.</p> <p>15a. Until the analysis. 15b. To prevent contamination.</p>	
<p>E. Spectrophotometer Warm-up</p>	<p>1. Plug the power cord on the instrument into a wall outlet.</p> <p>2. Turn on the power toggle switch.</p> <p>3. Allow the instrument to warm up for 45 minutes. While it is warming up, proceed with Section F.</p>	<p>1a. 115 volts, 50/60 cycle.</p> <p>2a. The indicator light will go on. 2b. Do not turn on the lamp toggle switch. It is not used when making flame photometry measurements.</p> <p>3a. With older instruments, 2 hours or longer would be a better warm-up time.</p>	

EFFLUENT MONITORING PROCEDURE: Determination of Potassium Using Flame Photometry

OPERATING PROCEDURES	STEP SEQUENCE	- INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>F. Preparation of Standards</p>	<ol style="list-style-type: none"> 1. While the instrument is warming up, prepare the potassium standards. 2. Assemble six 100 ml volumetric flasks. 3. Mark them 0, 2, 4, 6, 8, and 10 if you use the intermediate potassium solution (B.5., 0.1 mg/ml), or 0.0, 0.2, 0.4, 0.6, 0.8, and 1.0 if you use the standard potassium solution (B.6., 0.01 mg/ml). 4. Pipet 2.0 ml of the intermediate potassium solution into the flask marked 2, and 2.0 ml of the standard potassium solution into the flask marked 0.2. 5. Pipet 4.0 ml of the intermediate potassium solution into the flask marked 4, and 4.0 ml of the standard potassium solution into the flask marked 0.4. 6. Prepare the 6, 8, and 10, and 0.6, 0.8, and 1.0 flasks in a similar manner. 	<ol style="list-style-type: none"> 1a. Two standard potassium solutions were prepared. One concentration was 0.01 mg/ml (the standard potassium solution (B.6.)), and the other was 0.1 mg/ml (the intermediate potassium solution (B.5.)). Experience, or trial and error, will determine which potassium solution you use. 3a. If you <u>do not know</u> which potassium solution to use, then <u>prepare</u> two sets of standards with six more 100 ml volumetric flasks and a second 10 ml graduated pipet. If you <u>do know</u> the expected potassium concentration in the sample, and therefore, which potassium solution to use, only one set of standards is needed. 4a. Use one 10 ml graduated pipet for the intermediate solution, and a second 10 ml graduated pipet for the standard solution. 5a. Use the appropriate pipet. 6a. Use the appropriate pipet. 	

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>F. Preparation of Standards (continued)</p>	<p>7. Add deionized water to the mark in each of the six or twelve flasks.</p> <p>8. Mix thoroughly.</p>	<p>8a. The concentrations are: 0.0, 2.0, 4.0, 6.0, 8.0, and 10.0 mg/liter in the one set of standards, and 0.0, 0.2, 0.4, 0.6, 0.8, and 1.0 mg/liter in the other.</p>	
<p>G. Analysis</p> <p>1. Standards and Sample</p> <p>2. Lighting the Flame</p>	<p>1. Mark the 5 ml beakers (six or twelve) in the same way that the 100 ml volumetric flasks are marked.</p> <p>2. Pour the standards into the appropriate 5 ml beakers.</p> <p>3. Shake the sample thoroughly.</p> <p>4. Fill another 5 ml beaker with sample.</p> <p>5. Fill another 5 ml beaker with deionized water.</p> <p>1. Raise the door on the right side of the burner housing.</p>	<p>1a. Either glass or plastic disposable. 1b. Used for aspirating solutions.</p> <p>2a. Fill the beakers about three-fourths full. 2b. Caution: Throughout the remainder of this procedure, be careful not to get your fingers into any of the solutions. This could cause an error in the result.</p> <p>3a. Even if it was filtered earlier.</p> <p>4a. About three-fourths full.</p> <p>5a. About three-fourths full.</p>	

EFFLUENT MONITORING PROCEDURE: Determination of Potassium Using Flame Photometry

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>G. Analysis (continued)</p>	<p>2. Raise the sample positioner control between the instrument and the burner housing to the vertical position as far as it will go.</p> <p>3. Close the oxygen and fuel (hydrogen) valves on top of the control panel.</p> <p>4. Close the low pressure valve on the regulators attached to the hydrogen and oxygen cylinders.</p> <p>5. Close the fuel (hydrogen) needle valve on the right side of the control panel.</p> <p>6. Open the fuel (hydrogen) needle valve on the right side of the control panel.</p> <p>7. Open the main valve on the oxygen cylinder.</p> <p>8. Open the low pressure valve on the oxygen cylinder.</p> <p>9. Set the pressure for 30 pounds per square inch.</p>	<p>2a. Do not force the sample positioner control.</p> <p>3a. By turning them counter clockwise until they just begin to feel loose; i.e., until they can be wiggled very slightly left and right.</p> <p>4a. By turning the handles counter-clockwise until they just begin to feel loose; i.e., until they can be wiggled very slightly left and right.</p> <p>5a. By turning it clockwise with a screwdriver as far as it will go.</p> <p>5b. Do not force the screwdriver.</p> <p>6a. By turning it counter clockwise with a screwdriver.</p> <p>6b. About 1/8 turn.</p> <p>7a. By turning the handle counter clockwise two full turns.</p> <p>7b. For safety, <u>oxygen</u> is always turned on <u>first</u>, and turned off <u>last</u>.</p> <p>8a. By turning the handle clockwise until the needle moves away from zero.</p> <p>9a. By turning the valve clockwise.</p>	

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>G. Analysis (continued)</p>	<p>10. Open the oxygen valve on the control panel.</p>	<p>10a. By turning the handle clockwise until the needle moves away from zero.</p>	
	<p>11. Set the pressure for 13 pounds per square inch.</p>	<p>11a. By turning the handle clockwise. 11b. The actual oxygen pressure should be written on a tag attached to the atomizer burner. 11c. The optimum pressure may be checked if no tag is attached to the burner. 11d. Attach a small diameter piece of rubber or plastic tubing about 1 foot long to the brass tube extending from the bottom of the atomizer burner. The rubber or plastic tubing should fit the brass tube snugly. Place the other end of the rubber or plastic tubing in a 10 ml graduated cylinder. Fill it to the .0 line with deionized water. When the oxygen is flowing at the proper rate, the deionized water should be aspirated at the rate of 1.5 - 2.0 ml per minute. This rate corresponds to about thirteen pounds pressure per square inch.</p>	
	<p>12. Remove the tubing and cylinder.</p>		
	<p>13. Open the low pressure valve on the hydrogen cylinder.</p>	<p>13a. By turning the handle clockwise until the needle moves away from zero.</p>	
	<p>14. Set the pressure for 10 pounds per square inch.</p>	<p>14a. By turning the valve clockwise.</p>	
	<p>15. Open the fuel (hydrogen) valve on the control panel.</p>	<p>15a. By turning the handle clockwise.</p>	
<p>16. Set the pressure for 4.5 pounds per square inch.</p>			

EFFLUENT MONITORING PROCEDURE: Determination of Potassium Using Flame Photometry

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
G. Analysis (continued)	<p>17. Cautiously bring a lighted match to the tip of the atomizer burner.</p> <p>18. Close the door on the right side of the burner housing.</p> <p>19. Place the beaker containing the highest concentration standard (10 mg/liter or 1.0 mg/liter) into the holder.</p> <p>20. Swing the beaker into the burner housing compartment using the sample positioner control.</p> <p>21. Turn the sensitivity knob to position 1.</p> <p>22. Make the needle read 0% transmittance using the dark current knob.</p>	<p>17a. The oxygen-hydrogen mixture will ignite with a loud pop and hissing sound.</p> <p>18a. Caution: hot gases are escaping from the top of the coil. Do not place your hand or any other part of your body over the coil while the gases are burning.</p> <p>19a. See figure 3.</p> <p>19b. If two sets of standards are being used, take readings on the 10, 8, 6, 4, 2, and 0 mg/liter solutions <u>first</u>, and <u>then</u> the 1.0, 0.8, 0.6, 0.4, 0.2, and 0.0 mg/liter solutions. Do not <u>mix</u> the two sets of standards.</p> <p>20a. A distinct change in sound will be heard as the solution is aspirated into the burner.</p> <p>20b. For the remainder of this procedure, if the sound caused by the aspiration of a solution changes while the <u>solution is being aspirated</u>, it probably means that all of the liquid in the beaker has been aspirated, and the beaker is empty.</p> <p>20c. Swing the beaker out of the burner housing, refill it with the appropriate solution, and swing it back into the burner housing.</p> <p>20d. While the solution is aspirating, check that the hydrogen pressure is still 10 pounds per square inch and that the oxygen pressure is still at the optimum valve determined above. If one or both of the pressures have changed, reset them.</p> <p>22a. In this procedure, all readings will be made on the transmittance portion of the scale, not the absorbance portion.</p>	

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
G. Analysis (continued)	23. Turn the sensitivity knob to position 4. 24. Repeat step 22. 25. Adjust the fuel needle valve so the tip of the flame comes to the top of the burner housing compartment, not to the top of the coil, on top of the compartment.	25a. Use a screwdriver. 25b. Caution: extremely hot gases are escaping through the coil. 25c. Adjustment of this needle valve setting is not necessary each time the instrument is used, unless the setting has been changed for some reason.	
3. Aspiration of Standards and Sample	1. Turn the fltr-shtr-open knob to the open position. 2. Turn the wavelength knob very slowly clockwise. 3. Record this wavelength and proceed to step 14.	1a. The shutter is open in this position. 1b. In the sltr position, the shutter is closed. 2a. Turn the knob continuously. Do not turn the knob in steps. 2b. One of three things will happen, 2c., 2d., or 2e. 2c. <u>is the condition you are trying to achieve.</u> 2c. First: the needle will move to the left, and at some % transmittance between 50 and 100, will suddenly swing back to the right. In this case, proceed to step 3. 2d. Second: the needle will swing back to the right as in 2c, but at some % transmittance between 0 and 50. In this case, proceed to step 4. 2e. Thirdly: the needle will swing back to the right, but at some point greater than 100% transmittance; or, it will not swing back to the right, it will stay to the left. In this case, proceed to step 9. 3a. Use the example data sheet(s); see pages 29 & 30. 3b. Do not change this wavelength until the procedure instructs you to change it.	

246

247

EFFLUENT MONITORING PROCEDURE: Determination of Potassium Using Flame Photometry:

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATION ^c	TRAINING GUIDE NOTES
<p>G. Analysis (continued)</p>	<p>4. Close the shutter.</p> <p>5. Increase the reading on the slit knob by 0.05 millimeter.</p> <p>6. Turn the wavelength back to 760 nanometers.</p> <p>7. Repeat steps 1 and 2.</p> <p>8. Record this wavelength and proceed to step 14.</p> <p>9. Close the shutter.</p> <p>10. Decrease the reading on the slit knob by 0.05 millimeter.</p> <p>11. Turn the wavelength back to 760 nanometers.</p> <p>12. Repeat steps 1 and 2.</p>	<p>3c. This wavelength should be 768 nanometers. If it is too far away from this valve (less than 763, or more than 773), the calibration of the instrument should be checked using the instrument manual.</p> <p>5a. It is graduated in tenths of a millimeter between 0.0 and 1.5 millimeters.</p> <p>7a. If condition 2d still exists, repeat steps 4 through 7 until the needle moves as described in 2c.</p> <p>8a. Use the example data sheet(s): see pages 29 & 30. 8b. Do not change this wavelength until the procedure instructs you to change it.</p> <p>10a. It is graduated in tenths of a millimeter between 0.0 and 1.5 millimeters.</p> <p>12a. If condition 2e still exists, repeat steps 9 through 11 until the needle moves as described in 2c.</p>	

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
G. Analysis (continued)	<p>13. Record this wavelength and proceed to step 14.</p> <p>14. Turn the slit knob so as to make the needle read 100% transmittance.</p> <p>15. Close the shutter.</p> <p>16. Record a reading of 100% transmittance for the 10 or 1.0 mg/liter solution.</p> <p>17. Swing the beaker out of the burner housing compartment.</p> <p>18. Replace it with the 8 or 0.8 mg/liter solution.</p> <p>19. Swing it into the burner housing compartment.</p> <p>20. Open the shutter.</p> <p>21. Record the % transmittance reading.</p> <p>22. Close the shutter.</p>	<p>13a. Use the example data sheet(s); see pages 29 & 30.</p> <p>13b. Do not change this wavelength until the procedure instructs you to change it.</p> <p>15a. Do not change this setting until the procedure instructs you to change it.</p> <p>16a. Use the example data sheet(s); see pages 29 & 30.</p> <p>16b. This first set of readings you will get on the one or two sets of standards and sample are called peak readings.</p> <p>21a. Use the example data sheet(s); see pages 29 & 30.</p>	

250

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>G. Analysis (continued)</p>	<p>23. Repeat steps 17 through 22 for the 6, 4, 2, and 0 mg/liter solutions and the sample, or the 0.6, 0.4, 0.2, and 0.0 mg/liter solutions and the sample.</p> <p>24. Swing the beaker of sample out of the burner housing compartment.</p> <p>25. Replace it with the beaker of deionized water.</p> <p>26. Swing it into the burner housing compartment.</p> <p>27. Turn the dark current knob so the needle reads 0% transmittance.</p> <p>28. Replace the beaker of deionized water with the beaker of the 10 or 1.0 mg/liter solution.</p> <p>29. Open the shutter.</p> <p>30. Turn the wavelength knob very slowly clockwise.</p>	<p>23a. Use the example data sheet(s); see pages 29 & 30.</p> <p>28a. It may need refilling.</p> <p>29a. The needle should show 100% transmittance; it may be "off" by plus or minus 1 or 2% because of the dark current adjustment. Do not change the slit setting.</p> <p>30a. Turn the knob continuously. Do not turn the knob, stop, and start again.</p> <p>30b. The needle will move to the right.</p> <p>30c. At some low % transmittance value, perhaps less than 1, or even 0, the needle will move no farther to the right.</p>	

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>G. Analysis (continued)</p>	<p>31. Record this wavelength.</p> <p>32. Record this low % transmittance for the 10 or 1.0 mg/liter solution.</p> <p>33. Close the shutter.</p> <p>34. Swing the beaker out of the burner housing compartment.</p> <p>35. Replace it with the 8 or 0.8 mg/liter solution.</p> <p>36. Swing it into the burner housing.</p> <p>37. Open the shutter.</p> <p>38. Record the % transmittance reading.</p> <p>39. Close the shutter.</p> <p>40. Repeat steps 34 through 39 to obtain readings for the 6, 4, 2, and 0 mg/liter solutions and the sample, or the 0.6, 0.4, 0.2, and 0.0 mg/liter solutions, and the sample.</p>	<p>30d. If there is some doubt as to when the needle stops moving to the right, turn the wavelength knob about one-quarter turn counter clockwise and repeat step 30.</p> <p>31a. Use the example data sheet(s); see pages 29 & 30.</p> <p>31b. It is sometimes called the background wavelength.</p> <p>31c. Do not change this wavelength setting for the remainder of the procedure.</p> <p>32a. Use the example data sheet(s); see pages 29 & 30. This second set of readings you will get on the one or two sets of standards and sample are called background readings.</p> <p>38a. Use the example data sheet(s); see pages 29 & 30.</p>	<p></p>

254

255

EFFLUENT MONITORING PROCEDURE: Determination of Potassium Using Flame Photometry

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERAT.	ADJUSTMENTS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>G. Analysis (continued)</p> <p>256</p>	<ol style="list-style-type: none"> 1. Replace the beaker of sample with the beaker of deionized water. 42. Allow the deionized water to aspirate for 10 seconds. 43. Swing the beaker of deionized water out of the burner housing. 44. Close the main valve on the hydrogen cylinder. 45. When the hydrogen pressure registers L on the control panel hydrogen (fuel) gauge, close the main valve on the oxygen cylinder. 46. Turn the sensitivity knob to the standby position. 47. Turn off the power toggle switch. 48. Empty all of the small beakers used. 49. Rinse the beakers thoroughly with tap water. 50. Rinse the beakers thoroughly with deionized water. 	<ol style="list-style-type: none"> 44a. By turning clockwise all the way. 45a. By turning clockwise all the way. 48a. Don't forget the one still in the instrument. 48b. If the beakers are of the disposable type, discard them and proceed to H. If they are glass, do steps 49, 50, and 51. 		<p>257</p>

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
G. Analysis (continued)	51. Allow the beakers to drain dry, and then proceed to Section H.		
H. Calculations	<p>1. For all of the standards, subtract the % transmittance obtained at the background wavelength from the % transmittance obtained at the peak wavelength.</p> <p>2. On "regular" graph paper prepare a calibration graph using the % transmittance differences along the vertical axis, and the corresponding concentrations (10, 8, 6, 4, 2, 0, or 1.0, 0.8, 0.6, 0.4, 0.2, 0.0 mg/liter) on the horizontal axis.</p> <p>3. Repeat step 1 for the sample.</p> <p>4. Determine the potassium concentration in the sample.</p> <p>5. Record the result in mg/liter.</p>	<p>1a. Example: 10.0 mg/liter solution</p> <p>100 = % transmittance at the peak wavelength</p> <p>1 = % transmittance at the background wavelength</p> <p>99 = % transmittance difference</p> <p>2a. If you have used both sets of standards (the intermediate and standard potassium solutions), prepare two calibration graphs. Pages 31 and 32 are sheets of "regular" graph paper.</p> <p>2b. "Regular" graph paper is marked off in squares of equal size.</p> <p>2c. EMP CH:IN.cg.EMP.1a.1.77 may be referred to for specific instructions in calibration graph preparation and use.</p> <p>2d. In the case of the intermediate potassium solution, the calibration graph may have a slight bend.</p> <p>4a. Using the calibration graph.</p> <p>5a. Use the example data sheet(s); see pages 29 & 30.</p>	

258

259

EFFLUENT MONITORING PROCEDURE: Determination of Potassium Using Flame Photometry

EXAMPLE
DATA SHEET

Concentration of Potassium Standards in mg/liter	% Transmittance at the Peak Wavelength; Peak Readings	% Transmittance at the Background Wavelength; Background Readings	% Transmittance Differences; Values in 2nd Column Minus Values in 3rd Col.
_____*	_____	_____	_____
_____	_____	_____	_____
_____	_____	_____	_____
_____	_____	_____	_____
_____	_____	_____	_____
_____	_____	_____	_____
Sample	_____	_____	_____

Concentration of potassium in the sample in mg/liter _____

Peak Wavelength _____ nm

Background Wavelength _____ nm

* The six values in this column will be 0, 2, 4, 6, 8, and 10, or, 0.0, 0.2, 0.4, 0.6, 0.8, and 1.0. If both sets of standards were used, use a second data sheet exactly like this one; see page 30.


EFFLUENT MONITORING PROCEDURE: Determination of Potassium Using Flame Photometry


EXAMPLE
 DATA SHEET
 (See the paragraph at the bottom of Page 29)

<u>Concentration of Potassium Standards in mg/liter</u>	<u>%Transmittance at the Peak Wavelength; Peak Readings</u>	<u>%Transmittance at the Background Wavelength; Background Readings</u>	<u>%Transmittance Differences; Values in 2nd Column Minus Values in 3rd Col.</u>
_____	_____	_____	_____
_____	_____	_____	_____
_____	_____	_____	_____
_____	_____	_____	_____
_____	_____	_____	_____
_____	_____	_____	_____
_____	_____	_____	_____
_____	_____	_____	_____
Sample	_____	_____	_____

Concentration of potassium in the sample in mg/liter _____

Peak Wavelength _____ nm
 Background Wavelength _____ nm

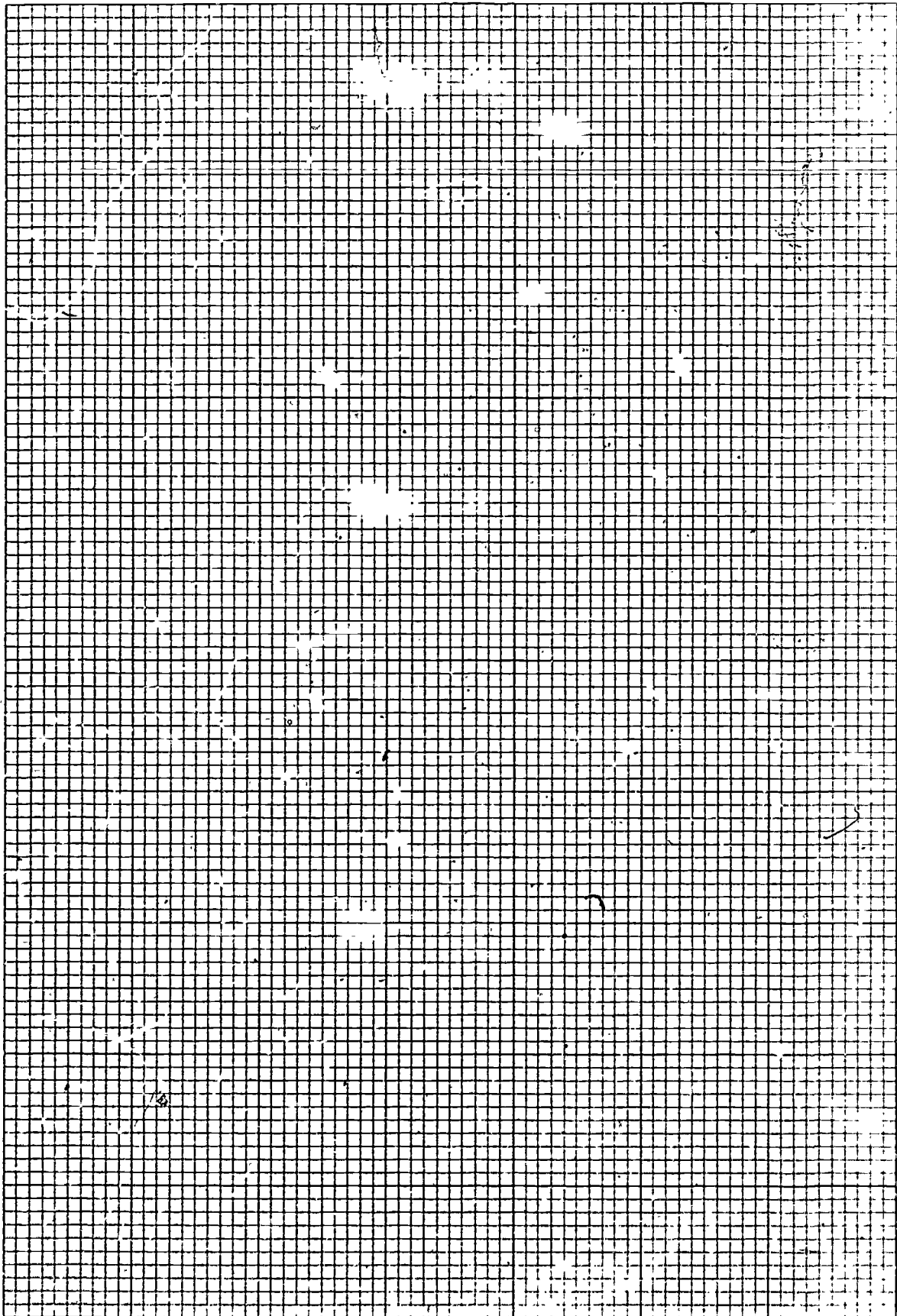

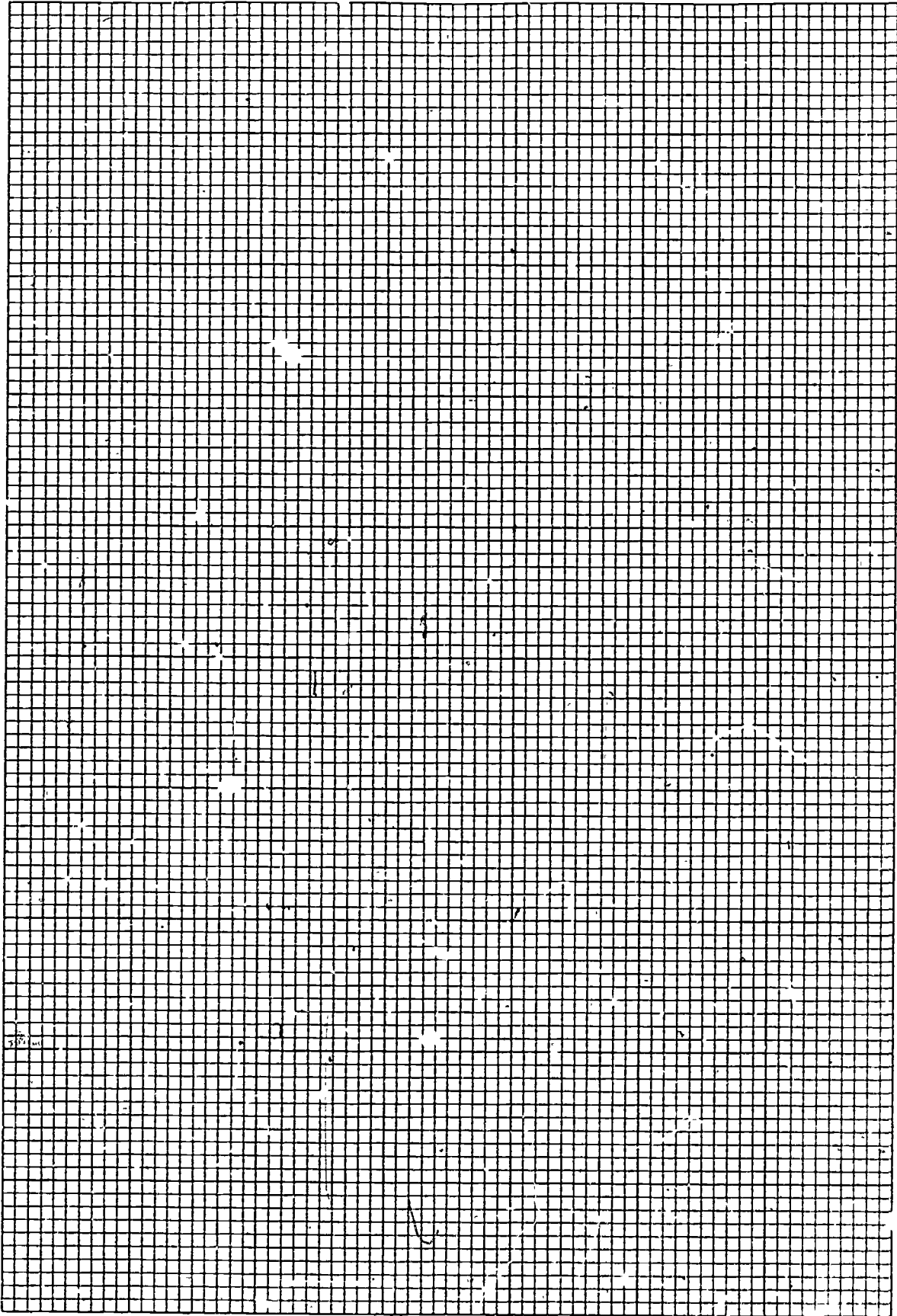

EFFLUENT MONITORING PROCEDURE: Determination of Potassium Using Flame Photometry

TRAINING GUIDE

<u>SECTION</u>	<u>TOPIC</u>
I	Introduction
II	Educational Concepts - Mathematics
III	Educational Concepts - Science
IV	Educational Concepts - Communications
V*	Field and Laboratory Equipment
VI	Field and Laboratory Reagents
VII	Field and Laboratory Analysis
VIII	Safety
IX	Records & Reports

* Training guide materials are presented here under the headings marked *.

EFFLUENT MONITORING PROCEDURE: Determination of Potassium Using Flame Photometry

FIELD AND LABORATORY EQUIPMENT

Section V

	TRAINING GUIDE NOTE	REFERENCES/RESOURCES.
A.1.1	<p>If the glassware is especially dirty and cannot be cleaned with ordinary detergents, chromic acid cleaning may be required.</p> <ol style="list-style-type: none"> 1. Pour 35 ml of distilled water in a 250 ml beaker. 2. Add about 1/8 teaspoon (simply estimate this quantity) of sodium dichromate, $\text{Na}_2\text{Cr}_2\text{O}_7$, to the water. 3. Swirl the beaker until the sodium dichromate has dissolved. 4. Keep repeating steps 2 and 3 until no more sodium dichromate will dissolve. 5. Pour the solution into a 2 liter beaker. 6. Slowly pour 1 liter of concentrated sulfuric acid, H_2SO_4, into the 2 liter beaker. <p>Caution: Use eyeglasses and protective clothing.</p> <ol style="list-style-type: none"> 7. Stir the mixture thoroughly. 8. Store it in a glass stoppered bottle. 9. The cleaning solution should be at a temperature of about 50°C when it is used. 10. It may therefore be necessary to warm the cleaning solution. 11. When using the warm cleaning solution, fill the piece of glassware with the solution. 12. Allow it to soak for 2-3 minutes (or longer). 13. Pour the cleaning solution back into the storage bottle. 14. Rinse the piece of glassware ten times with tap water. 15. The cleaning solution may be reused until it turns green. 16. It should then be discarded. 	<p>Standard Methods, 14th ed., P. 336, section 2c.2)</p>

**A PROTOTYPE FOR DEVELOPMENT OF
ROUTINE OPERATIONAL PROCEDURES**

for the

DETERMINATION OF SODIUM USING FLAME PHOTOMETRY

as applied in

**WASTEWATER TREATMENT FACILITIES
and in the
MONITORING OF EFFLUENT WASTEWATERS**

Developed by the

**National Training and Operational Technology Center
Municipal Operations and Training Division
Office of Water Program Operations
U.S. Environmental Protection Agency**

EFFLUENT MONITORING PROCEDURE: Determination of Sodium Using Flame Photometry

This operational procedure was developed by:

NAME Charles R. Feldmann

ADDRESS EPA, OWPO, NTOTC, Cincinnati, Ohio 45268

POSITION Chemist-Instructor

EDUCATION AND TECHNICAL BACKGROUND

B.S. - Chemistry

M.S. - Chemistry

1-1/2 years Industrial Chemist

4 years additional Graduate School

4 years college Chemistry Instructor

1-1/2 years CHEW - Air Pollution Program, Chemist

10 years DI - EPA, Chemist-Instructor

EFFLUENT MONITORING PROCEDURE: Determination of Sodium Using Flame Photometry

1. Analysis Objective:

The learner will determine the sodium content of a sewage effluent.

2. Brief Description of Analysis:

The learner will prepare a calibration graph using the percent transmittance and the concentration of a series of standard sodium solutions. The percent transmittance of a sample of sewage effluent will next be determined, and the sodium concentration found by using the calibration graph.

The work will be done using a Beckman Model B spectrophotometer with flame photometry attachments. Although this procedure has been prepared for use with a specific instrument, other equivalent commercially available instruments may of course be used. Mention of a particular brand name does not constitute endorsement by the U.S. Environmental Protection Agency.

3. Applicability of this Procedure:

a. Range of Concentration:

Concentration levels as low as 0.1 mg/liter can be detected in a good flame photometer. Although the cited reference (below) does not specifically state the highest sodium concentration which may be determined, a limit of 10 mg/liter may be inferred. Therefore, sample dilution may be required.

b. Pretreatment of Samples:

The digestion procedure referenced in the Federal Register, Wednesday, December 1, 1976, Part II, table I, footnote 15 (Methods for Chemical Analysis of Water & Wastes, 1974, page 83, par. 4.1.4.) is given in this effluent monitoring procedure.

c. Treatment of Interferences in the Sample:

When the following ratios are exceeded, an interference in the determination of sodium is present, and the cited reference must be consulted.

<u>Ratio</u>	<u>Value of the Ratio</u>
potassium/sodium	5/1
calcium/sodium	10/1
magnesium/sodium	100/1

Chloride, sulfate, and bicarbonate in relatively large amounts can cause radiation interference.

This procedure was taken from the Standard Methods, 14th ed., pg. 250, Method 320 A, and from the Beckman Instrument Manual.

EFFLUENT MONITORING PROCEDURE: Determination of Sodium Using Flame Photometry

Equipment and Supply Requirements

A. Capital Equipment:

1. Beckman Model B. spectrophotometer with flame photometer attachments and instrument manual
2. Source of distilled water
3. Source of deionized water
4. Oven, for use at 140°C.
5. Analytical balance, 100 g capacity

B. Reusable Supplies:

1. Trip balance, 100 g capacity
2. Pencil or pen
3. Twelve inch ruler
4. Graduated cylinder, 50 ml (for glassware cleaning)
5. Beaker, 150 ml
6. Beaker, 250 ml (for glassware cleaning)
7. Beaker, 2 liter (for glassware cleaning)
8. Laboratory apron
9. Safety glasses
10. Glass stoppered bottle, 1 liter (for storage of cleaning solution)
11. Three polyethylene bottles, 200 ml
12. One or two polyethylene bottles, 1 liter. (See page 7-13, steps 9 and 10.)
13. Graduated cylinder, 250 ml
14. Erlenmeyer flask, 125 ml
15. Erlenmeyer flask, 500 ml
16. Graduated cylinder, 100 ml
17. Graduated cylinder, 10 ml
18. Plastic weighing boat, about 2 inches square
19. Small spatula
20. One or two volumetric flasks, 1 liter (See page 7-13, 9a.)
21. Pipet, 100 ml
22. Two plastic squeeze bottles
23. Funnel, about 60 mm diameter
24. One piece Whatman number 40 (or equivalent) filter paper (to fit the funnel)
25. One ring (to support the funnel)
26. Six, or twelve, volumetric flasks; 100 ml (see page 7-17, 3a.)
27. One pipet bulb
28. One or two graduated pipets, 10 ml (See page 7-17, 3a.)
29. Eight, or sixteen, small beakers (depending on whether one or both sets of standards are prepared); either glass (supplied with the flame photometer), or disposable plastic, and used to hold solutions for aspiration
30. Two pressure regulators (one for hydrogen, one for oxygen)
31. Wrenches for use in attaching pressure regulators to gas cylinders
32. Wrench for use in making hose connections
33. Two clamps (for safely anchoring gas cylinders)
34. Screwdriver, medium size
35. One ring stand
36. Hot plate

EFFLUENT MONITORING PROCEDURE: Determination of Sodium Using Flame Photometry

B. Reusable Supplies (Continued)

37. Thermometer, 100°C
38. One piece of rubber or plastic tubing to fit onto the tube extending from the bottom of the atomizer burner, 1 ft.
39. Asbestos gloves or crucible tongs
40. One piece of pH sensitive paper (to measure a pH of 2)

C. Consumable Supplies:

1. Sodium chloride, NaCl, 3 g.
2. One piece graph paper; divided into squares of equal size
3. Sodium dichromate, Na₂Cr₂O₇ (for cleaning glassware)
4. Concentrated sulfuric acid, H₂SO₄ (for cleaning glassware)
5. Soap (for cleaning glassware)
6. Brush (for cleaning glassware)
7. Brush (for cleaning balance)
8. Paper towels
9. Matches
10. Concentrated nitric acid, HNO₃
11. Cylinder of hydrogen gas (for use with the flame photometer)
12. Cylinder of oxygen gas (for use with the flame photometer)
13. Concentrated hydrochloric acid, HCl
14. Distilled concentrated nitric acid, HNO₃*

All reagents should be of high quality. Different chemical manufacturers may have different ways of indicating a high quality reagent. While no endorsement of one chemical manufacturer over another is intended, the following are some designations used in four chemical catalogs to indicate high quality reagents.

<u>Catalog</u>	<u>Designations</u>
Thomas	Reagent, ACS, Chemically (Pure (CP))
Matheson, Coleman & Bell	Reagent, ACS
Curtin Matheson Scientific, Inc.	Primary Standard, ACS, AR
Fisher	Certified, ACS

* This reagent is used for acidifying the sample at the time of collection. It must be of higher purity than Reagent, ACS, etc. J. T. Baker "Ultrex" nitric acid is an example of a higher purity acid.

Beckman Model B spectrophotometer
with flame photometry attachments

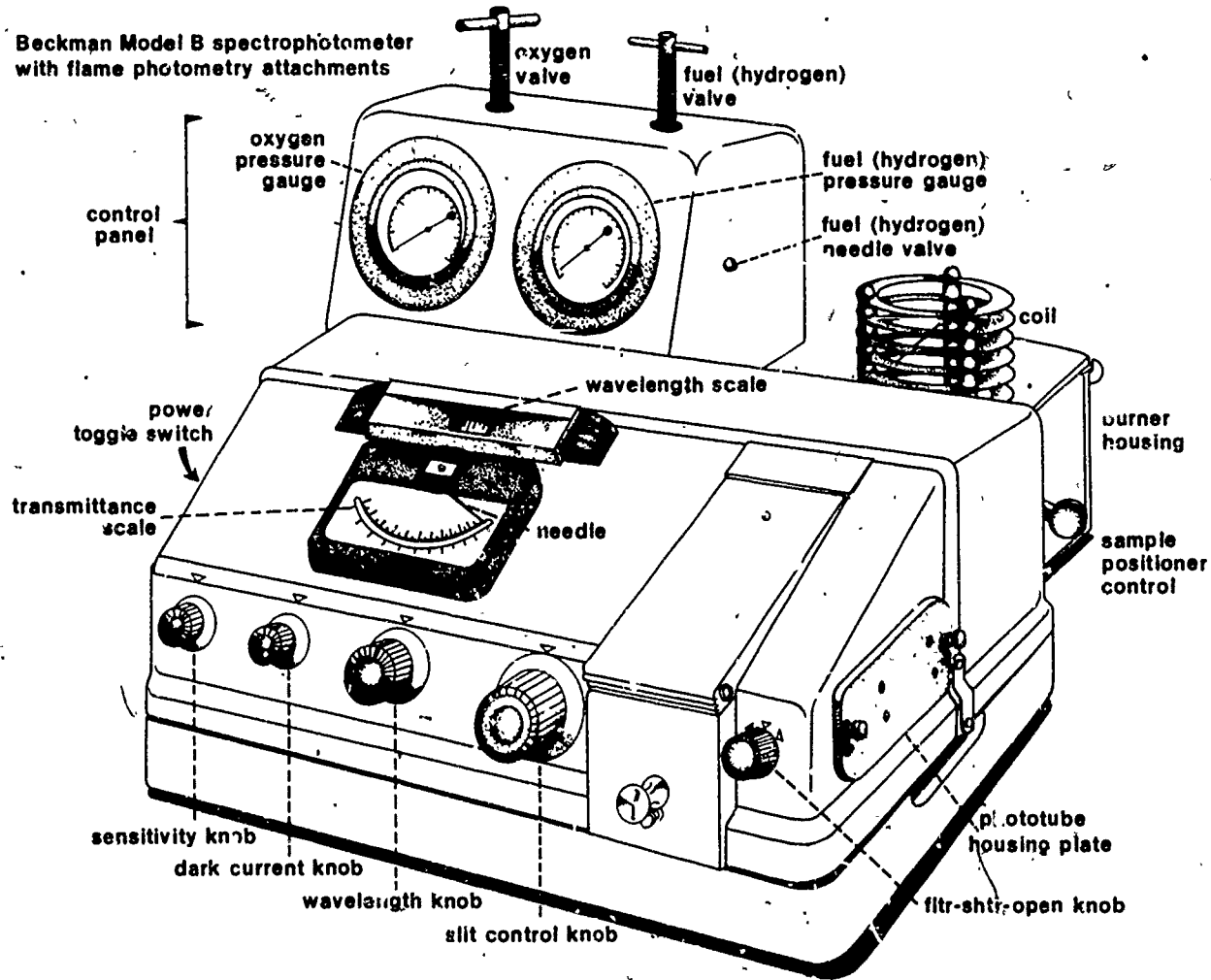


Figure 1

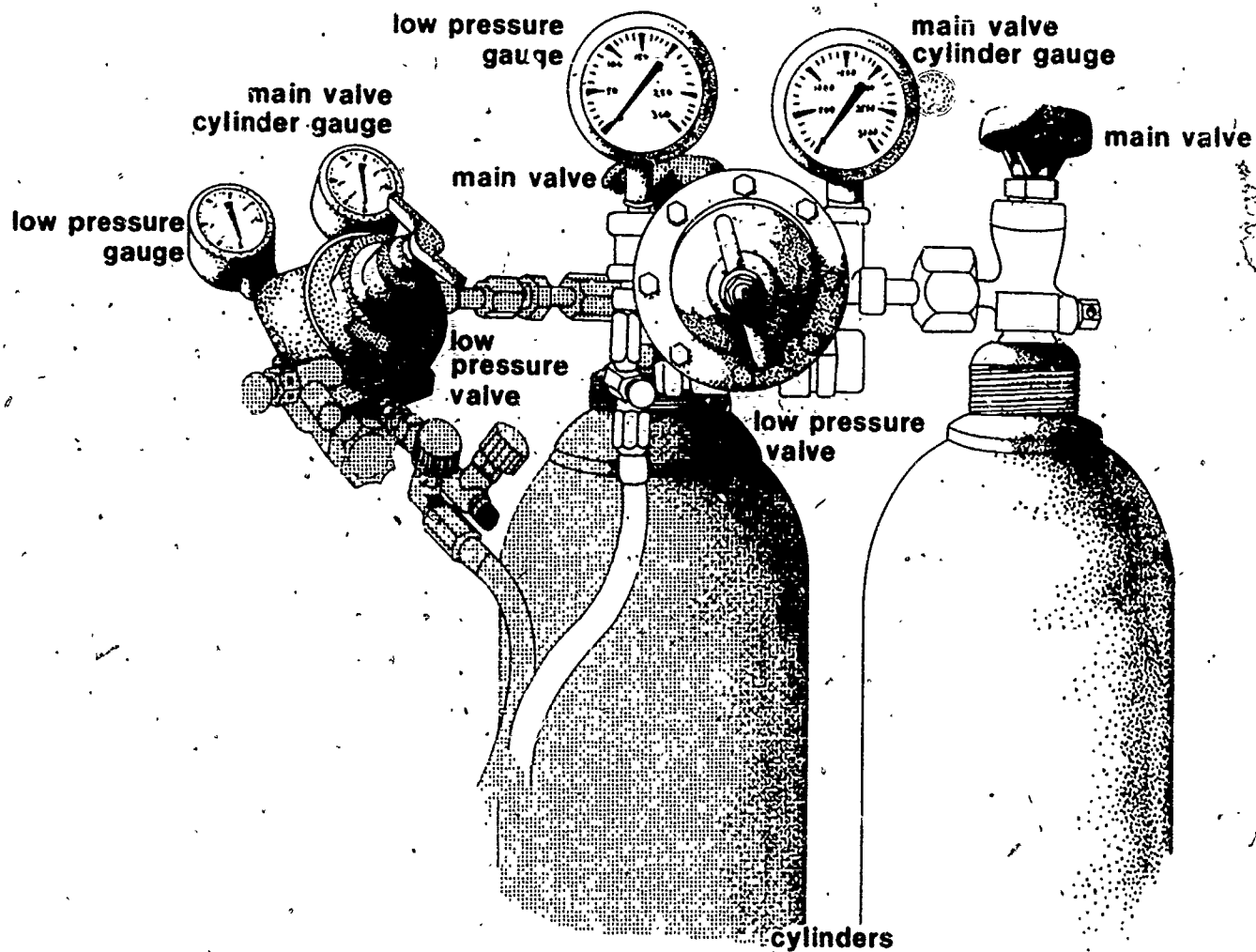


Figure 2

273

274

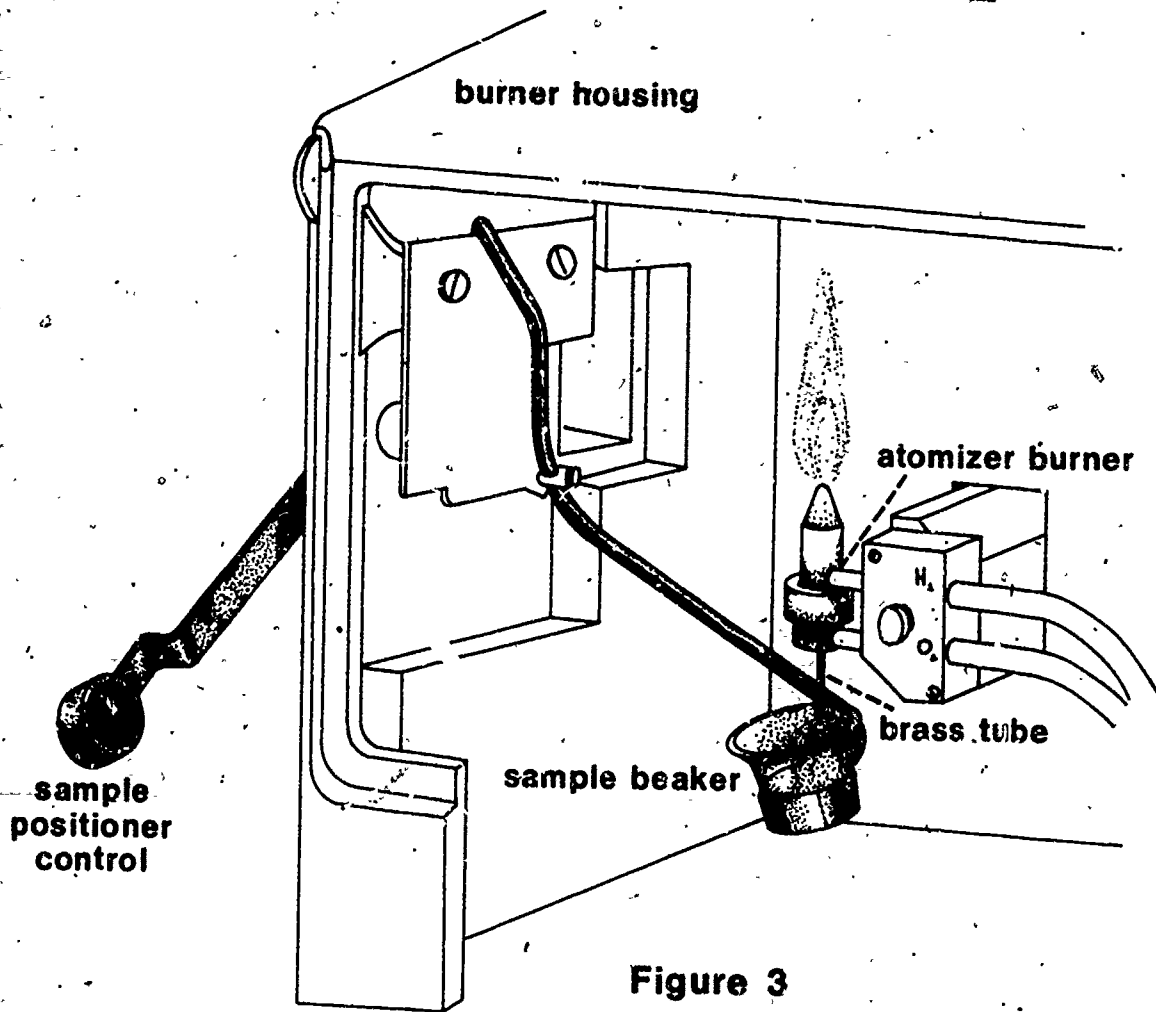


Figure 3

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
A. Equipment Preparation			
1. Cleaning of Glassware	<ol style="list-style-type: none"> 1. Clean all glassware and polyethylene bottles. 2. Rinse with nitric acid solution. 3. Rinse with tap water 4. Rinse with deionized water. 	<ol style="list-style-type: none"> 2a. Preparation of the acid is described in B.1. 3a. Ten times. 4a. Five times. 	V.A.1.1
2. Balance Preparation	<ol style="list-style-type: none"> 1. Check the analytical balance for cleanliness and proper operation. 	<ol style="list-style-type: none"> 1a. Consult the manufacturer's manual if the balance does not operate properly. 	
Spectrophotometer	<ol style="list-style-type: none"> 1. Attach the oxygen and hydrogen cylinders to the Beckman Model B spectrophotometer with flame photometry attachments. 2. Turn the wavelength knob to a reading of 580 nanometers. 3. Turn the slit control knob to a reading of 0.4 mm. 4. Turn the sensitivity knob to the standby position. 5. Turn the filter-shutter-open knob to the shutter position. 	<ol style="list-style-type: none"> 1a. Consult the manufacturer's manual for specific instructions. 1b. For the remainder of this procedure, the words instrument, or flame photometer, will be used to mean this particular instrument. 5a. This is the closed position. 	

276

277

EFFLUENT MONITORING PROCEDURE: Determination of Sodium Using Flame Photometry

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAI GUIDE NOTES
<p>A. Equipment Preparation (Continued)</p> <p>4. Phototube</p>	<ol style="list-style-type: none"> 1. Remove the phototube housing plate (see Figure 1). 2. Check the condition of the desiccant in the phototube compartment. 	<ol style="list-style-type: none"> 1a. The blue sensitive phototube should be used for sodium determination. 1b. Refer to figures 1, 2, & 3 when other parts of the instrument are mentioned in this procedure. 2a. It may need drying at 103-105°C for 1 hour, followed by cooling in a desiccator for 30 min. 2b. The spectrophotometer will be turned on in E. 	
<p>B. Reagent Preparation</p> <p>1. Nitric Acid, HNO₃, Solution, 1 + 15^v</p>	<ol style="list-style-type: none"> 1. Measure 150 ml of distilled water. 2. Pour it into a 500 ml Erlenmeyer flask. 3. Measure 10 ml of concentrated nitric acid, HNO₃. 4. Pour the nitric acid slowly into the distilled water. 5. Swirl the flask. 6. Store the solution in a polyethylene bottle. 	<ol style="list-style-type: none"> 1a. Use a 250 ml graduated cylinder. 3a. Use a 10 ml graduated cylinder. 5a. To mix the acid and water. 6a. This solution is used for rinsing glassware and polyethylene bottles. 6b. Larger amounts may be prepared using the same proportion of acid and water. 	

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
B. Reagent Preparation (continued)			
2. Nitric Acid, HNO ₃ , Distilled	1. See 1a.	1a. This reagent is used undiluted for acidifying the sample at the time of collection.	
3. Hydrochloric Acid, HCl, Distilled, 1 + 1	1. Measure 2.5 ml of distilled hydrochloric acid, HCl.	1a. Use a 10 ml graduated cylinder. 1b. Five ml of the 1 + 1 mixture are needed for each sample. Therefore, larger volumes may be prepared as necessary.	
	2. Pour it into a 125 ml Erlenmeyer flask.	2a. Use a larger flask if larger amounts of the mixture are being prepared.	
	3. Measure 2.5 ml of deionized water.	3a. Use the 10 ml graduated cylinder from 1a.	
	4. Pour it into the same flask.		
	5. Swirl the flask.	5a. To mix the acid and water.	
	6. Store the solution in a polyethylene bottle.		
4. Stock Sodium Solution	1. Weigh 3 g. of sodium chloride, NaCl.	1a. Use a trip balance. 1b. This is slightly more than is needed.	
	2. Dry it at 140°C.	2a. For 1 hour.	
	3. Cool it in a desiccator.	3a. For 30 minutes.	
	4. Weigh 2.542 g. of the sodium chloride, NaCl.	4a. Use an analytical balance.	
	5. Transfer it to a 1 liter volumetric flask.	5a. Use deionized water in the plastic squeeze bottle.	

280

281

EFFLUENT MONITORING PROCEDURE: Determination of Sodium Using Flame Photometry

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>B. Reagent Preparation (continued)</p>	<p>6. Dissolve the salt in de-ionized water in the flask.</p> <p>7. Dilute to the 1 liter mark.</p> <p>8. Mix thoroughly.</p> <p>9. If the expected sodium concentration in the sample is between 1.0 and 10.0 mg/liter, prepare the intermediate sodium solution (B.5 below).</p> <p>10. If the expected sodium concentration is between 0.1 and 1.0 mg/liter, prepare the standard sodium solution (B.6 below).</p>	<p>7a. With deionized water.</p> <p>8a. The concentration of this solution is 1.0 mg of sodium per 1.0 ml of solution.</p> <p>9a. If there is doubt about the expected sodium concentration, both the intermediate and standard sodium solutions should be prepared, along with two calibration graphs.</p>	
<p>5. Intermediate Sodium Solution</p>	<p>1. Pipet 100.0 ml of the stock sodium solution (B.4.) into a 1 liter volumetric flask</p> <p>2. Dilute to the 1 liter mark.</p> <p>3. Mix thoroughly.</p>	<p>2a. With deionized water.</p> <p>3a. The concentration of this solution is 0.1 mg of sodium per 1.0 ml of solution.</p>	

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>B. Reagent Preparation (continued)</p> <p>6. Standard Sodium Solution</p>	<ol style="list-style-type: none"> 1. Pipet 100.0 ml of the intermediate sodium solution (B.5.) into a 1 liter volumetric flask. 2. Dilute to the 1 liter mark. 3. Mix thoroughly. 4. Transfer the sodium solutions you have prepared to polyethylene bottles. 	<ol style="list-style-type: none"> 2a. With deionized water. 3a. The concentration of this solution is 0.01 mg of sodium per 1.0 ml of solution. 	
<p>C. Sample Collection</p>	<ol style="list-style-type: none"> 1. Collect the sample in a polyethylene bottle. 2. Acidify the sample with distilled nitric acid. HNO_3. 3. Gently swirl the bottle. 4. Check the pH of the sample. 	<ol style="list-style-type: none"> 1a. Although only a few ml of sample are actually needed, a larger volume, such as 200 ml, is more convenient to work with. 2a. Use 5 ml of acid/liter of sample; 1.0 ml for 200 ml of sample. 2b. Measure the acid in a 10 ml graduated cylinder. 3a. To mix the acid and sample. 4a. Use pH sensitive paper. 4b. It must be less than 2. If it is not, add a few more drops of acid, swirl the bottle, and recheck the pH. Repeat the procedure until the pH is less than 2. 4c. Do not store the sample for more than 6 months before completing the analysis. 	

EFFLUENT MONITORING PROCEDURE: Determination of Sodium Using Flame Photometry

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
D. Sample Digestion	<ol style="list-style-type: none"> 1. Shake the sample container. 2. Measure 100 ml of the acidified sample. 3. Pour it into a 250 ml Erlenmeyer flask. 4. Measure 5 ml of 1 + 1 distilled hydrochloric acid, HCl. 5. Pour it into the flask. 6. Swirl the flask. 7. Put the flask on a hot-plate. 8. Turn on the hotplate and adjust the temperature setting so the sample comes to a temperature of 95°C. 9. After the sample has come to 95°C, continue the heating for 15 min. 10. Assemble a filtration apparatus. 	<ol style="list-style-type: none"> 1a. To ensure that it is homogeneous. 2a. Use a 100 ml graduated cylinder. 3a. If any solids remain in the cylinder, rinse them into the flask with a few ml of deionized water. 4a. Use a 10 ml graduated cylinder. 6a. To mix the acid and sample. 8a. Check the temperature of the sample with a thermometer. 8b. Do not allow the temperature of the sample to go above 95°C. 9a. During the 15 minute period, proceed with Step 10. 10a. A ring stand, a funnel (approximately 60 mm diameter), a ring to hold the funnel, and a piece of medium porosity filter paper (such as Whatman number 40). 10b. Use a clean 100 ml graduated cylinder as receiver. 	

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>D. Sample Digestion (continued)</p>	<p>11. After the 15 min. heating period, remove the flask from the hotplate.</p> <p>12. Allow it to cool at room temperature for about 15 min.</p> <p>13. Filter the sample through the filter paper into the cylinder.</p> <p>14. Wash the filter paper with small portions (a few ml) of deionized water, and catch the washings in the cylinder.</p> <p>15. Cover the cylinder with an inverted 150 ml beaker.</p>	<p>11a. Be cautious of the hot flask. Use a glove or tongs.</p> <p>13a. Solids in the sample would clog the burner on the flame photometer.</p> <p>14a. When the volume in the cylinder is 100 ml, remove the cylinder from under the funnel.</p> <p>15a. Until the analysis. 15b. To prevent contamination.</p>	
<p>E. Spectrophotometer Warm-up</p>	<p>1. Plug the power cord on the instrument into a wall outlet.</p> <p>2. Turn on the power toggle switch.</p> <p>3. Allow the instrument to warm up for 45 minutes. While it is warming up, proceed with Section F.</p>	<p>1a. 115 volts, 50/60 cycle.</p> <p>2a. The indicator light will go on. 2b. Do not turn on the lamp toggle switch. It is not used when making flame photometry measurements.</p> <p>3a. With older instruments, 2 hours or longer would be a better warm-up time.</p>	

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>F. Preparation of Standards (continued)</p>	<p>7. Add deionized water to the mark in each of the six or twelve flasks.</p> <p>8. Mix thoroughly.</p>	<p>8a. The concentrations are: 0.0, 2.0, 4.0, 6.0, 8.0, and 10.0 mg/liter in the one set of standards, and 0.0, 0.2, 0.4, 0.6, 0.8, and 1.0 mg/liter in the other.</p>	
<p>G. Analysis</p> <p>1. Standards and Sample</p> <p>2. Lighting the Flame</p>	<p>1. Mark the 5 ml beakers (six or twelve) in the same way that the 100 ml volumetric flasks are marked.</p> <p>2. Pour the standards into the appropriate 5 ml beakers.</p> <p>3. Shake the sample thoroughly.</p> <p>4. Fill another 5 ml beaker with sample.</p> <p>5. Fill another 5 ml beaker with deionized water.</p> <p>1. Raise the door on the right side of the burner housing.</p>	<p>1a. Either glass or plastic disposable.</p> <p>1b. Used for aspirating solutions.</p> <p>2a. Fill the beakers about three-fourths full.</p> <p>2b. Caution: Throughout the remainder of this procedure, be careful not to get your fingers into any of the solutions. This could cause an error in the result.</p> <p>3a. Even if it was filtered earlier.</p> <p>4a. About three-fourths full.</p> <p>5a. About three-fourths full.</p>	

292

293

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>F. Preparation of Standards (continued)</p>	<p>7. Add deionized water to the mark in each of the six or twelve flasks.</p> <p>8. Mix thoroughly.</p>	<p>8a. The concentrations are: 0.0, 2.0, 4.0, 6.0, 8.0, and 10.0 mg/liter in the one set of standards, and 0.0, 0.2, 0.4, 0.6, 0.8, and 1.0 mg/liter in the other.</p>	
<p>G. Analysis</p> <p>1. Standards and Sample</p> <p>2. Lighting the Flame</p>	<p>1. Mark the 5 ml beakers (six or twelve) in the same way that the 100 ml volumetric flasks are marked.</p> <p>2. Pour the standards into the appropriate 5 ml beakers.</p> <p>3. Shake the sample thoroughly.</p> <p>4. Fill another 5 ml beaker with sample.</p> <p>5. Fill another 5 ml beaker with deionized water.</p> <p>1. Raise the door on the right side of the burner housing.</p>	<p>1a. Either glass or plastic disposable.</p> <p>1b. Used for aspirating solutions.</p> <p>2a. Fill the beakers about three-fourths full.</p> <p>2b. Caution: Throughout the remainder of this procedure, be careful not to get your fingers into any of the solutions. This could cause an error in the result.</p> <p>3a. Even if it was filtered earlier.</p> <p>4a. About three-fourths full.</p> <p>5a. About three-fourths full.</p>	

292

293

EFFLUENT MONITORING PROCEDURE: Determination of Sodium Using Flame Photometry

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>G. Analysis (continued)</p>	<ol style="list-style-type: none"> 2. Raise the sample positioner control between the instrument and the burner housing to the vertical position as far as it will go. 3. Close the oxygen and fuel (hydrogen) valves on top of the control panel. 4. Close the low pressure valve on the regulators attached to the hydrogen and oxygen cylinders. 5. Close the fuel (hydrogen) needle valve on the right side of the control panel. 6. Open the fuel (hydrogen) needle valve on the right side of the control panel. 7. Open the main valve on the oxygen cylinder. 8. Open the low pressure valve on the oxygen cylinder. 9. Set the pressure for 30 pounds per square inch. 	<ol style="list-style-type: none"> 2a. Do not force the sample positioner control. 3a. By turning them counter clockwise until they just begin to feel loose; i.e., until they can be wiggled very slightly left and right. 4a. By turning the handles counter clockwise until they just begin to feel loose; i.e., until they can be wiggled very slightly left and right. 5a. By turning it clockwise with a screwdriver as far as it will go. 5b. Do not force the screwdriver. 6a. By turning it counter clockwise with a screwdriver. 6b. About 1/8 turn. 7a. By turning the handle counter clockwise two full turns. 7b. For safety, <u>oxygen</u> is always turned on <u>first</u>, and turned off <u>last</u>. 8a. By turning the handle clockwise until the needle moves away from zero. 9a. By turning the valve clockwise. 	

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>G. Analysis (continued)</p>	<p>10. Open the oxygen valve on the control panel.</p>	<p>10a. By turning the handle clockwise until the needle moves away from zero.</p>	
	<p>11. Set the pressure for 13 pounds per square inch.</p>	<p>11a. By turning the handle clockwise. 11b. The actual pressure should be written on a tag attached to the atomizer burner. 11c. The optimum pressure may be checked if no tag is attached to the burner. 11d. Attach a small diameter piece of rubber or plastic tubing about 1 foot long to the brass tube extending from the bottom of the atomizer burner. The rubber or plastic tubing should fit the brass tube snugly. Place the other end of the rubber or plastic tubing in a 10 ml graduated cylinder. Fill it to the 10.0 line with deionized water. When the oxygen is flowing at the proper rate, the deionized water should be aspirated at the rate of 1.5 - 2.0 ml per minute. This rate corresponds to about thirteen pounds pressure per square inch.</p>	
	<p>12. Remove the tubing and cylinder.</p>		
	<p>13. Open the low pressure valve on the hydrogen cylinder.</p>	<p>13a. By turning the handle clockwise until the needle moves away from zero.</p>	
	<p>14. Set the pressure for 10 pounds per square inch.</p>	<p>14a. By turning the valve clockwise.</p>	
	<p>15. Open the fuel (hydrogen) valve on the control panel.</p>	<p>15a. By turning the handle clockwise.</p>	
	<p>16. Set the pressure for 4.5 pounds per square inch.</p>		

296

297

EFFLUENT MONITORING PROCEDURE: Determination of Sodium Using Flame Photometry

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>G. Analysis (continued)</p>	<p>17. Cautiously bring a lighted match to the tip of the atomizer burner.</p> <p>18. Close the door on the right side of the burner housing.</p> <p>19. Place the beaker containing the highest concentration standard (10 mg/liter or 1.0 mg/liter) into the holder.</p> <p>20. Swing the beaker into the burner housing compartment using the sample positioner control.</p> <p>21. Turn the sensitivity knob to position 1.</p> <p>22. Make the needle read 0% transmittance using the dark current knob.</p>	<p>17a. The oxygen-hydrogen mixture will ignite with a loud pop and hissing sound.</p> <p>18a. Caution: hot gases are escaping from the top of the coil. Do not place your hand or any other part of your body over the coil while the gases are burning.</p> <p>19a. See figure 3.</p> <p>19b. If two sets of standards are being used, take readings on the 10, 8, 6, 4, 2, and 0 mg/liter solutions <u>first</u>, and <u>then</u> the 1.0, 0.8, 0.6, 0.4, 0.2, and 0.0 mg/liter solutions. Do not <u>mix</u> the two sets of standards.</p> <p>20a. A distinct change in sound will be heard as the solution is aspirated into the burner.</p> <p>20b. For the remainder of this procedure, if the sound caused by the aspiration of a solution changes <u>while the solution is being aspirated</u>, it probably means that all of the liquid in the beaker has been aspirated, and the beaker is empty.</p> <p>20c. Swing the beaker out of the burner housing, refill it with the appropriate solution, and swing it back into the burner housing.</p> <p>20d. While the solution is aspirating, check that the hydrogen pressure is still 10 pounds per square inch and that the oxygen pressure is still at the optimum valve determined above. If one or both of the pressures have changed, reset them.</p> <p>22a. In this procedure, all readings will be made on the transmittance portion of the scale, not the absorbance portion.</p>	



OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>G. Analysis (continued)</p> <p>3. Aspiration of Standards and Sample</p> <p>300</p>	<p>23. Turn the sensitivity knob to position 4.</p> <p>24. Repeat step 22.</p> <p>25. Adjust the fuel needle valve so the tip of the flame comes to the top of the burner housing compartment, not to the top of the coil on top of the compartment.</p> <p>1. Turn the filter/shutter knob to the open position.</p> <p>2. Turn the wavelength knob very slowly clockwise.</p> <p>3. Record this wavelength and proceed to step 14.</p>	<p>25a. Use a screwdriver.</p> <p>25b. Caution: extremely hot gases are escaping through the coil.</p> <p>25c. Adjustment of this needle valve setting is not necessary each time the instrument is used, unless the setting has been changed for some reason.</p> <p>1a. The shutter is open in this position.</p> <p>1b. In the shutter position, the shutter is closed.</p> <p>2a. Turn the knob continuously. Do not turn the knob in steps.</p> <p>2b. One of three things will happen, 2c., 2d., or 2e. <u>2c. is the condition you are trying to achieve.</u></p> <p>2c. First: the needle will move to the left, and at some % transmittance between 50 and 100, will suddenly swing back to the right. In this case, proceed to step 3.</p> <p>2d. Second: the needle will swing back to the right as in 2c, but at some % transmittance between 0 and 50. In this case, proceed to step 4.</p> <p>2e. Thirdly: the needle will swing back to the right, but at some point greater than 100% transmittance; or, it will not swing back to the right, it will stay to the left. In this case, proceed to step 9.</p> <p>3a. Use the example data sheet(s); see pages 29 & 30.</p> <p>3b. Do not change this wavelength until the procedure instructs you to change it.</p>	<p>301</p>

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
G. Analysis (continued)		3c. This wavelength should be 589 nanometers. If it is too far away from this value (less than 584, or more than 594), the calibration of the instrument should be checked using the instrument manual.	
	<p>4. Close the shutter.</p> <p>5. Increase the reading on the slit knob by 0.05 millimeter.</p> <p>6. Turn the wavelength back to 580 nanometers.</p> <p>7. Repeat steps 1 and 2.</p> <p>8. Record this wavelength and proceed to step 14.</p> <p>9. Close the shutter.</p> <p>10. Decrease the reading on the slit knob by 0.05 millimeter.</p> <p>11. Turn the wavelength back to 580 nanometers.</p> <p>12. Repeat steps 1 and 2.</p>	<p>5a. It is graduated in tenths of a millimeter between 0.0 and 1.5 millimeters.</p> <p>7a. If condition 2d still exists, repeat steps 4 through 7 until the needle moves as described in 2c.</p> <p>8a. Use the example data sheet(s); see pages 29 & 30.</p> <p>8b. Do not change this wavelength until the procedure instructs you to change it.</p> <p>10a. It is graduated in tenths of a millimeter between 0.0 and 1.5 millimeters.</p> <p>12a. If condition 2e still exists, repeat steps 9 through 11 until the needle moves as described in 2c.</p>	

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>G. Analysis (continued)</p>	<p>13. Record this wavelength and proceed to step 14.</p> <p>14. Turn the slit knob so as to make the needle read 100% transmittance.</p> <p>15. Close the shutter.</p> <p>16. Record a reading of 100% transmittance for the 10 or 1.0 mg/liter solution.</p> <p>17. Swing the beaker out of the burner housing compartment.</p> <p>18. Replace it with the 8 or 0.8 mg/liter solution.</p> <p>19. Swing it into the burner housing compartment.</p> <p>20. Open the shutter.</p> <p>21. Record the % transmittance reading.</p> <p>22. Close the shutter.</p>	<p>13a. Use the example data sheet(s); see pages 29 & 30.</p> <p>13b. Do not change this wavelength until the procedure instructs you to change it.</p> <p>15a. Do not change this setting until the procedure instructs you to change it.</p> <p>16a. Use the example data sheet(s); see pages 29 & 30.</p> <p>16b. This first set of readings you will get on the one or two sets of standards and sample are called peak readings.</p> <p>21a. Use the example data sheet(s); see pages 29 & 30.</p>	

304

305

EFFLUENT MONITORING PROCEDURE: Determination of Sodium Using Flame Photometry

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
G: Analysis (continued)	23. Repeat steps 17 through 22 for the 6, 4, 2, and 0 mg/liter solutions and the sample, or the 0.6, 0.4, 0.2, and 0.0 mg/liter solutions and the sample.	23a. Use the example data sheet(s); see pages 29 & 30.	
	<p>24. Swing the beaker of sample out of the burner housing compartment,</p> <p>25. Replace it with the beaker of deionized water.</p> <p>26. Swing it into the burner housing compartment.</p> <p>27. Turn the dark current knob so the needle reads 0% transmittance.</p> <p>28. Replace the beaker of deionized water with the beaker of the 10 or 1.0 mg/liter solution.</p> <p>29. Open the shutter.</p> <p>30. Turn the wavelength knob very slowly clockwise.</p>	<p>28a. It may need refilling.</p> <p>29a. The needle should show 100% transmittance; it may be "off" by plus or minus 1 or 2% because of the dark current adjustment. Do not change the slit setting.</p> <p>30a. Turn the knob continuously. Do not turn the knob, stop, and start again.</p> <p>30b. The needle will move to the right.</p> <p>30c. At some low % transmittance value, perhaps less than 1, or even 0, the needle will move no farther to the right.</p>	

306

307

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>G. Analysis (continued)</p>	<p>30d. Record this wavelength.</p> <p>31. Record this low % transmittance for the 10 of 1.0 mg/liter solution.</p> <p>32. Close the shutter.</p> <p>33. Swing the beaker out of the burner housing compartment.</p> <p>34. Replace it with the 8 or 0.8 mg/liter solution.</p> <p>35. Swing it into the burner housing.</p> <p>36. Open the shutter.</p> <p>37. Record the % transmittance reading.</p> <p>38. Close the shutter.</p> <p>39. Repeat steps 34 through 39 to obtain readings for the 6, 4, 2, and 0 mg/liter solutions and the sample, or the 0.6, 0.4, 0.2, and 0.0 mg/liter solutions, and the sample.</p>	<p>30d. If there is some doubt as to when the needle stops moving to the right, turn the wavelength knob about one-quarter turn, counter clockwise and repeat step 30.</p> <p>31a. Use the example data sheet(s); see pages 29 & 30.</p> <p>31b. It is sometimes called the background wavelength.</p> <p>31c. Do not change this wavelength setting for the remainder of the procedure.</p> <p>32a. Use the example data sheet(s); see pages 29 & 30. This second set of readings you will get on the one or two sets of standards and sample are called background readings.</p> <p>38a. Use the example data sheet(s); see pages 29 & 30.</p>	

308

309

EFFLUENT MONITORING PROCEDURE: Determination of Sodium Using Flame Photometry

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
<p>G. Analysis (continued)</p>	<p>41. Replace the beaker of sample with the beaker of deionized water.</p> <p>42. Allow the deionized water to aspirate for 10 seconds.</p> <p>43. Swing the beaker of deionized water out of the burner housing.</p> <p>44. Close the main valve on the hydrogen cylinder.</p> <p>45. When the hydrogen pressure registers 0 on the control panel hydrogen (fuel) gauge, close the main valve on the oxygen cylinder.</p> <p>46. Turn the sensitivity knob to the standby position.</p> <p>47. Turn off the power toggle switch.</p> <p>48. Empty all of the small beakers used.</p> <p>49. Rinse the beakers thoroughly with tap water.</p> <p>50. Rinse the beakers thoroughly with deionized water.</p>	<p>44a. By turning clockwise all the way.</p> <p>45a. By turning clockwise all the way.</p> <p>48a. Don't forget the one still in the instrument.</p> <p>48b. If the beakers are of the disposable type, discard them and proceed to H. If they are glass, do steps 49, 50, and 51.</p>	

OPERATING PROCEDURES	STEP SEQUENCE	INFORMATION/OPERATING GOALS/SPECIFICATIONS	TRAINING GUIDE NOTES
G. Analysis (continued)	51. Allow the beakers to drain dry, and then proceed to Section H.		
H. Calculations	<ol style="list-style-type: none"> 1. For all of the standards, subtract the % transmittance obtained at the background wavelength from the % transmittance obtained at the peak wavelength. 2. On "regular" graph paper prepare a calibration graph using the % transmittance differences along the vertical axis, and the corresponding concentrations (10, 8, 6, 4, 2, 0, or 1.0, 0.8, 0.6, 0.4, 0.2, 0.0 mg/liter) on the horizontal axis. 3. Repeat step 1 for the sample. 4. Determine the sodium concentration in the sample. 5. Record the result in mg/liter. 	<ol style="list-style-type: none"> 1a. Example: 10.0 mg/liter solution <ul style="list-style-type: none"> 100 = % transmittance at the peak wavelength 1 = % transmittance at the background wavelength 99 = % transmittance difference 2a. If you have used both sets of standards (the intermediate and standard sodium solutions), prepare two calibration graphs. Pages 31 and 32 are sheets of "regular" graph paper. 2b. "Regular" graph paper is marked off in squares of equal size. 2c. EMP CH.IN.cg.EMP.1a.1.77 may be referred to for specific instructions in calibration graph preparation and use. 2d. In the case of the intermediate sodium solution, the calibration graph may have a slight bend. 4a. Using the calibration graph. 5a. Use the example data sheet(s); see pages 29 & 30. 	

312

313

EFFLUENT MONITORING PROCEDURE: Determination of Sodium Using Flame Photometry

**EXAMPLE
DATA SHEET**

<u>Concentration of Sodium Standards in mg/liter</u>	<u>% Transmittance at the Peak Wavelength; Peak Readings</u>	<u>% Transmittance at the Background Wavelength; Background Readings</u>	<u>% Transmittance Differences; Values in 2nd Column Minus Values in 3rd Col.</u>
_____*	_____	_____	_____
_____	_____	_____	_____
_____	_____	_____	_____
_____	_____	_____	_____
_____	_____	_____	_____
_____	_____	_____	_____

Sample

Concentration of sodium in the sample in mg/liter _____

Peak Wavelength _____ nm

Background Wavelength _____ nm

* The six values in this column will be 0, 2, 4, 6, 8, and 10, or, 0.0, 0.2, 0.4, 0.6, 0.8, and 1.0. If both sets of standards were used, use a second data sheet exactly like this one; see page 30.



EFFLUENT MONITORING PROCEDURE: Determination of Sodium Using Flame Photometry

EXAMPLE DATA SHEET
 (See the paragraph at the bottom of Page 29)

<u>Concentration of Sodium Standards in mg/liter</u>	<u>%Transmittance at the Peak Wavelength; Peak Readings</u>	<u>%Transmittance at the Background Wavelength; Background Readings</u>	<u>%Transmittance Differences; Values in 2nd Column Minus Values in 3rd Col.</u>
--	---	---	--

_____	_____	_____	_____
_____	_____	_____	_____
_____	_____	_____	_____
_____	_____	_____	_____
_____	_____	_____	_____
_____	_____	_____	_____
_____	_____	_____	_____
_____	_____	_____	_____


Sample


Concentration of sodium in the sample in mg/liter _____

Peak Wavelength _____ nm

Background Wavelength _____ nm



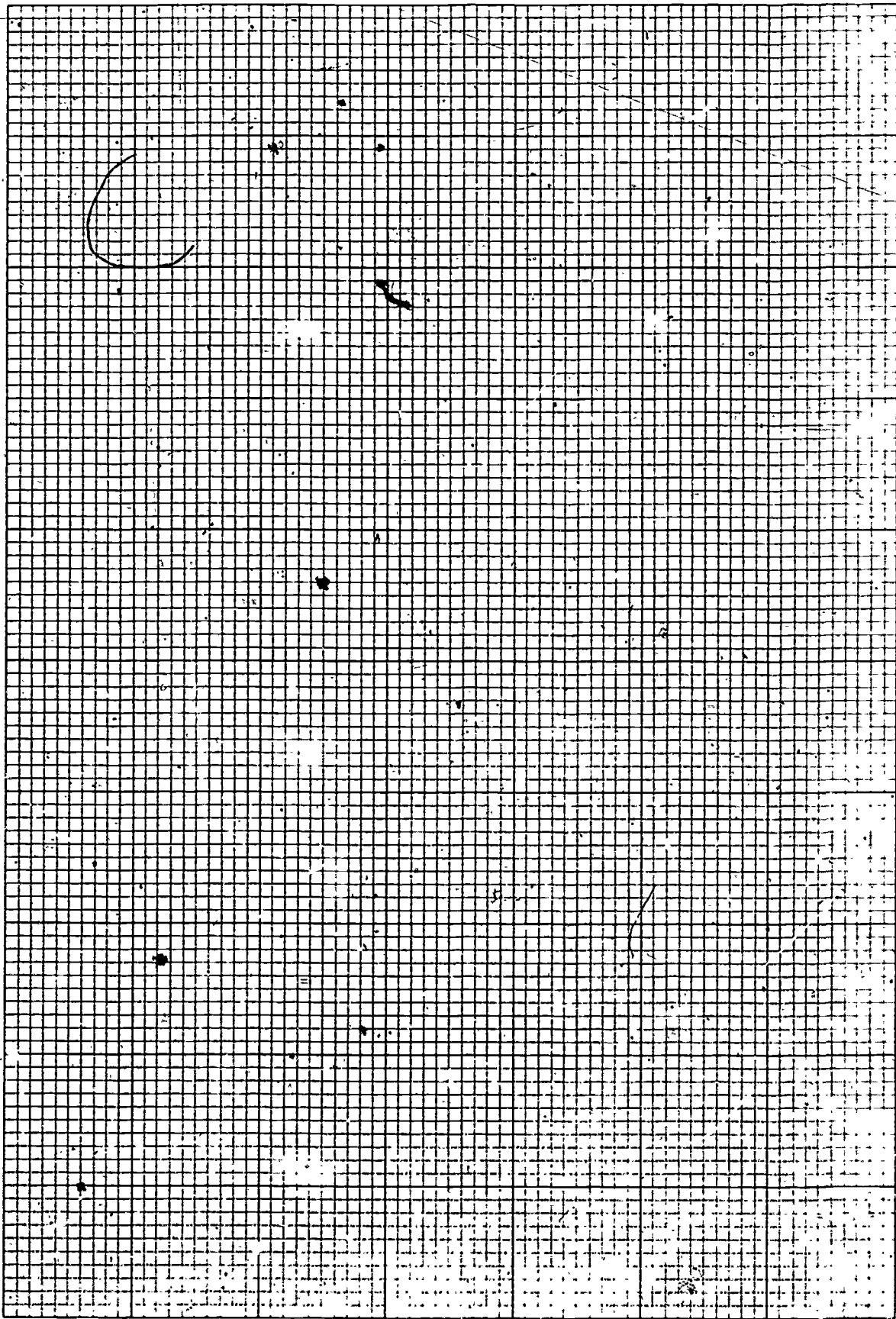

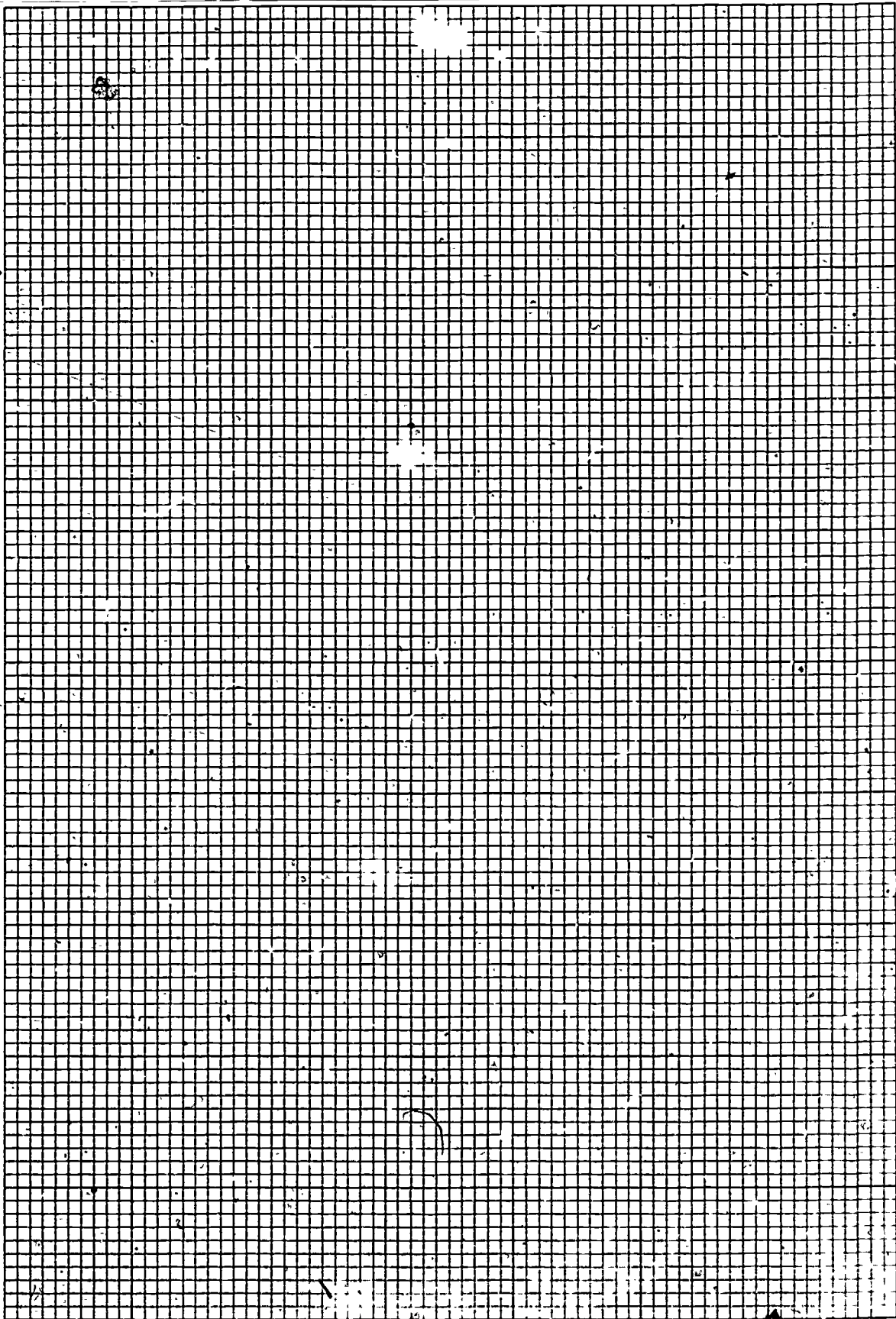

EFFLUENT MONITORING PROCEDURE: Determination of Sodium Using Flame Photometry

TRAINING GUIDE

<u>SECTION</u>	<u>TOPIC</u>
I	Introduction
II	Educational Concepts - Mathematics
III	Educational Concepts - Science
IV	Educational Concepts - Communications
V*	Field and Laboratory Equipment
VI	Field and Laboratory Reagents
VII	Field and Laboratory Analysis
VIII	Safety
IX	Records & Reports

* Training guide materials are presented here under the headings marked *.

FIELD AND LABORATORY EQUIPMENT

Section V

TRAINING GUIDE NOTE

REFERENCES/RESOURCES

A.1.1

If the glassware is especially dirty and cannot be cleaned with ordinary detergents, chromic acid cleaning may be required.

1. Pour 35 ml of distilled water in a 250 ml beaker.
 2. Add about 1/8 teaspoon (simply estimate this quantity) of sodium dichromate, $\text{Na}_2\text{Cr}_2\text{O}_7$, to the water.
 3. Swirl the beaker until the sodium dichromate has dissolved.
 4. Keep repeating steps 2 and 3 until no more sodium dichromate will dissolve.
 5. Pour the solution into a 2 liter beaker.
 6. Slowly pour 1 liter of concentrated sulfuric acid, H_2SO_4 , into the 2 liter beaker.
- Caution: Use eyeglasses and protective clothing.
7. Stir the mixture thoroughly.
 8. Store it in a glass stoppered bottle.
 9. The cleaning solution should be at a temperature of about 50°C when it is used.
 10. It may therefore be necessary to warm the cleaning solution.
 11. When using the warm cleaning solution, fill the piece of glassware with the solution.
 12. Allow it to soak for 2-3 minutes (or longer).
 13. Pour the cleaning solution back into the storage bottle.
 14. Rinse the piece of glassware ten times with tap water.
 15. The cleaning solution may be reused until it turns green.
 16. It should then be discarded.

Standard Methods, 14th ed., P. 356, section 2c.2)